



# SampleTrack<sup>TM</sup>

# SampleTrack Manual User Guide

Version 002



The information in this manual may be altered without notice.

BRUKER BIOSPIN accepts no responsibility for actions taken as a result of use of this manual. BRUKER BIOSPIN accepts no liability for any mistakes contained in the manual, leading to coincidental damage, whether during installation or operation of the instrument. Unauthorized reproduction of manual contents, without written permission from the publishers, or translation into another language, either in full or in part, is forbidden.

This manual was written by

#### Gilbert Cyppel / Stanley Niles

Special thanks to Birk Schütz who developed the Web interface.

© August 12, 2005: Bruker Biospin GmbH

Rheinstetten, Germany

P/N: Z31731 DWG-Nr.: 1438002

# Contents

	Contents	3
1	Copyright Notices	. <b>. 9</b> 9
2	Warnings and Notes	-
<b>-</b> 2.1	Warnings and Notes	
2.1 2.2	Contact for Additional Technical Assistance	
3	Introduction to Bruker LIMS	13
3.1	Laboratory Data Management Today	13
3.2	Bruker Laboratory Information and Management System	
3.3	The SampleTrack™ Interfaces	
3.4	SampleTrack™ as Data Reference	
3.5	Daily Routine Laboratory Functions	
3.6 3.7	Communication Overview SampleTrack™ Workflow Architecture	
3.7 3.8	Integrative Bioinformatics Solution - A Look into the Future	
4	Introduction to the SampleTrack™ Client?	21
4.1	SampleTrack™ Login	21
4.2	SampleTrack™ Overview	22
4.3	Sample Orders Module	
4.4	Details of a Preparation	
4.5	Sample Order Experiment Details	
4.6	Details Window Buttons	
4.7 4.8	Sample Order Window Buttons Sample Order Status Information	
4.8	Filters & Views	
1.0	The Right to Define Filters	
5	Functions of the SampleTrack™ Client Module	33
5.1	Submit Sample Orders	33
5.2	Manual Preparation	35
5.3	Proton Exchange	37
5.4	Finish Manual Experiments	
5.5	Delay Experiments	
5.6	Experiments Waiting (Delayed)	
5.7	Experiments Waiting for HD Exchange	
5.8	Order Approval	
5.9 5.10	E-mail Notification	
5.10	Environment Setup	43

6	Working with the SampleTrack™ Client45
6.1	Sample Order Details Window45
6.2	Entering the Sample and Client ID 45
6.3	Define a Method and Enter the Journal Number
6.4	Select the Preparation Workstep 47
6.5	Adding Experiments 48
6.6	Adding D20 for Proton Exchange 49
6.7	Copy an Existing Experiment 49
6.8	Starting a Preparation and Measurement 51
	Standard Method51
	Manual Submit in the Sample Order Details Window51
6.9	Rerunning a Faulty or Finished Experiment 52
6.10	Adding an Experiment to a Finished Sample Order 53
6.11	Filter and Search Tools
	Filter Tools
	Search Tools55
7	Interaction with Analytical Applications57
7.1	Interaction with TOPSPIN®57
7.2	Interaction with AMIX™58
7.3	Interaction with Structure Editors 59
7.4	Interaction with other Windows® Applications 59
7.5	Creating a New Report Within the Details Window
7.6	Application Management 61
	Application Type62
	Selecting Applications in the Sample Order Window 63
	Selecting Applications in the Detail Window
8	EasyDialog65
8.1	Introduction to EasyDialog
8.2	Opening EasyDialog
8.3	EasyDialog Login Window
8.4	The EasyDialog Window
8.5	Entering Data in the EasyDialog Window
8.6	EasyDialog Confirmation Message67
9	SampleTrack™ Web Client69
	-
9.1	WEB Client - Introduction
9.2	Viewing Sample Orders
	Sample Orders – View
	Structure and Web Spectra Viewer
	Search
	Extended Search74 Instrument Order View75
	Matrix Orders
9.3	Enter Orders Using the EasyWeb Dialog
9.3 9.4	Applications (only for Windows® Systems)
JT	
10	Interaction with Automation Robots

10.1	SampleRail™	
10.2	TECAN® - Pooling Module	79
	TECAN® Requirements	80
	TECAN® Plugin Installation	80
	TECAN® Matrix Settings	
	TECAN® Pooling Settings	
	TECAN® Instrument Setup	
10.3	Protein Screening / Module Pooling	
10.5		
	TECAN® Generating Samples	
	TECAN® Starting the Automation Process	83
11	Environment Setup	85
11.1	SampleTrack™ Environment Setup	85
11.2	Inserting a New User in the Companies Setup Module .	
11.3	Inserting a New Company	
11.3	• · ·	
	Entering a New Department	
11.5	Entering a New User	
11.6	Deleting a User, Department, or Company	
11.7	Inserting a New Instrument	
11.8	Inserting a New Method Type	
11.9	Inserting a New Solvent	90
11.10	Inserting a Proton Exchange Solvent	91
11.11	Inserting a New Preparation Method	92
11.12	Inserting a Solvent Mixing Method	93
11.13	Inserting New Experiments	94
11.14		
11.14	Experiment Substitution for Imported Files	
11.14	Experiment Substitution for Imported Files	95
<b>12</b>	The Object Manager	
	The Object Manager	97
12	The Object Manager	<b> 97</b> 97
<b>12</b> 12.1 12.2	The Object Manager Introduction to the Object Manager Opening the Object Manager	<b>97</b> 97 98
<b>12</b> 12.1 12.2 12.3	The Object Manager Introduction to the Object Manager Opening the Object Manager The Object Manager Main Menu	97 97 98 99
<b>12</b> 12.1 12.2 12.3 12.4	The Object Manager Introduction to the Object Manager Opening the Object Manager The Object Manager Main Menu The Group Tree and Content Sub-windows	97 97 98 99 100
<b>12</b> 12.1 12.2 12.3 12.4 12.5	The Object Manager Introduction to the Object Manager Opening the Object Manager The Object Manager Main Menu The Group Tree and Content Sub-windows Viewing Group Details in the Content Window	97 97 98 99 100 101
<b>12</b> 12.1 12.2 12.3 12.4 12.5 12.6	The Object Manager         Introduction to the Object Manager         Opening the Object Manager         The Object Manager Main Menu         The Group Tree and Content Sub-windows         Viewing Group Details in the Content Window         Client Profile Example	97 97 98 99 100 101 101
<b>12</b> 12.1 12.2 12.3 12.4 12.5 12.6 12.7	The Object Manager         Introduction to the Object Manager         Opening the Object Manager         The Object Manager Main Menu         The Group Tree and Content Sub-windows         Viewing Group Details in the Content Window         Client Profile Example         User Classification and Client Profiles	97 97 98 99 100 101 101 102
<b>12</b> 12.1 12.2 12.3 12.4 12.5 12.6	The Object Manager         Introduction to the Object Manager         Opening the Object Manager         The Object Manager Main Menu         The Group Tree and Content Sub-windows         Viewing Group Details in the Content Window         Client Profile Example	97 97 98 99 100 101 101 102
<b>12</b> 12.1 12.2 12.3 12.4 12.5 12.6 12.7	The Object Manager         Introduction to the Object Manager         Opening the Object Manager         The Object Manager Main Menu         The Group Tree and Content Sub-windows         Viewing Group Details in the Content Window         Client Profile Example         User Classification and Client Profiles	97 97 98 99 100 101 101 102 103
<b>12</b> 12.1 12.2 12.3 12.4 12.5 12.6 12.7 12.8 <b>13</b>	The Object Manager         Introduction to the Object Manager         Opening the Object Manager         The Object Manager Main Menu         The Group Tree and Content Sub-windows         Viewing Group Details in the Content Window         Client Profile Example         User Classification and Client Profiles         Adding a User to an Existing Group	97 97 98 99 100 101 101 102 103
<b>12</b> 12.1 12.2 12.3 12.4 12.5 12.6 12.7 12.8 <b>13</b>	The Object Manager         Introduction to the Object Manager         Opening the Object Manager         The Object Manager Main Menu         The Group Tree and Content Sub-windows         Viewing Group Details in the Content Window         Client Profile Example         User Classification and Client Profiles         Adding a User to an Existing Group         Example: Create a User Group.         Creating a New Group	97 97 98 99 100 101 101 103 105
<b>12</b> 12.1 12.2 12.3 12.4 12.5 12.6 12.7 12.8 <b>13</b> 13.1 13.2	The Object Manager         Introduction to the Object Manager         Opening the Object Manager         The Object Manager Main Menu         The Group Tree and Content Sub-windows         Viewing Group Details in the Content Window         Client Profile Example         User Classification and Client Profiles         Adding a User to an Existing Group         Creating a New Group         Adding a Client Profile	97 97 98 99 100 101 101 102 103 105 105
<ul> <li>12.1</li> <li>12.2</li> <li>12.3</li> <li>12.4</li> <li>12.5</li> <li>12.6</li> <li>12.7</li> <li>12.8</li> <li>13.1</li> <li>13.2</li> <li>13.3</li> </ul>	The Object Manager         Introduction to the Object Manager         Opening the Object Manager         The Object Manager Main Menu         The Group Tree and Content Sub-windows         Viewing Group Details in the Content Window         Client Profile Example         User Classification and Client Profiles         Adding a User to an Existing Group         Creating a New Group         Adding a Client Profile         Adding Table Defaults to the New Group	
<ul> <li>12.1</li> <li>12.2</li> <li>12.3</li> <li>12.4</li> <li>12.5</li> <li>12.6</li> <li>12.7</li> <li>12.8</li> <li>13.1</li> <li>13.2</li> <li>13.3</li> <li>13.4</li> </ul>	The Object Manager         Introduction to the Object Manager         Opening the Object Manager         The Object Manager Main Menu         The Group Tree and Content Sub-windows         Viewing Group Details in the Content Window         Client Profile Example         User Classification and Client Profiles         Adding a User to an Existing Group         Creating a New Group         Adding a Client Profile         Adding a New Group         Inserting a New Group EasyDialog Configuration	97 97 98 99 100 101 101 102 103 105 105 105 106 108
<ul> <li>12.1</li> <li>12.2</li> <li>12.3</li> <li>12.4</li> <li>12.5</li> <li>12.6</li> <li>12.7</li> <li>12.8</li> <li>13.1</li> <li>13.2</li> <li>13.3</li> <li>13.4</li> <li>13.5</li> </ul>	The Object Manager         Introduction to the Object Manager         Opening the Object Manager         The Object Manager Main Menu         The Group Tree and Content Sub-windows         Viewing Group Details in the Content Window         Client Profile Example         User Classification and Client Profiles         Adding a User to an Existing Group         Creating a New Group         Adding Table Defaults to the New Group         Inserting a New Filter Method	
<ul> <li>12.1</li> <li>12.2</li> <li>12.3</li> <li>12.4</li> <li>12.5</li> <li>12.6</li> <li>12.7</li> <li>12.8</li> <li>13.1</li> <li>13.2</li> <li>13.3</li> <li>13.4</li> </ul>	The Object Manager         Introduction to the Object Manager         Opening the Object Manager         The Object Manager Main Menu         The Group Tree and Content Sub-windows         Viewing Group Details in the Content Window         Client Profile Example         User Classification and Client Profiles         Adding a User to an Existing Group         Creating a New Group         Adding a Client Profile         Adding a New Group         Inserting a New Group EasyDialog Configuration	
<ul> <li>12.1</li> <li>12.2</li> <li>12.3</li> <li>12.4</li> <li>12.5</li> <li>12.6</li> <li>12.7</li> <li>12.8</li> <li>13.1</li> <li>13.2</li> <li>13.3</li> <li>13.4</li> <li>13.5</li> </ul>	The Object Manager         Introduction to the Object Manager         Opening the Object Manager         The Object Manager Main Menu         The Group Tree and Content Sub-windows         Viewing Group Details in the Content Window         Client Profile Example         User Classification and Client Profiles         Adding a User to an Existing Group         Creating a New Group         Adding Table Defaults to the New Group         Inserting a New Filter Method	97 97 98 99 100 101 101 102 103 103 105 105 105 105 106 108 109 109
<ul> <li>12.1</li> <li>12.2</li> <li>12.3</li> <li>12.4</li> <li>12.5</li> <li>12.6</li> <li>12.7</li> <li>12.8</li> <li>13.1</li> <li>13.2</li> <li>13.3</li> <li>13.4</li> <li>13.5</li> <li>13.6</li> </ul>	The Object Manager         Introduction to the Object Manager         Opening the Object Manager         The Object Manager Main Menu         The Object Manager Main Menu         The Group Tree and Content Sub-windows         Viewing Group Details in the Content Window         Client Profile Example         User Classification and Client Profiles         Adding a User to an Existing Group         Creating a New Group         Adding a Client Profile         Adding Table Defaults to the New Group         Inserting a New Group EasyDialog Configuration         Inserting a New Group Name	
<ul> <li>12.1</li> <li>12.2</li> <li>12.3</li> <li>12.4</li> <li>12.5</li> <li>12.6</li> <li>12.7</li> <li>12.8</li> <li>13.1</li> <li>13.2</li> <li>13.3</li> <li>13.4</li> <li>13.5</li> <li>13.6</li> <li>14</li> </ul>	The Object Manager         Introduction to the Object Manager         Opening the Object Manager         The Object Manager Main Menu         The Object Manager Main Menu         The Group Tree and Content Sub-windows         Viewing Group Details in the Content Window         Client Profile Example         User Classification and Client Profiles         Adding a User to an Existing Group         Creating a New Group         Adding Table Defaults to the New Group         Inserting a New Group EasyDialog Configuration         Inserting a New Group Name         Handling the Tube ID         The Tube ID and TubeID.txt File	
<ul> <li>12.1</li> <li>12.2</li> <li>12.3</li> <li>12.4</li> <li>12.5</li> <li>12.6</li> <li>12.7</li> <li>12.8</li> <li>13.1</li> <li>13.2</li> <li>13.3</li> <li>13.4</li> <li>13.5</li> <li>13.6</li> <li>14.1</li> <li>14.2</li> </ul>	The Object Manager         Introduction to the Object Manager         Opening the Object Manager         The Object Manager Main Menu         The Group Tree and Content Sub-windows         Viewing Group Details in the Content Window         Client Profile Example         User Classification and Client Profiles         Adding a User to an Existing Group         Creating a New Group         Adding a Client Profile         Adding a Client Profile         Adding a New Group         Inserting a New Group EasyDialog Configuration         Inserting a New Group Name         Inserting a New Group Name         Inserting a New Group EasyDialog Configuration         Inserting a New Group Name         Inserting a New Group Name	
<ul> <li>12.1</li> <li>12.2</li> <li>12.3</li> <li>12.4</li> <li>12.5</li> <li>12.6</li> <li>12.7</li> <li>12.8</li> <li>13.1</li> <li>13.2</li> <li>13.3</li> <li>13.4</li> <li>13.5</li> <li>13.6</li> <li>14.1</li> </ul>	The Object Manager         Introduction to the Object Manager         Opening the Object Manager         The Object Manager Main Menu         The Object Manager Main Menu         The Group Tree and Content Sub-windows         Viewing Group Details in the Content Window         Client Profile Example         User Classification and Client Profiles         Adding a User to an Existing Group         Creating a New Group         Adding Table Defaults to the New Group         Inserting a New Group EasyDialog Configuration         Inserting a New Group Name         Handling the Tube ID         The Tube ID and TubeID.txt File	

15	System Parameters	. 113
15.1	SampleTrack™ Client Parameters	113
15.2	EasyDialog Parameters	
16	Shortcuts in SampleTrack™	. 115
17	SampleTrack™ Tables	. 117
17.1	SampleTrack™ Base Tables	117
	Introduction to SampleTrack Base Tables	. 117
	The EasyDialog Fields Overview	. 117
	The SAMPLE_ORDER Table	
	The METHODS Table	
	The NMR_PREPARATION_WS Table	
	The Experiment Table	
17.2	VCU Request Tables	
	Introduction to VCU REQUEST Tables	
	The VCU Sample Order Table	
	The VCU Preparation Table	
	The Experiment Table (VCU_REQUEST_ALL)	. 122
18	EasyDialog Configuration	.123
18.1	Opening the EasyDialog Configuration Window	123
18.2	Opening a Form / Creating a New Form	123
18.3	The Configuration Window	124
18.4	Drag & Drop Objects	
18.5	How to Drag & Drop Objects	126
18.6	Grids	
18.7	The Property Editor	
18.8	The Fields Overview Window	
18.9	Fields Overview Required Values	
18.10	Default Values in the Fields Overview Window	
18.11	The STATUS Field	
18.12	Configuring a Status	
18.13	The Administrator Toolbar	
18.14	Example: Creating a New Input Form	
	Creating a New Form: Opening the Form	
	Creating a New Form: Establishing Group Boxes	
	Creating a New Form: Adding the First Object	
	Creating a New Form: Adjusting the Group Box	
	Creating a New Form: The Completed Group Box	
	Creating a New Form: Adding the Other Group Boxes	
	Creating a New Form: Adding Buttons	
40.45	Saving an EasyDialog Configuration	
18.15	EasyDialog Parameters	135
19	E-Mail Support	137
19.1	E-Mail Support	137
19.2	Installation of E-mail support	138
	System Requirements	. 138

	E-mail Internal Processes138
	E-mail Interface139
	E-mail Trigger Setup139
	E-mail Parameters140
	E-mail Interface Trigger Table141
	E-mail Address Registering142
20	Data Exchange & Directory Structure143
20.1	Introduction to Data Exchange 143
20.2	SampleTrack <sup>™</sup> bstsdata Directory Structure
21	JCAMP Definition File145
21.1	JCAMP Definition File Introduction145
21.2	Communication File with JCAMP Format 146
22	Script Library159
22.1	Introducing Scripts for Automation159
22.2	Common Variables and Functions 159
	Variables with Instrument Type: Interface
	Columns on Tab 'Interface File Settings'160
22.3	Common Terms for Script Processing 161
	Functions for Script Processing161
	Functions for File Import/Export162
22.4	Functions for Tables 164
22.5	Lookup Functions Database 167
22.6	Field / Value Evaluation168
22.7	Show Messages 168
22.8	Function: Send Emails 169
22.9	Special Functions 170
22.10	Function: Archiving 171
22.11	Obsolete Functions 172
22.12	Alphabetical List of Function and Procedures 173
23	System Requirements177
23.1	General Hint 177
23.2	Server Requirements 177
23.3	Client PC Operation System Requirements 178
23.4	Spectrometer Software Requirements 178
23.5	Directory Sharing with UNIX Instrument Computers 178
24	Oracle® 9 Server and Client Installation179
24.1	ORACLE® 9i Database Server Installation
	The Path in the Environment System
24.2	ORACLE® 9i Client Installation 186
25	Oracle® 8 Server and Client Installation189
25.1	Oracle® 8.1.7. Database Server on a Windows PC 189
25.2	Installation Method 189
25.3	Installing Oracle 8.1.7. Client on a Windows PC 195

26	Client/Server Network Connection	.199
26.1	Net8 Assistant	199
27	SampleTrack™ System and Client Installation	203
27.1	SampleTrack™ System vs. SampleTrack™ Client	203
27.2	Installation of the SampleTrack™ System	204
27.3	SampleTrack™ Patches	206
27.4	Running SampleTrack™ for the First Time	207
27.5	SampleTrack <sup>™</sup> Program Icons	213
27.6	ST Config - A Program for the Administrator	214
27.7	Database Connection Error After New Installation	215
27.8	Installation of Instruments	216
27.9	Installation of BRUKER NMR Spectrometer Using B-ACS	217
	1. Installation of a Dummy Instrument from a Template.	
	2. Make a Copy of the Dummy Instrument "SPECT"	
	3. Make this Instrument Available	. 220
27.10	SampleTrack <sup>™</sup> Client Installation	221
28	SampleTrack™ Web Client Installation	225
20	-	
28.1	Introduction / System Requirements	
28.2	Installation	225
	Installation of Program Files	
	Activation of the Web Server	. 226
	Secure Socket Layer (SSL-HTTP)	. 227
28.3	Configuration	227
	User Identification	. 227
	Group or Individual User Settings	. 228
	Navigation Board	. 228
	Layout	. 229
	Custom-page Wizard	. 229
	HTML Packages	. 230
	Rights	. 230

# **Copyright Notices**



#### Copyright © 2004 Bruker Biospin GmbH

All rights reserved. No part of this publication may be reproduced in any form without the written permission of Bruker Biospin GmbH.

SampleTrack<sup>™</sup> is a trademark of Bruker Biospin GmbH.

ICON-NMR<sup>™</sup> is a trademark of Bruker Biospin GmbH.

AMIX<sup>™</sup> is a trademark of Bruker Biospin GmbH.

Topspin® is a registered trademark of Bruker Biospin GmbH.

XWIN-PLOT<sup>™</sup> is a trademark of Bruker Biospin GmbH.

ProteinScape<sup>™</sup> is a trademark of Bruker BioSciences .

Biacore® systems are registered trademarks of Biacore AB.

Oracle® is a registered trademark of Oracle Corporation and/or its affiliates.

All Borland® and Borland® brands and product names are trademarks or registered trademarks of Borland.

Gilson product names are trademarked products or registered trademarks of Gilson, Inc.

 $\ensuremath{\mathsf{Tecan}}\xspace{\mathbbmath{\mathbb{R}}}$  product names are trademarked products or registered trademarks of Tecan Group Ltd.

Crystal Reports and Crystal Enterprise are trademarks or registered trademarks of Business Objects SA or its affiliated companies in the United States and other countries.

Apache is a trademark of The Apache Software Foundation, and is used with permission

Excel® is a registered trademark of Microsoft Corporation.

Word® is a registered trademark of Microsoft Corporation.

Microsoft® and Windows® and Microsoft® Internet Explorer are registered trademarks of Microsoft Corporation.

# **Copyright Notices**

Windows® Me is known officially as: " Microsoft® Millenium Edition operating system".

Windows® 2000 is known officially as: " Microsoft® Windows® 2000 operating system".

MDL® ISIS/Draw is a trademark of MDL Information Systems, Inc.

All product names and marks mentioned in this document are trademarks or registered trademarks and property of their respective owners.

# Warnings and Notes

There are two types of information notices used in this manual. These notices highlight important information or warn the user of a potentially dangerous situation. The following notices will have the same level of importance throughout this manual.



WARNING: Indicates the possibility of severe personal injury, loss of life or equipment damage if the instructions are not followed.

### **Contact for Additional Technical Assistance**

**BRUKER BioSpin GMBH** 

am Silberstreifen D-76287 Rheinstetten

Germany

Phone:

FAX:

Email:

Internet:

For further technical assistance please do not hesitate to contact your nearest BRUKER dealer or contact us directly at:

User Guide Version 002

**BRUKER BIOSPIN** 

or www.bruker-biospin.de

sampletrack@bruker-biospin.de

+ 49 721 5161 0

+ 49 721 5171 01

www.sampletrack.de

2.2

2.1



# Warnings and Notes

# Introduction to Bruker LIMS

3.1

### Laboratory Data Management Today

The steadily increasing number of samples in analytical spectroscopy means that laboratories have to deal with a massive data throughput. This is created on the one side by spectral data and analytical reports and on the other side by data archival, referencing and distribution. This situation, along with GLP requirements, FDA regulations, and ISO 9000 certification, means that the instrumentation in analytical laboratories have to be linked to an existing corporate LIMS. Increasingly, many chemical, pharmaceutical and biotech companies not only have their QA/QC processes integrated via LIMS, but also their R & D departments, where typically BRUKER NMR, FT-IR, and mass spectrometers are to be found. Therefore, many industrial users are now seeking to obtain a LIMS connection for BRUKER spectrometers. Traditional LIMS programs are laid out more for chromatography and for the very widespread, less costly analytical instrumentation. They do not usually include NMR spectroscopy. SampleTrack<sup>™</sup> has been designed to fill this gap!

<u>See also: SampleTrack Planning Questionnaire (http://www.bruker-bio-spin.de/NMR/automat/samtrac/samtrac6.html)</u>

or the SampleTrack<sup>™</sup> homepage ((http://www.sampletrack.de/)



Figure 3.1. Laboratory Data Management - LIMS

# Bruker Laboratory Information and Management System

SampleTrack<sup>™</sup> is a Laboratory Information and Management System with a standardized interface for BRUKER spectrometers. It is a software package that provides the link between the analytical laboratory, the synthesis laboratory and the information processing department. More importantly, it is a software tool for the laboratory network that coordinates all automation steps in analytical measurements. It monitors the activity on the laboratory floor, and provides chronological precise and detailed information on each step along a sample's path. Sample-Track<sup>™</sup> provides complete sample tracking for the NMR laboratory within a modern research environment.




Figure 3.2. SampleTrack<sup>™</sup> - Worksteps

### The SampleTrack<sup>™</sup> Interfaces

SampleTrack<sup>™</sup> offers a configurable, easy to use interface with all of the features necessary for the administration of day-to-day routine operations in laboratories, such as: sample log-in, bar code printing, report generation, lists of sample submitter's, and lists of sample preparation methods and requested experiments. These features, combined with the underlying database, provide a powerful organizational tool for laboratory managers.

SampleTrack<sup>™</sup> is designed to optimize sample tracking, beginning at the submitter's site and continuing through to the final return of results. Information management and hardware control are carried out concurrently as the sample proceeds from the synthesis laboratory through the analytical procedure. Information on the sample is constantly updated and made available to the laboratory personnel as well as to the submitter. Laboratory data is always on-line, active and accessible.

The typical environment of SampleTrack<sup>™</sup> is within the NMR laboratory. All worksteps can be processed individually from SampleTrack<sup>™</sup> or in conjunction with an external LIMS.

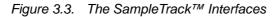



- SampleTrack<sup>™</sup> client administration tool <u>"SampleTrack<sup>™</sup> Environment Set-up" on page 85</u>
- <u>"Enter Orders Using the EasyWeb Dialog" on page 76</u> for easy order setup in the web
- EasyDialog for easy order setup <u>"Introduction to EasyDialog" on page 65</u>
- "WEB Client Introduction" on page 69
- <u>"Structure and Web Spectra Viewer" on page 72</u>

### SampleTrack<sup>™</sup> as Data Reference

SampleTrack<sup>™</sup> consists of a straight forward Client/Server architecture and intuitive user interface within the Microsoft Windows<sup>®</sup> operating environment. Document usage is easily accomplished through drag & drop selection, and the Client Tool provides easy access to all of your related data (e.g. spectra, structure diagrams, analytical reports etc.)

The SampleTrack<sup>™</sup> integrated automation allows you to perform standard and composite experiments without any limitations. And you will have access to your related data even after the archiving process.

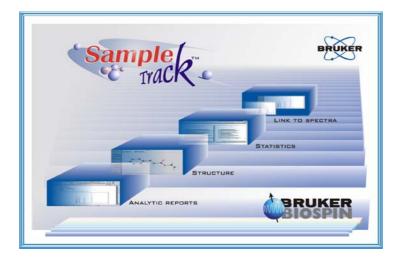


Figure 3.4. SampleTrack<sup>™</sup> as Data Reference

## **Daily Routine Laboratory Functions**

In its role as a local management system, SampleTrack<sup>™</sup> provides availability for all the daily routine laboratory functions, such as:

- Order and Report generation
- Supervision of the work routine/production
- Management of all clients
- Preparation methods
- The available and permissible experiments
- Barcode printing
- Order verification
- Archiving of all associated data

Figure 3.5. Daily Routine Laboratory Functions



### **Communication Overview**

Orders can be setup with EasyDialog, with the EasyWeb Dialog, within the SampleTrack<sup>™</sup> Client, or imported by an external LIMS system. When the status of the manual or automated preparation is set to "finished", the orders are sent to one of the existing spectrometers, and after measurement, reported back.

<complex-block>

Figure 3.6. Communication Overview

## SampleTrack<sup>™</sup> Workflow Architecture

In addition to sample tracking and procedures control, an important feature of SampleTrack<sup>™</sup> is the complete recording and documentation of the steps within the procedures. The sample data, path and status, as well as preparation, analysis, spectra transfer and storage are recorded in the local database on the server. This is an essential feature for laboratories operating within GLP regulations.

When no company LIMS is available, laboratory managers can take advantage of the recording and reporting features of SampleTrack<sup>™</sup> to document their laboratory studies, in order to utilize them for GLP documentation.



Figure 3.7. SampleTrack™ Workflow Architecture

### Integrative Bioinformatics Solution - A Look into the Future

Unique Proteomics RIMS<sup>™</sup> integrates multiple information-rich methods (MS/ SPR/NMR/XRAY) for increased research productivity in proteomics.

#### Proteomics RIMS<sup>™</sup> combines

- Bruker BioSpin's SampleTrack<sup>™</sup> LIMS (Laboratory Information Management System)
- Bruker BioSciences' ProteinScape<sup>™</sup> mass spectrometry software
- Bruker AXS X-ray crystallography technologies
- Biacore®'s surface plasmon resonance (SPR) technology for detection and monitoring of biomolecular binding
- Bruker BioSpin's NMR

Dr. Herbert Thiele, Director of Bioinformatics at Bruker Daltonics. "Proteomics RIMS is the logical consequence of Bruker BioSciences and Bruker BioSpin companies providing the entire range of systems solutions for protein MS, x-ray crystallography and NMR. The synergy of their combined and integrated usage is a major step towards increased productivity in drug discovery and target validation."

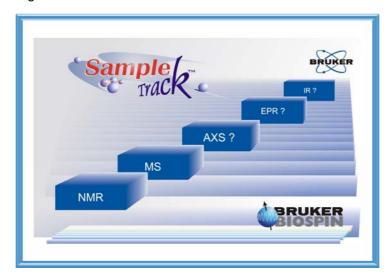


Figure 3.8. A Look into the Future

# Introduction to the SampleTrack™ Client?

### SampleTrack<sup>™</sup> Login

When first starting SampleTrack<sup>™</sup> Client you will be required to login by entering the appropriate login name and password. The login date and time will be saved at this time.


Login Screen for SampleTRACK Client

Figure 4.1. Login Window

4.1

#### SampleTrack<sup>™</sup> Overview

On the left side of the main SampleTrack<sup>™</sup> Client window is a SampleTrack<sup>™</sup> Overview window displayed.

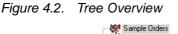




As shown in the figure, this window depicts a tree overview of the various displays that are available. With this overview window, as is customary with most Windows® based applications, the details that are available for each area are displayed when a module is selected. For example, selecting Sample Orders will display all of the current sample overviews that are available.

4.3

The Sample Orders module provides a good overview of the complete order, without having to change to other parts of the program. All global information, such as Sample ID, Client ID, Instrument, etc., as well as the status of the complete order, are displayed in the Sample Orders heading.

This module also allows you to add, copy and delete sample orders, and to view experiments. You must only click on the corresponding buttons to implement these routines.

Figure 4.3. The Main Window

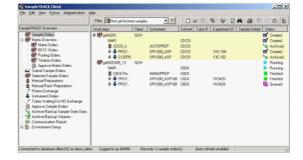

In the NMR method Sample Orders window (shown above), all the worksteps are clearly portrayed within the window. The assignment of individual experiments to the respective preparation steps can be accomplished through inserting appropriate values.

The details of the actual orders are displayed simultaneously in the Sample-Track<sup>™</sup> overview window.

#### 4.4

# ٥

For viewing the details of the individual preparation for a Sample order, doubleclick on the Sample ID line of the menu. This will result in the display of the sample order details window, as shown below.

\_ | D | × Filter 😡 gd 🖳 🔽 🕞 🛍 🦉 🖏 🗳 🖓 🖬 🎒 🚳 ۵ 🎼 Solvent Tube ID Exp Workstep: Ø 😻 gah EST 8 H 4 P H B A D 8 0 0 ٤. ments Start Date 29.04.2002 14: . End Date 29.04.2002 14 --\* Archi मन

Figure 4.4. Double Click to See the Sample Order Details

If you wish to view details of an individual experiment, e.g. a preparation, double click on the line of the corresponding preparation workstep. This will display the Preparation Setup window as shown in the figures that follow.

Pier


Figure 4.5. Preparation Details

Solvent Settings Preparation	Solution Method
Solvent Quantity	USONIC  Solution Duration
Dient Solvent	60 Second(s) Solution Temperature
Concentration in %	310 Kelvin
	Cloning Number of Clones
	Clone Quantity
	ug

ŧ

You can also view the details of an experiment from the Sample Order Details window by double-clicking on the corresponding workstep.

Figure 4.6. Experiment Details

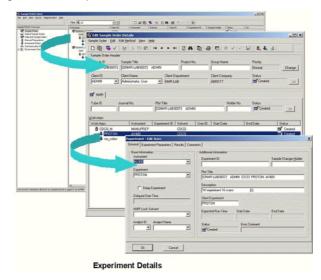

On the right edge of the sample order input window you will find **Buttons**, which allow you to quickly add a Preparation, an Experiment or a Proton Exchange.

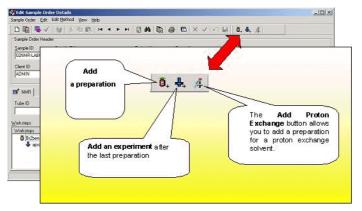



Figure 4.7. Add Preparation / Experiments / Proton Exchange Buttons

The Add Preparation button allows you to add a new preparation. The Add Experiment button allows you to add a new selected experiment The Add Proton Exchange button allows you to add a preparation for a proton exchange solvent.

### Sample Order Window Buttons

4.7

On top of the Sample Order window you will find **Buttons**, which allow you to quickly select the individual functions for order adjustment and modifications.

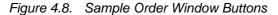



#### **Create New Sample Order Button**

D

With the Create New Sample Order you can add a complete new order. Pressing this button results in display of a blank Sample Order Details window.

#### **Edit Item Button**

1

The Edit Item button allows you to edit a workstep or the details of an order.

#### **Copy Sample Order Button**

With the Copy Sample Order button you may copy an order to form a new one or generate a set of new orders.

#### Submit Sample Order Button

The Submit Sample Order button is used to begin the measurement, or to restart an experiment with a **Finished** or **Error** status.

#### **Delete Sample Orders Button**

in the

With the Delete Sample Orders button you can delete all of the selected sample orders.

#### **Reload or Refresh Button**

t.

Reloads the data from the database. This can be automized in the menu Extras.

#### View all spectra Button

**b** 

With this button all spectra from one sample order will be displayed.

#### **View Experiment Button**

With the View Experiment button you have the possibility of viewing and manipulating the associated spectra of the experiment with an experiment viewer.

#### **View Next Experiment Button**

,

With the View Next Experiment button you can jump to the next existing dataset for viewing or manipulation.

#### **Expand / Contract Button**

8:

With the Expand/Contract Button items can be expanded or contracted.

Items expanded

Figure 4.9. Items Contracted

Filter 🛐 CY	x 💽 🗹 🔽 🖌 🖉			GA @	) 🖽 o 🖆 🎼 🗕 🛶		
Warksteps	Chert.	Instrument	Solvent	Tube ID	Experiment ID	Sample Holder	Status
Cr_020709002	CY						E Created
🗱 cy_05_07_2002_01	CY						E Finished
👹 cy_05_07_2002_2	CY						E Created
W cy_05_07_2002_3	CY			5678			Created
Cy_05_07_2002_4	CY						Created
👹 cy_05_07_2002_5	CY						E Created
😻 cy_05_07_2002_6	CY						E Created
👹 cy_05_07_2002_7	CY						Created
W cy_05_07_2002_8	CY						Created

Figure 4.10. Items Expanded

Fiker 🕵 CY		- Y	口运用	S 3		日回臣	-		
Worksteps	Client	Instrument	Solvent	Tube ID	Experiment ID	Sample Holder	St	atus	
W cy_020709002	CY	a in the second	1.	to the sheet			· 💷	Created	1
NMR			DMSO				1	Created	
DM50_M		MANUPREP	DMS0				m	Created	
C134PT		DPX300_ASP	DMSO		cy_020709002		1	Created	
/ DMSO_M		MANUPREP	DMS0+D2D				. m²	Created	
C13APT		DPX300_ASP	DM50		cy_020709002		100	Created	
10 2002 05 07 2002 01	CY						1	Finished	
NMR			Acetan				1	Finished	
CDCI3_M		MANUPREP	Aceton				- 🔳	Finished	
PROTON		DPX300_ASP188	Aceton		1H3618		- 📕	Finished	
CDCI3_M		MANUPREP	Aceton+D20				- 12	Finished	
PROTON		DPX309_ASP188	Aceton		1H3626		1	Finished	
W cy_05_07_2002_2	CY						m	Created	
NMR			Aceton					Created	
CDCI3_M		MANUPREP	Aceton					Finished	
PROTON		DPX300_A5P188	Aceton		1H3519		1	Finished	
COCB_M		MANUPREP	Aceton+D20				10	Created	18
PROTON		DPX300_ASP188	Aceton		1H3628		<u> </u>	Created	
W cy_05_07_2002_3	CY						<b></b>	Created	
NMB			Aceton	5678				Created	

In the Sample Orders window it is also possible to receive the complete status information for the selected orders.

Figure 4.11. Status Information

	Faer Tala		- V	口學的專	1 E A & 0	00 1		
	Worksteps	Ciert	Instaument	Solvent Tube ID	Esperiment ID Sample	Holder Sta	hus .	
Sangle Deckers	ABC123#4277	ADMN				1	Approved	
E 👹 Matix Overview	NMB			D20		-	Approved	
Matix Order:	TECAN LPD		TECAN	028		1-2	Approved	17
BEST Orden	# # PROTON		DPX300_SAMP.	D20	ABC123#4277	$i \sim$	openand	
Pooling Enders	ABC123#4278	ADMN					a second	
- W Tittetion Orders	NMR			D20		1	Addressed	
Approve Matrix Di	C TECAN_LPD		TECAN	D20	1	1	proved.	
Selected Sample Orders	0 🖶 PROTON		DPX300_SAMP	D28	ABC123#4278	1	restrict	
Manual Preparations	ey 05_07_2002_01	CY			and the second second	5	Created	
A Marual Rack Prepara	NMR			Aceton		5	Created	
- Proton Exchange	CDD3_M		MANUPREP	Aceton		<b>1</b>	Created	
A Instrument Orders	PROTON		DPX300_ASP188	Aceton	1H3618		Created	
7 Tubes Waiting For HE	Wey_05_07_2002_2	CY				i 🖬	Created	
Approve Sample Bide	NMR			Aceton		N 🖬	Created	
Archive/Backup Sam	CDC3_M		MANUPREP	Aceton		/ s'	Created	
- Archive/Backup Volu	PROTON		DPX300_ASP188	Aceton	1H3619		Created	
- Communication Report	W cy_06_07_2002_3	CY.					Consted	
注意 Environment Setup	NMR			Aceton		<b></b>	Created	
	COC3_M		MANUPREP	Aceton			Created	
	- PROTON		DPX300_ASP188	Aceton	1H3820		Causted	
	Wey_05_07_2002_4	CY					Created	
	NMR			Aceton			Created	
	CDC3_M		MANUPREP	Aceton			Created	
• •	PROTON		DPX300_ASP188	Aceton	1H3621	<b>5</b>	Consted	

Figure 4.12. List of all Status Symbols

10		Created				
20	65	Verified/Submitted				
25		Waiting				
30	Fair	Queued				
40	1	Started/R unning				
50		Paused				
60		Finished				
70	~	Approved				
80	*=	Archived				
85	6	Archived on CD				
95	0	Instrument Error				
99	X	Canceled				

## Filters & Views

Specific client or department fields and their views are managed by the filter & view assistant.

Figure 4.13. Filter & View Assistant Button

V

**Basic filters** select only those orders which the user or their department are principally allowed to see.


Within this personal data pool user specific filters can be set.

🖉 Filter & View Managem				<u>_   ×</u>
Standard Filters   Base Filters	View Fields Available Fields	N		
Availat ields		<del>ር</del> ጎ		
Field Title	Table Name			
☑ ♦ Sample ID	Sample Orders	Composite Fields		
☑ ◆ Client ID	Sample Orders	Field Title		
🔽 🔷 Status	Sample Orders	Client ID		
🔽 🔷 Sample Title	Sample Orders			
🗹 🔷 Project Number	Sample Orders	Composite Field	Info	
🗹 🔷 Mod Date	Sample Orders	1. Level Field		Variant (HTML)
IP ◇ IP	Sample Orders	Client ID (San	nole Orders)	
🗹 🔿 Comment	Sample Orders			
Not CP	Sample Orders	2. Level Field		Variant (HTML)
🗹 🗘 Poison	Sample Orders			• •
Back	Sample Orders			
Smell	Sample Orders	3. Level Field		Variant (HTML)
🗹 🔷 Method Type Key	Methods			
🗹 🔷 Tube ID	Methods	4. Level Field		Variant (HTML)
Plot Title	Methods	4. Level 1 iolu		
Journal Number	Methods			
Status	Methods	5. Level Field		Variant (HTML)
	Preparations	<b>•</b>		
LIVI ♀ Pade	Methods		-	
Check Selected It	ems 🛛 🔚 Save Fields			
Uncheck Selected It	ems 🛛 😰 Refresh View			
			I	
			Ok	Apply Cancel

Figure 4.14. Filter and view management

The users might be allowed to define their standard filters but are not allowed to define basic filters and available fields. In order to use this option, the necessary user privileges must first be assigned by the Administrator in the Object Manager as shown below.

Figure 4.15. The Object Manager

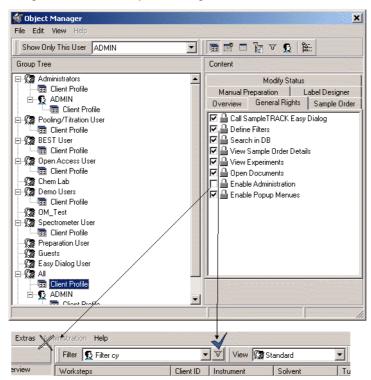

When the **Define Filter** box is checked and the **Enable Administration** box unchecked as in the example above, the user only has minimal rights to define their personal views and filters.

🖋 Filter & View Management			
Standard Filters Base Filters View F	Fields		
🕼 Manage Filters		Filter Name	
Available Filters			
Filter Name	Owner		
マ 01 Created Samples マ Filter 2	All Cyppel, Gilbert	🚺 🗖 Sample ID (Sample Orders) 🛛 🖛 💌	) and 💌
🗹 🐨 Filter JEA	Cyppel, Gilbert	Client ID (Sample Orders)	▼ ) and ▼
♥ ♥ Filter cy ■ ♥ 02 Verified Samples	Cyppel, Gilbert All	👘 🔲 Status (Sample Orders) 👘 🔳	• ) and •
□ ♥ 03 Queued Samples         □ ♥ 04 Prepared Samples         □ ♥ 05 Finished Samples (This         □ ♥ 07 Archived Samples (This         □ ♥ 08 Archived On CD (This M         □ ♥ 08 Archived On CD (This M         □ ♥ 09 Not Yet Finished Samples         □ ♥ 10 Not Yet Approved Sample         □ ♥ Filter 1         □ ♥ Filter 3         □ ♥ Running Samples for DFX3	All All All All All All Cyppel, Gilbert		Y ) and Y
		Ok	pply Cancel

Figure 4.16. Filter Management

When the **Define Filter** box and the **Enable Administration** box are checked, which is usually a privilege for the administrator, the entire functionality is available:

👹 Filter & View Management		
Standard Filters   Base Filters   View Fields   Available Fields		
Manage Groups 🕼 Manage Filters	Filter Name	Group Name
Available Filters		
Filter Name Owner		
v v Filter 1 All	🚺 🗖 Sample ID (Sample Orders)	) and •
√ ▼ 01 Created Samples All		
V ▼ 02 Verified Samples All	( 🔲 Client ID (Sample Orders) =	() and (
√ ▼ 03 Queued Samples All		
√ ▼ 04 Prepared Samples All	🚺 🔲 Status (Sample Orders) 📃 💻	r ) and 💌
√ ▼ 05 Finished Samples All		
V V 06 Approved Samples (This All		
√ ▼ 07 Archived Samples (This All	김 동물에서 가장을 물러 물건을 받았다. 것이	아님은 것은 것이 같은 것은 것은 것을 것 같아.
✓ ▼ 08 Archived On CD (This M All	김 한 것이 있는 것은 것은 것을 것을 것 같아.	
√ ▼ 09 Not Yet Finished Samples All	김 화장님은 것은 것이 아파 영화를 얻는 것을 했다.	
√ ▼ 10 Not Yet Approved Sampl All		
Groups		
Group Name Group		
Easy Dialog User 01		
Guests 02		
Show Users		
🗌 Uncheck All 🛛 Check All 🔄 Save Item 🛃 Refresh View		
		Ok Apply Cancel

Figure 4.17. Filter Management

For further information see <u>"Filter and Search Tools" on page 54</u> and <u>"Insert-ing a New Filter Method" on page 109</u>

# Functions of the SampleTrack™ Client Module

5.1

#### Submit Sample Orders

Depending on the workflow in the laboratory, the order must be handled by the official person or set automatically by the SampleTrack<sup>™</sup> system.

In the following scenario the order is verified manually. After creating an order using EasyDialog or within the client the status of the order remains **Created**.

Figure 5.1.	Scenario 1
i igui e ei ii	econario i

	Interaction	Interaction	Interaction
Status	Create 📕	Submit	Manual Preparation and Supply Tube ID
Sample Order	Created	Verified	Running
Method	Created	Verified	Running
Preparation	Created	Verified	Finished
Experiment	Created	Verified	Verified

In the next scenario the step **Submit Order** is skipped, whereas the next step is the **Manual Preparation**:

Figure 5.2. Scenario 2

	Interaction	Interaction
Status		Manual
Status	Create	Preparation and
	<b>•</b>	Supply Tube ID
Sample Order	Verified	Running
Method	Verified	Running
Preparation	Verified	Finished
Experiment	Verified	Verified

Another scenario could be a LIMS Import by an external LIMS workaround with prepared NMR samples:

	Interaction	Interaction	
Status	Sample Orders   by LIMS	Submit	
Sample Order	Created	Running	
Method	Created	Running	
Preparation	Finished	Finished	
Experiment	Created	Verified	

A completely automized workaround, without interaction, is possible when the orders are imported and immediately submitted by the system.

The area Submit Sample Orders on the left side provides a method of displaying an overview of the manually created and/or imported orders, and sample orders with an **Error** status.

Figure 5.4. Status Overview

	ignistration Belp						
	Filter All		I 🖸 🖉 🖬	1 🐴	9 B # 6 (	3 II ( ) III ( )	
stiple1RACK Overview	Worksteps	Diere IC	Indument	advert.	TubeID	Statun	
👹 Sample Diders	W cy 05 07 2002 01	CY	1			Created	
Matrix Overview	NMR		/	Aceton		Created	
Matrix Orders	CDCB M		MANUPPEP	Aceton		Created	
- BEST Ordero	PROTON		DPX308 ASP188	Aceton		Created	
Pooling Orders	cy 05 07 2002 2	CY	AND A DECK OF A DECK	a desired	103035200040319530	Created	100
- W Titration Daters	NMB		1	Acetan		Created	_
Approve Mahis Dr	6 CDCB M	1	MANJPREP	Aceton		Created	
Submit Sample Oxden	PROTON	1 .	DPX300_A5P188	Aceton.		Created	
- Manual Preparations	W 05 07 2002 3	188/				Created	
Marual Back Prepara	NMR	11		Aceton		Created	
- A Proton Exchange	CDCB M	~	MANUPREP	Aceton		Created	
A Instrument Orders	PROTON		DPX300_A5P188	Aceton		Created	
Tubes Waiting For HE	Way 05 07 2002 4	CY.				Created	
Approve Sample Dide	NMR			Aceton		Created	
Archive/Backup Sam	CDCB_M		MANUPREP	Aceton		Created	
Archive/Backup Volu	PROTON		DP/(300_ASP188	Aceton		Created	
😂 Communication Repor	ay 05 07 2002 5	CY				Created	
🙀: Environment Setup	NMR			Aceton		Canaled	88 T
	CDCB_M		MANUPREP	Aceton		Canadad	
	PROTON		DPX300_ASP188	Aceton		Canaled	
	W oy_05_07_2002_6	CY				Created	
	NMR			Aceton		Created	
	CDCI3_M		MANUPREP	Acaton		Cented	
×	PROTON		DPX300_ASP188	Aceton		Created	

The orders can be verified from here using the **Submit** button.

Figure 5.5. The Submit Button



Select All - use the right mouse click to mark all of the orders as Submitted

Figure 5.6. Select All


Select All	Strg+A
	 100 million (100 m

When a modification of an order is necessary, you can automatically move to the Sample Order Details window by double-clicking on the order.

## **Manual Preparation**

This module is used by the preparation work area. Orders must have a submitted/ verified status in order for the Sample ID to be reported here.

Figure 5.7. Tree - Manual Preparation Area

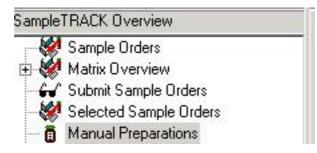

The goal of this module is that the preparer needs only to provide the minimum information necessary to confirm the implemented preparation and with it the generation of the necessary instrument order.

To display the respective preparation data of an order in the Manual Preparation window, the Sample ID should first be selected by using a barcode reader, by entering the ID number manually, or by selecting the ID number from a pull-down menu.

Sample Order Information	Preparation	Polution Cottings
<u>S</u> ample ID	Sample Settings	Solution Settings
03NMR-LAB09015	Sample Quantity	Solution Method
Sample Title	ug 🔽 Filter	
03NMR-LAB09015 234-342 ADMIN	Minimal Temperature Maximal Temperature	Solution Duration
	Kelvin Kelvin	Second(s)
Client ID		Solution Temperature
ADMIN	Solvent Settings	Kelvin
	Solvent Name	Proton Exchange
Tube ID	CDCI3	
	Solvent Quantity	
Einished Cancel	J W	
	Experiments	
	Experiment Instrument HD E	× Modify Experiment List
	lige aproton_schnell A∨400	Set Experiment Instrumer
		Set To Instrument

Figure 5.8. The Manual Preparation Window

If the solvent that has been selected is not applicable, a different solvent can be selected from this dialog window.

# Functions of the SampleTrack<sup>™</sup> Client Module

The instrument that is used for the experiment can also be selected or changed in this window.

Once the preparation has been performed and the barcode collar placed on the tube, the Tube ID from this collar must be entered in the Tube ID block in the Manual Preparation window.

After assigning a Tube ID, the order is **submitted** by pressing the **Finished** button.



Figure 5.9. B-BCE for Easy Tube ID Barcode Reading

An instrument often used for this purpose, the B-BCE, is connected with the computer and can automatically fulfill this procedure.

## Proton Exchange

This module has been created to give the preparer the possibility of searching for an order associated with a sample tube ID.

After the preparer adds the proton exchange solvent to the sample in the lab, the new order for the spectrometer is automatically generated by pressing the Finished button.



SampleTRACK Client			<u> ×</u>
File Edit View Extras A	Administration Help		
ampleTRACK Overview	Sample Order Information	Sample Title	Proton Exchange Settings Proton Exchange Solvent
Selected Sample Ord			
	Sample ID	Client ID	EnishedCancel
- 은 Communication Repol 한 않는 Environment Setup			
onnected to database dben:	152 as ki abbott	Logged in as ADMIN Records: 26 sample o	rder (slAuto refresh enable)

#### Finish Manual Experiments

When experiments from a sample order have been measured on an instrument which is **not connected** to the SampleTrack<sup>TM</sup> system it is necessary to manually set the status of the experiment to **Finished.** This is necessary to be able to use the Order Approval window to approve the orders and release the Tube ID for reuse.

Figure 5.11. The Finished Button

When you press the Finished button, all selected experiments are set to the **Finished** status. If all experiments in an order have been completed, the complete sample order is also set to **Finished**.

5.4

You are then able to use the **Order Approval** window to approve the orders and release the Tube ID.

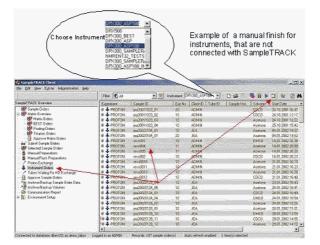


Figure 5.12. Example of a Manual Implemented Finish

#### **Delay Experiments**

In some cases it is desirable to delay the measurement from a specific experiment or experiment set. For example, to allow all C13 experiments to be run over the night or weekend in order to save the spectrometer time for proton experiments during the day.

Figure 5.13. Status Icon of Waiting / Delayed Experiments t

# 

The Delay Experiments window and its button allows you to set these experiments to the **Delayed** status for processing at a later time. The pull-down menu **Experiment Type** allows you to select various experiment types (nucleus). Experiments that have been delayed have a **Waiting** status.


Figure 5.14. Delay Experiments

S Samphi 3/4 ( CAH)	1m  10	e 07.2002 e 00.00	× ×						~(0) ×)
	fitze Date	a II	and shad been	PX300_ASP188 💌		50 b	📕 🥡 🕹	3 4	8 K
Elempiel PIALX, Oververv	Experiment	Sanple ID	ExpNo	Clevil TubelD	Sangle Hot	School	C StanDara		End +
Sample Cadersi	0 🕹 PROTON	#520011023_01	50	0.704.99		CDUB	24 10 2001		24.14
S - St Hats Dyarvay	0 🕹 PROTON	ins20011623_02	10	ACHIN		CDCI3	34,12,2001	217.00	24.1(5)
- Minis Gideo	u 🕹 enarou	#520011023_03	10	40MW		Acerone	24,15,2001	122200	24,11
-80 DEST Order:	V 🕹 PROTON	ma20011023_04	10	ADM N		Acesse	25.10.2001	15.4200	35.0
Pooling Orden	PROTORS & 0	jpa20020:34_01	10	JEA		Actives	04 0# 2082		0.10
- Of Lindian Diden	NO TORY & D	pa20020904_01	20	JEA		Applore	34.01.2082	4,52:00	04.0
Submit Sample Orders	и 🕹 Ряргоч	NPVD-021	10	ADMAN		CINED	14.0120021		14.0
Selected Sample Orders	0 🕹 FRDTON	sevo004	11	ADMN		Acesse	14/85.20021		14.0
B Manual Preparatory	0 🕹 FROTON	MN/0005	11	ADMIN		Applore	14 (2 2002)		14.0
- 5 Manual Rack Preparations	0 👆 FROTON	NV0002	10	ADMIN		CDCD	14 01 2012		14.0
/ Proton Exchange	0 🐥 Marun	DF33avee	10	ADMIN		Acetone	21 07 2202		21.0
Inthiner#Oider:	0 🕹 PROTON	HV/00011	10	ACM 32		Acaione	21.61 2002		21.0
7 Tubes Wating For HD Exchange	0 🕹 PROTON	NevroR013	10	326426		Actione	21.04.2002.1		210
Approve Somple Diden:	NOTORS 🐥 0	E1000/M	10 .	000430		CDCI3 ·	21.01.2002	184200	21.0

# Experiments Waiting (Delayed)

5.6

The **Start** the button within the Instrument Orders Window used to start the measurement of **Waiting (delayed)** status experiments.

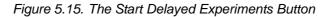

To start the experiment you must only highlight an experiment and press the **Start Experiment** button. If you have multiple experiments you can check the **Start Box** to the right of the desired experiments (or use **Select All** to select all of the experiments) and then press the **Start Experiment** button to begin the experiments.

You can also edit a selected sample order from this window, before starting the experiment, by pressing the **Edit Sample Order** button.


	Film 🕵 🕯	1	. 19	Indouneet D	PX300_ASP18	33 .	000	5 II I		1	3 4	0
ampial RADI: O verview	Expenser	Sarce	D.	ExpNo	Ciercito 1	TubelD	Sanple Hol	Schowd	514	Dist.		End
- 😸 Sanple Criters 🔹	0 🕹 PROTO		11023_07	SUST SALES	NING			CDCR			104200	260
Maria Gideo	0 4 PR010		11023_02 HIDBX_03	10 (2010)(9)(7)(2)	ADMIN ADMIN			CDCR Aceiore			12 17 00	24,11 9/28/16
DEST Dides	0 & PRDTO		11023_04	10	ADMIN			Acetore			194200	25.1
Profing Orden:	0 🕹 PROTO		201010101	iajaijas) ( <b>K</b> irak G	o gegeneel i	6446046076		Acoure			14 37.00	000
Approve Makis Orders	0 🕹 PROTO		20104_01	20 2029/2018/10/2	JEA U alforationes			Azeisee			14.52.00	040 949
Gul Subrit Sample Oxdero	0 - PROTO			10	ADMIN			Acatore			2010000	14.0
Selected Sample Under     Manual Perparations	0 🕹 PROFE			1004.040 <b>1010 :</b> 4446:	AINE			Acetera			302010	448
Manual Pack Preparations	0 🕹 PRDTD			10	AGMIN			CDCB			20:22:00	14.0
Proton Exchange	0 - PROTO			010001518-003	CREWNICS &			Azetore			101000	~210
Instrument Didets Tuber Wating For HD Exchange	0 4 PROFO			10 society?ppinioni	ADMIN COMIN			Azelore Azelore			16 22 00	21.0
Approve Sangle Orden	0 & PRDID			10	ADMIN .			CDUR			394800	21.0

Figure 5.16. Highlight and Start Delayed Experiments

#### **Experiments Waiting for HD Exchange**

This module is NMR specific. In this window all of the orders that are waiting for the addition of the solvent for the proton hydrogen/deuterium exchange can be viewed or printed out in report form. As with the Experiments Waiting (Delayed) you can also edit a sample order from this window.



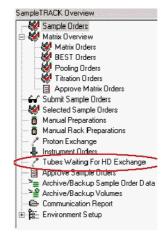

The HD Exchange Solvent window is opened in the detail window using the folowing button



Figure 5.18. The HD Exchange Window

	ADABY/ODAL
Sample Dider Header	nt lane Graphane
	beral Change
reparation - Add HD Exchange Solvent	×
General	
Base Information	Extended Information
Instrument	Start Date
MANUPREP	03.07.2002 14:48:40
Preparation Short Name	End Date
Solvent Name	Expected Bun Time
Aceton	
Add HD Exchange	김 중 이상입지만 동작하는 것이라. 전
	Status
	Created
Preparer ID Preparer Name	Error Comment
CY * Cyppel, Gibert *	
Carlo and a start of the start	and the second
Ok Cancel	신경 화장님 집 집을 알려야 한 것이 없다.

When the previous experiment within the order is finished, the exchange experiment can be found and handled in the **Tubes Waiting for Exchange** window.

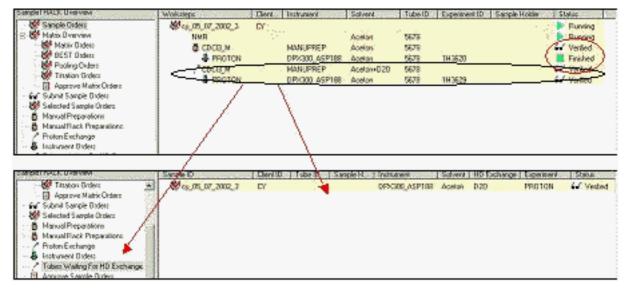


Figure 5.19. The Tubes Waiting for Exchange Window

#### Order Approval

The Order Approval module allows you to approve finished orders so that you can **reuse Tube ID barcode collars**. Individual reports can also be generated for viewing or printing by using corresponding filters.

With this module you may approve all finished orders at once (Select All / CTRL A) or you may select orders individually by clicking on the **Approved** box with your mouse.

To reverse the **Select All** process, press Refresh, which will remove the checkmarks.

Through this process the header status is set to approved and the Tube ID is removed from the valid data field releasing it for reuse.

Once the order has the **Approved** status, no other changes are possible, unless the order is not reset.


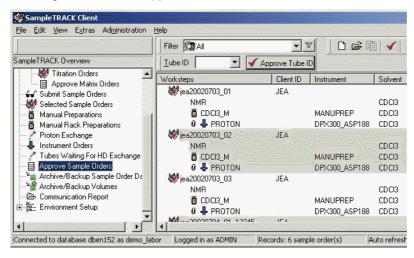


Figure 5.20. The Approve Window

The B-BCE can be used to easily approve an order and release the Tube ID.

Figure 5.21. B-BCE

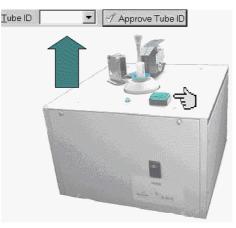

# E-mail Notification

5.9

It is also possible to provide automatic customer notification via E-mail , when the order is complete.




With this E-mail it is possible to send the spectra data as an attachment.


See also: "E-Mail Support" on page 137

#### Environment Setup

5.10

The Environment Setup depicts the area, where all of the adjustments and definitions for the complete system environment are implemented. Refer to the Sample-Track Environment for further information <u>"SampleTrack™ Environment Setup"</u> on page 85.

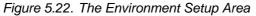

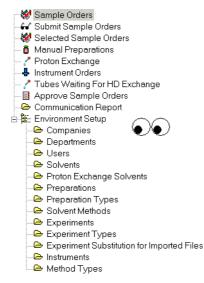

# Working with the SampleTrack™ Client

Sample Order Details Window

To set up a sample order move to the Sample Order window and press the Add Sample Order button.

This will open a new,	empty Sample Order Details window.
-----------------------	------------------------------------

Figure 6.1. SampleTrack Order

🖇 View Sample Order Details			
Sample Order Edit Edit Method View He	p		
	H - F - B B A D - 6 D	n i seen sen i 2000 📶 i noomaan	
Sample Order Header			
Sample ID Sample Title	Project Name Group	Name Priority	
1		Normal 333	Change
] ]	J J	Normal	Change
Client ID Client Name	I I I	Company Status	

Entering the Sample and Client ID

Enter your Sample ID and optionally a Project No. and Group ID. Using the TAB key or mouse move to the client ID field and enter the Client ID, or select it from the list of values from the pull-down menu.

6.2

6.1

	-  🗰   🗞 🕮 🖽   1			⊜  <b>⊡</b>  ×√ ∽			
ample Order Head	er Sample Title		Project Name	Graup Marca	Priority		
ample 10 cy_090702_01	test2		test		Normal	Change	
lient ID States and	Client Name	Client Dep		Client Company	Status	<u></u>	
ADMIN	Cyppel, Gilbert	NMBENH		BA	Created		
demo_open_acces demo_samplerai demo_tecan GDH GDE GUEST JEA KLE MBE REST TG THW	5						

Figure 6.2. Entering the Client ID

# Define a Method and Enter the Journal Number


6.3

To define a method like NMR, BEST, ... click the right mouse button:

Figure 6.3. Defining a New Method Using Right Mouse Button

New Method 
Method BEST-NMR Strg+1
Delete Method Strg+D
Method NMR Strg+2

or press Ctrl 1, Ctrl 2, ...

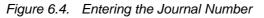

aksteps Underson Instrument Examinant ID Exhand Elect Data End Data Status	xksteps	
Industry Industry Example Contract Clark Data End Data Status		
Tonsteps industrian Experimental Survers Star Date End Date Status	Vorksteps Instrument Experiment ID Solvent Start Date	End Date Status

Optionally enter your journal number in the Journal No. block. As you move to this block the header of the order is automatically stored in the database.

When you save the order the Edit Sample Order Details window will appear.

	New Jonana and C. H. Charachi, C. Sta					
	Edit Method View I					and the second
D 💼 🖷 🗸	이 🗞   🐰 🏨 💼	H A 🕞 🖬 🔛	<b>m</b> 🖻 🥭 !	🖻   🗙 🖓   🏷	🕞 🛛 Ö+ 🖶 🌾	
ample Order Head	ler				ener Silsentinen lier	
ample ID	Sample Title	Proje	ect Name Gro	up Name	Priority	
cy_020709002	cy_020709002/CY/				Normal	Change
Client ID	Client Name	Client Departmer	nt Clie	nt Company	Status	
CY 🔻	Cyppel, Gilbert					the second s
	Journal No.	NMRENH     Plot Title     [cy_020709002/CY/]	BA	Holder No	Status	>>) 
- NMR   ube ID orksteps Vorksteps	Journal No.	Plot Title 	BA Solvent	Holder No	Status	>> Status

Figure 6.5. The Edit Sample Order Details Window

#### Select the Preparation Workstep

Move to the Worksteps block and press the Add Preparation button to enter the necessary preparation. A Preparation Setup window will be opened, whereby you can select the desired preparation from the pull-down list of available preparations located under the Preparation block.

#### **0**+

Figure 6.6. Select a Preparation

Preparation - Add Preparation	×
General Sample Preparation Pooling/Titration	Comments
Base Information Instrument MANUPREP	Extended Information Start Date
Preparation Short Name	End Date
Solvent Name Acetic Aceton Acetone ACND20	Expected Run Time
Aethanol Br2ben Buffer pH7 CDCl3 CD2Cl2 CD3CN CH3CN+D20 Cl2ben C6D6 C6F6	Status Created Error Comment

6.4

Upon selection of a preparation, all predefined values will be displayed. You may modify them as needed, or accept the values by pressing the OK button.

The preparation entry is now complete and will be displayed as a workstep in your order. The Sample Title and Plot Title is automatically created, using your pre-defined function, and entered in the database.

Edit Sample Or Sample Order Edit	der Details Edit <u>M</u> ethod <u>Vi</u> ew <u>H</u> e	lp			<u>.</u>	<u>-     ×</u>
	a national sector and the		M D 3		🕞   ō, 4, 4	
Sample Order Hea	der					
<u>S</u> ample ID	Sample Title	Projec	t Name	Group Name	Priority	
cy_020709002	cy_020709002/CY/				Normal	
Client ID	Client Name	Client Department		Client Company	Status	
CY .	<ul> <li>Cyppel, Gilbert</li> </ul>	MMRENH	and a second	BA	Created >>	
Tube 1D	Journal No. cy_AHA03481	Plot Title 		Holder No	Status Status	
<u>W</u> orksteps Worksteps	Instrument	Experiment ID	Solvent	Start Date	End Date Status	
DMSO_M	MANUPR		DMSO	Stait Date	End Date Status	

Figure 6.7. Order Details

#### Adding Experiments

**₽**+

Your order is now ready for you to define the experiments that are planned. Press the button Add Experiment.

A window for inserting new experiments will be opened.

nstrument DPX300_ASP	Experiment ID		Sample Changer <u>H</u> older
Experiment	Plot Title		
AL27ND B112G C13APT C13CPD C13CPD02 C13CPD5N C13CPD5N C13CPd5N	Description Client Experiment		
C13DEPT135 C13DEPT135p C13DEPT45 C13DEPT90		art Date	End Date
C13GD C13HUMP C13IG C13MULT	Status	Error Comment	

Figure 6.8. Add Experiment

Select the desired instrument and experiment from the respective pull-down menus and change the Analyst blocks if necessary. To commit the new experiment press the OK button. The experiment will then be added and displayed as the next workstep in your sample order. If required, you may enter additional experiments in this same manner.

#### Adding D20 for Proton Exchange

6.6

If you want to add D2O for a proton exchange experiment to your order, press the Add Proton Exchange button located in the Sample Order Details window. A small Preparation Setup window will appear, allowing you to add the D2O by placing a check in the Add D2O box. Pressing the OK button will return you to the Sample Order Details window, while simultaneously adding the D2O workstep (a workstep with D2O added is shown in the following figure).

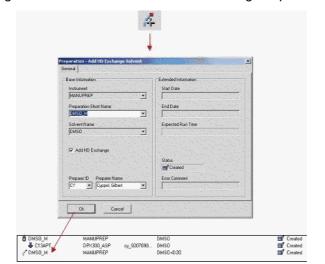


Figure 6.9. Add D2O for a Proton Exchange Experiment

Add the associated proton experiment as described above (see Adding Experiments) or use the standard Window's copying procedure to copy an existing proton experiment.

# Copy an Existing Experiment

6.7

To copy an existing experiment:

- Press and hold down the Control key on your keyboard.
- Select the gray area to the left of the existing proton experiment with your mouse, and drag the experiment into the gray area under the new preparation.
- Release the mouse button and Control key. The experiment should now be copied under the new preparation. The following figure is a demonstration of this process:

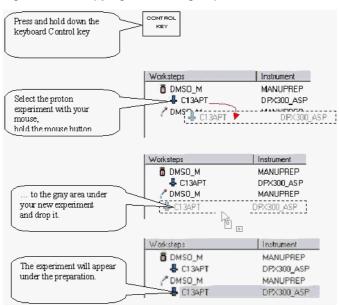


Figure 6.10. Copying an Existing Experiment

Figure 6.11. A Sample Order Created Using the Copy Function

]   √	8			\$ . <b>⊡</b> ` ×~√ . 	> 🖬 🛛 🗖₊ ♣₊ 🌾	
ample Order Head	er Statistics (Statistics)					
<u>S</u> ample ID	Sample Title	Pro	oject Name	Group Name	Priority	
cy_020709002	cy_020709002/CY/				Normal	Change
Client ID	Client Name	Client Departme	ent	Client Company	Status	
	1					
CY <u> </u>	Journal No.	Plot Title [cy_020709002/CY/		BA Holder No	Status	
NMR       Inde ID     orksteps	Journal No.	Plot Title [cy_020709002/CY/		Holder No	Status	
Vorksteps	Journal No. [cy_AHA03481 Instrument	Plot Title cy_020709002/CY/			Status	>>>
NMR       Inde ID     orksteps	Journal No.	Plot Title 	DMSO	Holder No	Status	
Vorksteps	Journal No. [cy_AHA03481 ] Instrument MANUPRE	Plot Title (cy_020709002/CY/ Experiment ID :P :P :P :P :P	DMSO 12 DMSO DMSO+D2O	Holder No	Status	🛒 Created

#### Starting a Preparation and Measurement

There are two methods available for starting the preparation and/or measurement:

- Standard Method
- Manual Method in the Sample Order Detail Window

Standard Method	6.8.1

The standard method used to start the preparation and measurement is to use the Submit/Verify Sample Orders Tool located in the SampleTrack Overview window.

Figure 6.12. The Submit Window



With this tool all external LIMS imported or manually generated sample orders are viewed and submitted for measurement.

You can submit the sample orders individually with the submit button, or all together by pressing the Select All button (CTRL A) and the Submit button.



The relevant orders for the instrument will then be generated for all submitted sample orders.



Note: Delayed experiments will not be generated until the start time has elapsed.

Manual Submit in the Sample Order Details Window

6.8.2

6.8

As an alternative method it is also possible to start an individual order generation by using the **Submit Sample Order** button in the Sample Order Details window.


This method is effective only when you have a limited amount of orders.

## Rerunning a Faulty or Finished Experiment

Mark the order and select one of the following options:

- Reset complete Sample Order
- Reset Sample Order Experiments

#### Figure 6.13. Rerun an Experiment

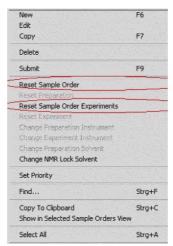


Resetting a distinguished experiment or a preparation is possible in the Sample Order Details window. Select the desired faulty or finished experiment or preparation by marking it, and use the right mouse button to select Reset Item (Ctrl R). This will set the status to Created.

Now resubmit the order.

In the case, where the whole order with preparation should be restarted, use Reset Sample Order. Submit the order to start the experiment.

#### Reset Item to Rerun the Complete Order

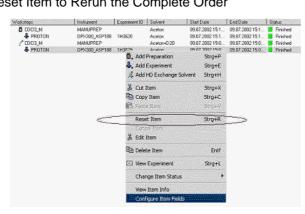

6.10

#### Adding an Experiment to a Finished Sample Order

At times it may be desirable to add an additional experiment to a finished sample order. To do so please follow the instructions below.

- Move to the Sample Order Details window of the relevant sample order by double-clicking on the order.
- Click once on the last experiment of the preparation in the Worksteps block and press the Add Experiment button. The Insert a new Experiment window will appear.
- Select the desired Instrument and Experiment data and update the Analyst blocks if required.
- When the parameters are acceptable confirm them by pressing the OK button.
- Start the measurement of your new experiment by pressing the Verify/Submit Sample Order button.

Figure 6.14. Adding an Additional Experiment

ase Information Instrument	Additional Information		
DFX300_ASF	Experiment ID		Sample Changer Holder
Experiment C134PT	Plot Title		-!
C Delay Experiment	Description		
	Allached Proton Test o	using imod pulse program	
Delayed Start Time	Client Experiment		
NMR Lock Solvent	Expected Bun Time	Start Date	End Date
Analyst ID Analyst Name	Status	Error Comment	
의 의 의	Created		
Ok ( Cancel			The second
steps	Experiment ID   Solvent	Start Date	End Date Sta
CDCI2 N MANUPREP	Aceton	09.07.2002.15/1	09.07.2002.15-1

#### Filter and Search Tools

#### Filter Tools

6.11.1

SampleTrack<sup>™</sup> runs in combination with the relational Oracle database, therefore, the displaying all of the stored sample orders is quite inefficient. The use of configurable filters reduces this inefficiency by allowing you to select only the samples that meet your search criteria.

Many of the window views within SampleTrack<sup>™</sup> allow the user to define various filters. The user specific filters can be stored with a user defined name and are available after each new log-in. To open the filter definition window, press the button located to the right of the filter bar in the respective window.

A Filter and Sort Sample Order window will appear, which will allow you to define a variety of filter settings, or select an existing filter.

ilter List		Filter Settings Enable		
Name	Owner	Sample ID	= <b>•</b> AA*	
🕼 Created samples	Spectron			
🕼 Verified samples	All	Client ID	= -	-
🕼 Started samples	All			
Prepared samples	All	Tube ID	= 💌	•
Finished samples	All			
Approved samples (this month)	All	🔽 Status	= 🗾 Finished	-
Error samples	All			Ŧ
Not yet finished samples	All	mounding		
PROTON Experiments	ADMIN	Date Entry	Today	-
		Sort Settings	Availabla Fields	
		Sorting Order	Available Fields	
			Up Sample Title Project Number	
		Sorting Order	Up Sample Title Project Number Group	
		Sorting Order	Up Sample Title Project Number Group Client Company	
(		Sorting Order	Up Sample Title Project Number Group Client Company Status	
		Sorting Order	Up Sample Title Project Number Group Client Company Status Date Tube ID	
	▶ <u>S</u> et Owner	Sorting Order	Up Sample Title Project Number Group Down Client Company Status Date	
		Sorting Order	Up Sample Title Project Number Group Down Client Company Status Date Tube ID	

Figure 6.15. Example of a Filter Definition

In the example above a filter with the following functions has been defined:

All sample ID's begin with **AA**, have a status **Finished** and are sorted in Client ID and Tube ID order.

When creating a new filter you should always use a meaningful name to ease identification at a later time.

Defining these special filters makes working with SampleTrack<sup>™</sup> both efficient and easy.

Although working with the filters search function makes SampleTrack<sup>™</sup> efficient and easy, it does not eliminate the desirability of a standard search function. To ease searching for a specific Sample ID, Tube ID etc., a standard search function has been implemented in SampleTrack<sup>™</sup>.

There are two separate search possibilities, one in the Client Tool window, and one in the Sample Order Details window.

The Client Tool window has a Search in View feature that will **search through the current list of data in the display window** for a specific parameter. To perform this search, press the Search in View button located at the top of the Client Tool window.

Figure 6.16. Search in View - Client Tool Window

Search For Sample Order	×
Search Sample Order with this parame	ters
In Field	Search <u>P</u> attern
Sample ID 💌	99030948
Options	Direction
🔽 🔟 hale text only	Forward     Search
☑ <u>C</u> ase sensitive	C Backward Close

The Sample Order Details window has a Search in DB feature that will **search through the Oracle® database** for a specific parameter that you specify. To perform this search, press the Search in DB button located at the top of the Sample Order Details window.



Search For Sample Order		×
Search Sample Order with this parame	eters	
In Field	Search <u>P</u> attern	
Sample ID 💌	99030948	
Options	Direction	
✓ Whole text only	<ul> <li>Forward</li> </ul>	<u>S</u> earch
Case sensitive	C Backward	Close
10 2-11 10 10 10 10 10 10 10 10 10 10 10 10 1		Liose

In both cases select the relevant search field from a list of values in the In Field pull-down menu and enter your search parameter in the Search Pattern field. If your search is to be case sensitive, check the Case Sensitive box. You can also indicate whether you want to search in ascending or descending order by selecting Forward or Backward in the Direction box.

Press Search to begin the search operation, or close to cancel and exit the window.

If the search program reaches the end of the list (when you are searching while in the Client Tool) without finding your search parameter, you will be prompted whether you want to restart from the beginning. This is significant if you have started your search in the middle of the list. In this case, press Yes to continue, or No to end the current search.



Attention! The search function corresponds directly with your selected filters. Therefore, you will not find information outside of your filtered data.

# Interaction with Analytical Applications

7.1

#### Interaction with TOPSPIN®

SampleTrack<sup>™</sup> allows you to directly call TOPSPIN® with your associated experiment data set. This results in easy verification of measurement results and structure elucidation.

Figure 7.1. The Paperclip Icon

0

The paperclip icon to the left of the experiment symbol shows existing experiment paths.

Doubleclick an experiment marked with the paperclip icon. After selecting TOP-SPIN® SampleTrack<sup>™</sup> transfers the information about the location of the data set to the program. TOPSPIN® then converts the data and displays it on the screen.

With the View next Experiment button, the program automatically switches to the next experiment and passes the associated data to TOPSPIN®.

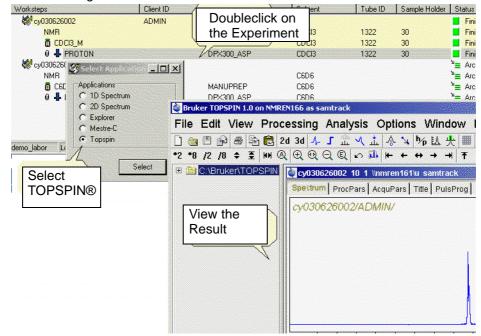


Figure 7.2. Interaction with TOPSPIN®

The result data path is stored in the database and can be seen in the experiment detail view.

Figure 7.3. The Results Data Path


xperiment - Edit Item	
General   Experiment Parameters   Results   Comments	
Result Data	
Disk Unit Partial Proc Path Dataset Name	Instrument Login Name Experiment Number
	су 10
Result Data Path	
C:\u/data/SamTrack/nmr/cy030430003/10	

#### Interaction with AMIX<sup>™</sup>

SampleTrack<sup>™</sup> allows you to directly call AMIX<sup>™</sup> with your associated experiment data set. This results in easy verification of measurement results and structure elucidation.

The paperclip icon to the left of the experiment symbol shows existing experiment paths.

0

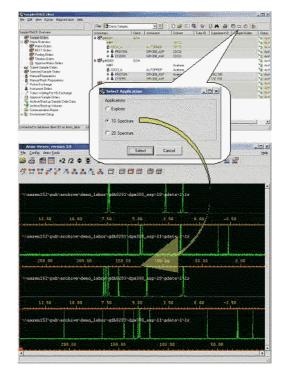

#### Figure 7.4. Interaction with AMIX™

#### Interaction with Structure Editors


7.3

SampleTrack<sup>™</sup> also allows you to call a structure editor e.g. MDL® ISIS/Draw<sup>™</sup> where you can enter the expected chemical structure results.

This feature is also activated in the Sample Order Details window by pressing the Documents button, and selecting Structure. The Structure Editor program will then be opened and the sample structure can be viewed. Usually the structure is entered when the sample order is created. The saved file will be linked with the sample order. The file name could be for example the laboratory journal number or the sample order number or a project number.

If desired a template file with customized settings can be opened as default.

Figure 7.5. MDL® ISIS/Draw<sup>™</sup> Template

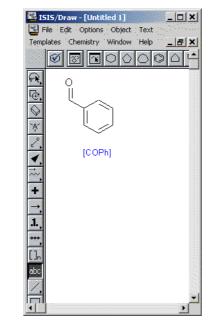

#### Interaction with other Windows® Applications

7.4

SampleTrack<sup>™</sup> allows you to link to other Microsoft Windows<sup>®</sup> applications (e.g. for reports) directly from the Sample Order Details window.

To create a new UNC (Universal Naming Convention) formatted link to an existing Microsoft Word® document, for example, open the Sample Order Details window and click on the **More** button.

Figure 7.6. The More Button.

 $( \rangle \rangle$ 

This will expand the viewing options to include a Document Files box. Open the Windows® Explorer, and drag the desired file to the Document Files box, and drop it in the box. This will create a UNC format link to the desired file.

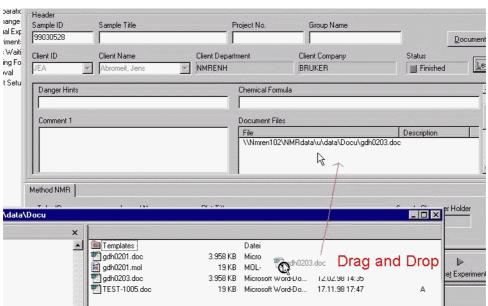


Figure 7.7. Creating a New UNC Formatted Link to a Document File

Once a new UNC formatted link has been created you must only double-click on this link to open e.g. Microsoft Word® to view the file.

#### Creating a New Report Within the Details Window

It is also possible to create a new report within the Sample Order Details window. To create a new report press the Documents button, then select New Document -Report as shown in the following figure.

Figure 7.8.	Creating a New Report
riguio r.o.	



This will open a report template file in your default word processor (e.g. Microsoft® Word). The file will automatically be saved as the pre-selected destination file.



This will work only if you have set the appropriate parameters in the Options - General Settings - Applications - Report menu in the SampleTrack<sup>™</sup> Client window (See next chapter).

7.5

)verview			•		
50455051 Tunner	Microsoft Word				
ple Orders			igen Forma <u>t</u> E <u>x</u> tra: T		::::::::::::::::::::::::::::::::::::
eparation:	Standard	and the second se	sNew Roman 💌 [12		
ual Experime		<u>a</u> (*)	1 🖻 🖻 🖉 🔊		🗐 ¶ 75% 💽 👰
eriments Is Waiting (E	+ 2 + 1 + 1 +	·	1 • 3 • 1 • 4 • 1 • 5 • 1 • 6 • 1 •	7 - 1 - 9 - 1 - 9 - 1 - 10 - 1 - 11 - 1 - 12 - 1 - 13 -	1 - 14 - 1 - 15 - 1 - 2 - 1 - 17 - 1 - 19 -
ting For HD		NMR-Report			
itSetup ດູ່			gdh0203		
÷			Günther Hiss Demo001		
. :		Result:	20110001		
. +					
م		The following	display represents		
- -		1		1	
-1-2					
1 ÷					
0 - -					
ę.					
Ę					R
12-1-1					45
- 181 – I			1		
\$				1 11	

Figure 7.9. An Example of a Report

#### **Application Management**

7.6

The definition of all applications in SampleTrack<sup>™</sup> can be found using the menu **Administration - Application management.** 

In the Application Management window, on the left side, all existing applications are displayed. To add or copy an application, use the right mouse button. On the right side of the window are the options **mixed**, **global** and **local**. These options allow you to display various kind of entries:

**Mixed** is a view of the entries in the global and local window.

Global entries are saved in the SampleTrack<sup>™</sup> database. They consist of data used by the most or all the users.

For example, the program executable path for 90% of the SampleTrack<sup>™</sup> users in one company could be:

C:\Programme\Microsoft Office\Office\WINWORD.EXE.

And the rest of the users could have a path like

C:\Program folders\Microsoft Office\Office\WINWORD.EXE.

In this example local entry is written, which is different from the global entries. The program reads the local entry first, and then the global entry.

If an additional program is used only by a few people: this program must be defined in the local viewer. These definings will be saved in the local registry of the PC.

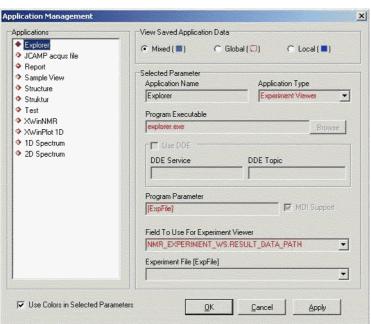


Figure 7.10. Administration - Application Management

#### Application Type

7.6.1

The Application type decides in which environment within SampleTrack<sup>™</sup> an application may be started.

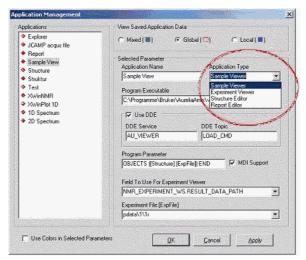


Figure 7.11. Defining the Application Type

- · Sample Viewer (one or more experiments in an order)
- Experiment Viewer (one or more experiments with different orders)
- Structure Editor (IMDL® ISIS/Draw™/MESTRE-C™ or others)
- Report Editor (Word®/Excel®)

In the **Sample Order View** the **Application Selection** window is called by simple double clicking the experiment (Experiments with the papeclip icon only!).

You can also mark one or more experiments in the **sample orders view** window and click the 'view the selected experiments' button.

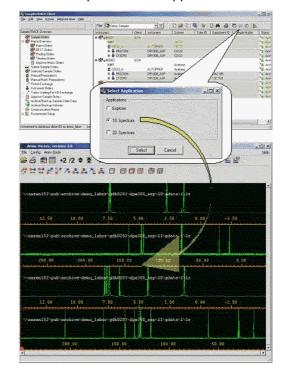


Figure 7.12. Example 1: Select Application

#### Selecting Applications in the Detail Window

7.6.3

The Applications defined as '**Experiment Viewer**' is selectable in the **order detail** window pressing the right mouse button as shown in the next figure.

Delete Item	Entf
🖸 View Experiment	Strg+L
Change Item Status	
View Item Info Configure Item Fields	
Select Application  Applications  Explore  D D Spectrum	5
C 2D Spectrum	

Figure 7.13. Example 2: Select Application

# EasyDialog

8.1

# Introduction to EasyDialog

EasyDialog is a program that provides administrator-definable forms that allow an open access user to quickly and easily add a sample order to SampleTrack<sup>™</sup>. EasyDialog is ideal for a user who needs to quickly add a routine sample order, or for an inexperienced user to easily begin productivity.

EasyDialog lacks the full features obtained by using the Add Sample Order in the client window, however, provides a productivity-enhancing feature for routine orders.

The administrator can configure EasyDialog for individual or groups, based on their specific requirements.

# **Opening EasyDialog**

To open the EasyDialog window press the Windows® Start button and select SampleTrack<sup>™</sup> - EasyDialog.

Figure 8.1. EasyDialog Icon



After a moment, a login window will appear, prompting you to login to EasyDialog.

#### EasyDialog Login Window

After you enter your login name and password, press the Login button to finish the start-up process. The EasyDialog Window will appear, similar to the one shown below. Remember that your specific dialog window may vary depending on how the administrator has configured it. If you have questions on specific features of your EasyDialog window, check with your administrator.

Note: If you have administrator privileges, you can also start EasyDialog by selecting File - Open EasyDialog from the SampleTrack<sup>™</sup> Client Tool. Refer to the EasyDialog Configuration help for details

"Opening the EasyDialog Configuration Window" on page 123.

8.2

8.3

Figure 8.2. EasyDialog Login

Figure 8.3. EasyDialog Example

# The EasyDialog Window

8.4

8.5

00001005			
99031005			
Client ID	Client Name	Client Department	Client Company
NIL	<ul> <li>Niles, Stan</li> </ul>	▼ NMRENH	BRUKER
<sup>o</sup> reparation			
C Chloroform		Tube ID	
			the second s
C DMSO			
C DMSO			
Experiments			Dink   Lenut
			Print Logout
Experiments	<u> </u>		Erint Logout
Experiments	nit D20		
Experiments 1.Experiment	nit D20		<u>Exit</u>

In the EasyDialog example above the sample order information is automatically loaded when the window is opened (this is an administrator-configurable option). To add a sample order you must only select or enter the required preparation and experiment parameters and press one of the corresponding buttons.

# Entering Data in the EasyDialog Window

When you enter the data as shown in the figure below, and then press **Save & Add**, the new sample order will be automatically added to the SampleTrack<sup>™</sup> database. A confirmation message will be displayed when this process is complete.


99031007							
Client ID Client Name NIL Niles, Stan			Client Department	Client Company			
		-	NMRENH	BRUKER			
Preparation							
Chloroform			Tube ID				
			1234				
			1234				
C DMSO			11234				
			1234				
			1234	Print	Logout		
Experiments	<u> </u>			Print	<u>L</u> ogout <u>E</u> xit		
Experiments 1.Experiment				Print ave & Add			
Experiments 1.Experiment PROTON							

Figure 8.4. Entering Data with EasyDialog

#### EasyDialog Confirmation Message

8.6

The EasyDialog window will remain open after the confirmation of a data entry so that you can add additional sample orders. In this example if you wanted to add only the one sample order you could press Save & Logout, which would add the order to the database, but would then exit EasyDialog.

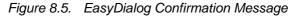



The above example is just one of many different features that an administrator can add to EasyDialog. EasyDialog can be configured for your own unique group or individual laboratory requirements. The EasyDialog field **tube ID** can be automatically filled out by the barcode reader from B-BCEThe tube ID is inserted automatically by pressing the green button on the B-BCE.

Figure 8.6. Automatic Reading of the Barcode Tube ID



# EasyDialog

# SampleTrack™ Web Client

#### WEB Client - Introduction

The **SampleTrack Web Client** allows connection to the SampleTrack database with a standard HTML browser from any PC, that has connection to the Internet or company Intranet – without additional installation of the SampleTrack Client software.

Its functionality does not cover the whole SampleTrack spectrum but is sufficient for most standard measurements. All information for a sample order can be queried and with the **EasyWeb Dialog** tool orders can be initiated the same way as EasyDialog does.

Within the **SampleTrack Web Client** services are available for structure and spectra viewing or as a quick evaluation tool. For a more intensive evaluation, an application wizard is available, that can call TOPSPIN®, AMIX<sup>™</sup> or other used applications directly.

SampleTrack Web Client is optimized for Microsoft® Internet Explorer version 5.0 or higher, but is compatible with common browsers like Netscape®, Mozilla<sup>™</sup> or Opera®.

Although Java simplifies the HTML handling, the SampleTrack Web Client does not use any Java applets.

All user and group profiles, group filters and views that were previously defined with the SampleTrack administration tool are integrated within the web tool. An additional configuration is not necessary.

All data sent through the Web can be secured by a secure connection with SSL.



For the SampleTrack Web Client installation see *"Installation" on page 225* 

#### Viewing Sample Orders

To view sample orders, login into the **SampleTrack Web Client** using the SampleTrack user name and password:

Figure 9.1. SampleTrack Web Client Login

BRUKER BRUSHIN Sample Frack Microsoft Intern	t Explorer zur Verfügung gestellt von Bruker IT- Dept.	E 🗖 🔀
Bolei Bearbeitor Greicht Eskenten Edites 2	· O · B Z S / *** 0 0+4 0+9	
Mintelli 🕘 Http://www.en166.cov.rech.bruker.de/sowplebrock		🛩 🛐 Wechseln zu
CHIOSPIN Sample K.	BRUKER	
Navigation + Login Procedure		
Contraction on Alexandria on Ale		
		and the second second second
	You are carrently not logged in to the RampleTrack System.	
	Please login:	
	Username: Jahnn	
	Password:	
	Logn	
1073		
E Funta		bitareat

In the left frame the following views are available:

- Sample Orders: Complete orders within the SampleTrack system.
- Instrument orders: Order information related to an instrument
- Matrix orders: Order information about matrix orders like BEST
- EasyWeb Dialog

Figure 9.2. Navigation Board

	SampleTrack - Lounge and Hews - Microsoft Internet Explorer sur Verfügung gestellt von Bruker IT. Dept.	
gate pearbonan Ara	di (sentes Less ) 0 • 0 − 3 2 0 / ± 4 0 0 + 5 0 • 5	8
Advanta 🕲 Http://www.	06 Instructulari dej teorglettaci	😁 💽 Wetherin to
CAROSPEN *	smith .	â
Nevigation	1 Lounge and News	
Views • Looppe and News • Sample Orders • Instrument Orders • Matrix Orders	Welcome	
Fasy Dislags + standard + standard = th 000	to	
Hiscalianeous • Druker Homopoos • ST SOAP Interface • Inter • Inter • Change Paceword • Logout	SampleTRACK <sup>TM</sup>	
laggeliten Admostation Gewengern deeldi te deenjates	Sobielle: Init Weithda	
Time Job Peace 8 1000 Michael - Lander Michael - Lander	o 1904 - sadak 10. milak **** ayos sent way	• MMM

#### Sample Orders – View

9.2.1

The **Sample Order View** lets you track the orders and view their results. You can select your favourite view by selecting the various filters in the upper window. These individually customizable selections let you view the fields that are of interest. The other selection window, which also has customizable filters, lets you select only the records that are of interest – personal orders, orders of a department or orders with a predifined status. Working with these filters and views are identical as those found in the SampleTrack client.

The number of records that can be displayed in one window can be entered in the upper window. Navigation through the pages is possible using the links "First, Prev, Next, Last". A link "Print this page" navigates to a printer friendly view.

	icht Eaueriten Kates 2	0.0		4.00	· & 🖬 ·	-			
Carlos ( Handar)	164. ner sen hubruken, dej som plebror	Storpel/eur	A STATE OF A STATE OF					💌 🛐 Wechse	
CAROSPER *	Sample &		1		BRUK	<u> </u>			
Enripelies	4 Sample Orders	Strate Contractor	and Party and a second	COLORING.					
Sears - Lounge and News - Sample Orders - Instrument Orders - Matrix Orders	n Haur 110 Nor Ver Agreend Samples an view Standard and Jampies per sage 12 - Amerikan Samples per sage 12 - Amerikan Sample								
ny Diskogs	and the second second second second	rat offrey Nexto La					2000/02/02/02/02		
Standard Standard with 020	Werksteps	El El	0 Independent	Solved	Tide ID	Start Date Ewd Date	Status Project Numbs	040707001/JEA/	
iscellaneous	A DICTOROUS				4454	7/8/2004 2:04:01 PM	Cueved		
Scelleneous Envirences	8 SEST_EVEN.gcp		01.50N	CDCR	4454	7/23/2004 2:14 34 FM	d Output		
ST SOAP arbertace		PROTON	SPECT	CDCID	4154		W vertice		
Help Info		Citron	SPECT	CDCIT	4154		Fr vented		
Change Pageword	We 040715003	ADMIN				7/13/2004 12:58:45 PM	Burrown St/Weerzere	010715003.Aprenation (benze	
Logist	CINHR .				1234	7/15/2004 12:58:45 PM	P Running		
	8 coco w		NUMEREP	coda	1234	7/15/2004 12:58:45 PM 7/15/2004 12:58:45 PM	Finished		
		PROTON	CP9G00_ASP188	CECO	1234		W Vermed		
		CLISCED 4	DPXD00_ASP108	0000	1234		64 Verified		
rapped in co	W 0+0715005	ADADA				7/15/2004 1:05:38 PM	P Running	040715003/ADM01/	
	CI NMR				2225	7/15/2004 1:08:39 FM	P Running		
Invested to down53 or lives. Salar	COCI3_M		MANUPTER	COCIS	2225	7/15/2004 1:05:39 PM 7/15/2004 1:05:39 PM	Finished		
122012201210000000000000000000000000000		PHOTON 4	DPX300_ASP	CDCD	2221		Queutd		
		C13CF0	DEX100_ASP	COCIS	2223		Queutd		
	<b>#</b> # 040723005	er .					Created	040724003/CY/	
	CI NHR			162262			Created		
	CDC3 Preparat		NUMPREP	COCID			Created		
		PROTON 4	DP9Q00_ASP	CDCD		Company of the second second second second second	Created		
	# 040722001	EA				7/22/2004 3(56)14 PM 117/22/2004 3(56)14 PM	Russing Lonvoerzere	040722001/JEA/Ethylbenzene	
	0 CDCD M		HMAPSEP	ceas		17/22/2004 3:49:19 PM 7/22/2004 3:49:19 PM	Friend		
		penno 🗳	DPICTOR ASP INTE		04072200		Comment.		
	A 0+0010001	MIL			- ALEN	a second s	d Created	040910001.9/08/	
	CINK				ST100121		Created		
	S INN CLEAN W	ater and	AUTOFREP	UNRIÓN	NST100123		a created		
	ABC123+5749	HEE					# verified	ABC12345743/MBE/	
	6 default Urine to	et nov oip	GILSON, MOE	Buffer pH	17		fil verified		
	0 0000 040	10333510002010	OPX100 REST	Define pil	0035306		W verified		

Figure 9.3. Sample Orders - View


The worksteps are presented in a tree view, as is within the SampleTrack client. By clicking on the sample ID, the method, the preparation or the experiment the program navigates to a window with more detailed information. In the **Detailed Sample Order** window information can be found on the complete order, e.g. methods, attached documents or comments.


#### Figure 9.4. Complete Order

				10		and the second		
Sample Order Details								
		View: Standa	ard View 💌				Close wind	
ample Information: > Advance	ed Information							
Status	₩ Verified							
Sample ID	A8C123#5749							
Sample Title	ABC123#5749/M	BE/						
Client ID	MBE							
Docu files								
no files available.								
lethods:								
BEST-NMR-2 W BEST	-NMR							
Tube ID: Journal Nun	nber: P	lot Title:	Sample Chang	er Holder:		Status & Veri		
/orksteps:								
Vorksteps	are por tester de la casa de	Instrument	Solvent	User ID	Start Date	End Date	Status	
default_Urine test mbe.gsp		GILSON_MBE	Buffer pH7	MBE			Werified	
CDCl3 stan		DPX300_BEST	Buffer pH7	MBE			Werified	
	PROTON	DPX300_BEST	Buffer pH7	MBE			Verifi	

The SampleTrack Web client provides structure and spectra viewing capabilities, including zooming in and out. Zoom In - Zoom Out

To view distinct areas of spectra click at the starting point, choose the area to be zoomed in (without clicking), and click to confirm. The area selected is also shown in the printer friendly view.

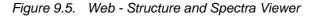

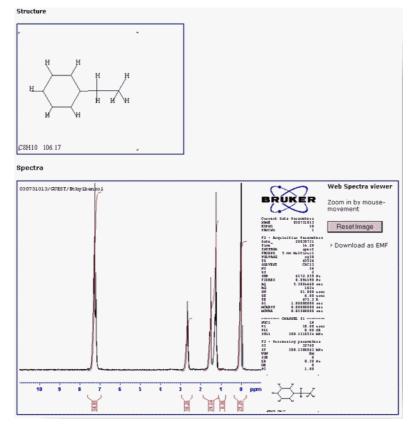

Click the upper left corner once and afterwards click the lower right corner to zoom in.


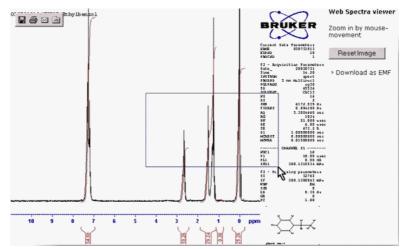


Figure 9.6. Web - Structure and Spectra Viewer, Zooming In

To review the original size just click the **Reset Image** button

Figure 9.7. Web - Structure and Spectra Viewer, Reset Image

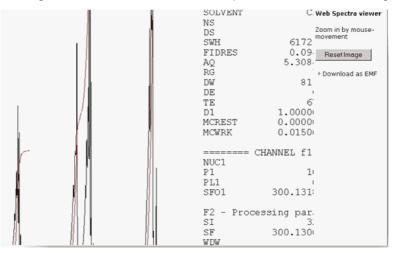


After finishing an experiment, the **Experiment Detail** window shows the spectra results, when XWIN-PLOT<sup>™</sup> has been triggered to generate the results as emffile.

The link "Search" provides the possibility of performing a manual search, e.g. for a sample ID. Select the kind of field information to search, such as sample ID, status or a date and the search criteria.

When "Whole text only" is checked, the program searches for an exact match.

When "Whole text only" is unchecked, the program looks for patterns that include the entered words, e.g. "BCD" finds "ABCD123".

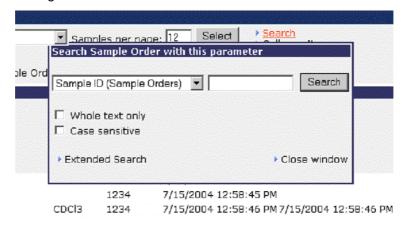

#### Figure 9.8. Search

#### **Extended Search**

9.2.4

With the link "Extended Search" the Web module (like SampleTrack client) provides more complexe queries over several database fields. Remember, that you can only find records within your predefined datapool, that were defined as your group filter. Changes here can only be made by the administrator in the Sample-Track client.



Datei Bearbeiten Ansicht	and the state of the	and the second second		n @Meden 3 2-3 5							
partners and partners in a subscription of the	and the second se	Constitution of the	and the second second	Constant from the property of the state of the	ŋ •	6312 1		22523			
Adresse (a) http://nmren166.r						1000		2.22.2			
Coogle -	<u> </u>	D MeD-24	sche •	Basel boovert	ptionen	-					
	Sample mac	k.				000		₽	RUK	<u>g</u> r	
Navigation	+ Sample	Orders	1202		and.	322				i de la constante	
Views  Counge and News  Sample Orders  Matrix Orders  Matrix Orders  Standard  Standard  Matrix D20	Goto page Workster	Exten	God ded So ed Filte				Samt Copy	oles pi	er page	n 12 Sele	ct > Searc > Collap > Print t
Miscellaneous		[	র 🗖	Status (Sample Orders)			Archived on CD	-	) *	end 💌	
<ul> <li>Bruker Homepage</li> <li>ST SOAP Interface</li> </ul>	<b>8</b> 0 040	[	<b>v</b>	Mod Date (Sample Orders)			This Month	•	) •	and •	м
<ul> <li>Help</li> <li>Info</li> </ul>	🛄 N	(	<b>v</b> .	Please Select					) .	and 💌	м
<ul> <li>Change Password</li> <li>Logout</li> </ul>		[	·г	Please Select			[		) •	and 💌	M7/
	§	[	Ψr	Please Select			-		) •	and T	
	8 04071	(	- r	Please Select	-	•			) -	and 💌	
Logged in as	🗋 N	(	-	Please Select					) •	and 💌	7/
Administrator	I CREAKE	-		Disease Salard						and w	

#### Instrument Order View

The **Instrument Order View** lets you view experiments that are to be measured in a certain instrument. The filters here are also customizable.


Figure 9.10. Instrument Orders

Retei Brerbeten Art			and the second		P * * * @ P+1	Salta Managara Salta ang			w 🛃 Wedler
Cooge-		Sucha - Ø B							
	sample k.	1111		1	BRUKE				
levisories	Fiber: 06 Not Ye		les 🕑 view	Standard	Instrument. DPX	Max. nest 9	Experiments Splect		* Search • Prizz this page
Sangla Criders     Sangla C	Сталист В расток Ф в касток Ф в сазова Ф в касток Ф в расток Ф в расток Ф	245025010 246073002 046713605 046713605 046723805 540723805 517200123	Clinet 10 Cr Abrets Abrets Cr MBE	Tebe (I). Samuk 2221 2231	Holder Geleni CDC3 CDC3 CDC3 CDC3 CDC3 CDC3 CDC3 CDC	8246 0004 (0)42004 203 58 90 (0)42004 203 58 10 (0)42004 203 51 88 90 (0)42004 203 51 88 90 (0)420004 407 13 89 (0)420004 407 13 89 (0)420004 407 13 89	4 - 4 5004 - 23318 PM (9042006 23318 PM (9042006 23718 PM (9042006 4718 PM (9042006 41118 PM (9042006 41118 PM (9042006 41318 PM)	M Verified Quarter Connect Connect Connect Connect M Connect M Connect	

**Matrix Orders** 

9.2.6



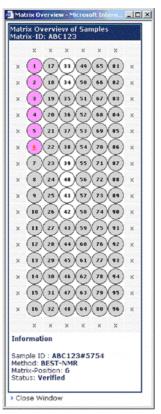

Special matrix orders, e.g. for BEST, can be found with this view.

Enter the matrix ID to find all orders with the leading matrix ID.

Using the link "Matrix-Overview" a graphical overview is displayed with a special color for each status of a matrix order.

The worksteps can be diplayed by clicking on the suborder.

# Enter Orders Using the EasyWeb Dialog

On the left navigation board one or several customized order entry forms can be found – **EasyWeb Dialogs**. Those customer interfaces are not connected to the **SampleTrack Client EasyDialog** and must be configured separately.

Figure 9.12. EasyWeb Dialog

↓ Zurück • → · ③ 🔄 付 ③ Adregse 🕘 http://nmren166.rmrenh.bru	ker.de/sampletrack/ProcessPage?first_click=10/fie-		Wechseln zu Lini
	🚯 Web-Suche 🔹 🦪 🗗 🗗 🖏	· · · · · · · · · · · · · · · · · · ·	
	le k	BRUKER	
Navigation + Star	ndard with D2O		
Views > Lounge and News > Sample Orders		Please enter your data	
Instrument Orders     Matrix Orders	Sample Order		
Easy Dialogs > Standard > Standard with D20	Sample ID * 040924003 (unique) Client ID * ADMIN	Tube ID *	
Miscellaneous Bruker Homepage ST SOAP Interface Helo	Solvents*	Experiments*	
Info     Change Password	C Chloroforme	E PROTON	
Logout	C DMSO	PROTON with D20	
	C C6D6 C MeOH	F C13	
Logged in as Administrator	1	Continue to next step	
Connected to dben152 as demo_labor			

The order can be entered on a single page or with a wizard. Only the minimal information is required, usually the method, the solvent and the experiment.


After the entry is made a confirmation window is displayed, where the order can be submitted or modified.

After the submission the new order is displayed.

9.4

# Applications (only for Windows® Systems)

The experiment results can be edited with additional applications. In principal it is not possible for security reasons - to start an application from the Web. Therefore the **SampleTrack Application-Wizard** is needed, which starts the selected program with the necessary parameters. Click the experiment Toolbox icon and select "Application" in the following menu.

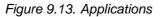



If the application wizard was not started before, a new window is displayed.

Click "Execute ST-Application-Wizard". When the program appears in the system border, click "Continue".

Figure 9.14. Application Wizard

🖉 Bruker - Install Applicaton Opener Serve 💶 🗙
Need SampleTrack Application Wizard
You need to execute the Sample-Track Application Wizard. Please
First: > Execute ST-Application-Wizard Second: > Continue
For the first installation use following settings: IP-Address of Server: 149.236.16.166 Port: 45222
> Cancel

To view a spectra, select one of the viewers from the **View Spectra** window. The program will open and display the spectra in a new window.

Figure 9.15. The View Spectra Window

Please select the program for showing the spectra:
Amix Viewer
C TopSpin
C Explore
Show Spectra Cancel

# Interaction with Automation Robots

SampleRail<sup>™</sup>

SampleRail<sup>™</sup> is part of a system under the management of SampleTrack<sup>™</sup>, that automatically prepares an NMR probe, inserts it into an NMR magnet, performs NMR experiments on the probe, and transports it back to the preparation system. The SampleRail<sup>™</sup> fulfills the transporting task from the preparation system into the NMR magnet and back while a TECAN® Genesis pipetting robot is used as the preparation system. The SampleRail<sup>™</sup> may also be used in a 'manual insertion mode' to insert/ eject single NMR probes manually into the NMR magnets.

#### **TECAN® - Pooling Module**

The Pooling Module is designed to create special sample orders. These orders contain several worksteps for automation using well plates. The first workstep is to create a preparation of different well plate substances in one tube. The second workstep encompasses the measurement of these substances. After finishing the measurements, the results are interpreted and new sample orders are generated, dependant on the results of the interpretation. Each generation of a new sample order is part of a level. Levels are used to visualize the progress of the automation process. Using the Pooling Module it is very easy to handle large amounts of samples and well plates. The Pooling Module is initially designed for protein screening but can also be used for other applications.

The SampleRail<sup>™</sup> is part of a system under the management of the Sample-Track<sup>™</sup> Laboratory Information Management System and ICON-NMR<sup>™</sup> automation software, that provides automatic preparation and mixing of multiple compounds, and subsequent NMR measurements.

A typical application example is the protein-ligand binding detection by NMR. This system automatically prepares an NMR sample, inserts it into an NMR magnet, performs NMR experiments on the sample, and transports it back to the preparation system.

The SampleRail<sup>™</sup> fulfills the transporting tasks from the preparation system into the NMR magnet and back, while a TECAN® GENESIS<sup>™</sup> pipetting robot is used as the preparation system. The SampleRail<sup>™</sup> may also be used in a **Manual Insertion Mode** to insert and eject single NMR probes manually into the NMR magnets.

10.1

10

#### **TECAN®** Requirements

SampleTrack version 2.35 standard installation.

Stmatrix.bpl and Stpooling.bpl must be installed.

Required SQL tables MATRIX\_TABLE and MATRIX\_ITEMS must be installed. These are installed automatically beginning with DB release V2.039.

Stmatrix.bpl and Stpooling.bpl must be loaded and configured.

Script library PoolingLib V5.x must be installed.

If TECAN® system is present, thescript library TecanLib V5.x must be installed.

Preparation instruments must support the pooling option.

#### **TECAN®** Plugin Installation

10.2.2

Matrix and pooling support are activated using the Configure Plugins window. To open the Configure Plugins window select: Menu->Administration->Configure Plugins. In this window you can add the necessary Stmatrix.bpl and Stpooling.bpl Plugins to SampleTrack Server and Client.

Figure 10.1. Install PlugIns

gIn Overview for SampleTRACK Version 2.44	
art Plugins   Local Packages&Plugins	
Server Plugins	Available Plugins
SampleTRACK Matrix Support Package SampleTRACK Proving Package	SampleTRACK Archiving Package SampleTRACK Barcode Prinhing Package SampleTRACK BEST Package SampleTRACK Easy Dialog Package SampleTRACK Easy Dialog Package SampleTRACK Marks Support Package
Client Plugins	SampleTBACK Pooling Package SampleTBACK Titration Package
SampleTRACK Archeirop Package SampleTRACK Matrix Support Package SampleTRACK BEST Package SampleTRACK Pooling Package SampleTRACK Tradin Package SampleTRACK Say Diabag Package Cathol Tradin	SampleTRACK Web Configuration Package
Easy Dialog Plugins	
Up Down	1
	Ok Cancel

Once the PlugIns have been installed, they will need to be configured. This involves two steps:

- Completing the matrix settings
- Completing the pooling settings

#### **TECAN® Matrix Settings**

10.2.3

The first step is to complete the matrix settings. To do this open the Menu->Administration->Options window and select the Matrix Settings tab. Within this tab sheet you can add (or delete) available matrix configurations, by assigning a name and the well plate design, and pressing the Add Matrix Type button. In this tab sheet you should add all the matrix types that will need to be available later for pooling. Matrix type names can be in various forms, based on any established naming conventions.

The following example is based on the classic format used in BEST-NMR:

WH12=hx12

Whereas:

W is for Well plate,

H is a size from A to H, and,

12 is a size from 1 to 12.

WH12 is the name assigned to the configuration

= is a separator

hx12 is the current well plate design

In most laboratories, libraries with well plates are available for protein screening. In order to take advantage of these libraries, an import interface has been designed to import these libraries automatically into the matrix database. To import a library, select an import file and check the 'Enabled' checkbox, which activates the automatic import for the communication server. This information will then be displayed in the information section of the server.

**TECAN®** Pooling Settings

10.2.4

The second step of the configuration process involves setting the default values for pooling. To do this open the Menu->Administration->Options window and select the Pooling Settings tab sheet. The default settings include the instruments and matrix settings. These values will remain fixed based on the actual instrument configuration. Depending on the kind of matrix type, different matrix organization can be useful. The matrix organization represents from which point to which point each well should be counted to within the well plate. The Pooling Volumes and Pooling Size are flexible values, and can be changed as required.

Figure 10.2. Pooling default values

nectivity Database Archiving Options	Matrix Settings   BEST Options   Pooling Opt	ions Titration Options Easy Dialog Options	
Addite Settings           Matrix Type           State           Matrix Type           Matrix Ty	Destination Matrix Settings Matrix Type [PRUKE_TECAN # [Caparication # 2 # 3 # 4 # 5 # 5 # 5 # 5 # 5 # 5 # 5 # 5	Instruments Preparation Instrument IECAN Spectrometer Instrument DPc000_SAMPLER Pooling Pooling Size 2 Fooling Size 2 Fooling Size Complete Differentien Sorpt Name GER_Sample Differentien Fiel Name Browne Browne	

In most laboratories, libraries with well plates are available for protein screening. In order to take advantage of these libraries, an import interface has been designed to import these libraries automatically into the matrix database. To import a library, select an import file and check the 'Enabled' checkbox, which activates the automatic import for the communication server. This information will then be displayed in the information section of the server.

When you check the 'Enabled' checkbox the automation module in the server is activated. This means that new samples are generated for all samples with hits. This data will be displayed in the beginning of the information section.

#### **TECAN®** Instrument Setup

10.2.5

10.3.1

Scripts must be installed and configured for both preparation and experiment instruments. With scripts, the Sample ID is the Matrix ID with a sequence. For file imports and exports the following scripts must be configured:

Instrument	Action	Scripts
Prep Instrument	Export	ASP_EXPORT
		Pooling_Export
Prep Instrument	Import	ASP_IMPORT
		TECAN_WORKAROUND (if prep instrument is
		TECAN system)
Exp Instrument	Export	AVANCE_Export
		Pooling_Export
Exp Instrument	Import	AVANCE_Import
LIMS Instrument	Import	GEN_SAMPLE_ID
		LIMS_Import_SO_Pooling_Defaults
		LIMS_Import_SO_Pooling

The script GEN\_SAMPLE\_ID creates an unique identifier for samples, with the MATRIX\_ID as the prefix and MATRIX\_SEQ as the sequence.

Protein Screening / Module Pooling	10.3
------------------------------------	------

#### **TECAN®** Generating Samples

The dialog for sample setup is started when you press the 'New' button in the Pooling Orders Window. Here you can enter the well plate for pooling, and select the wells that are to be used by clicking on them with your mouse. After entering the Project Title, Sample Title and Experiment Name, press the Generate Order button to generate the sample orders. Worksteps with tubes will be created based on the pooling size. This level is also referred to as level 0. If Matrix ID specific in-

formation is imported from the communication server, the Ligand's are displayed in the Ligand list in the table BSTS\_REGISTRY.


- 0 > na Mai Matrix ID POOL124 • GEN Sample ID From ▼ Ad ▼ B4 TECAN • Field Po A01 A02 A03 A04 A05 A06 A07 A08 A09 A10 A11 A12 R01 ◀ Exp Solv) ent Name 1 ent Name • Generate Orders Close


Figure 10.3. Protein Screening / Module 'Pooling'

**TECAN® Starting the Automation Process** 

10.3.2

Before a new automation process can be started, all currently running automation processes must be stopped. This means the Tecan® Genesis System must be reset to its initialization state, and ICON-NMR<sup>™</sup> must be reset. Following this step the Pooling Setup window will be closed and the updated sample list for this Matrix ID should be visible.

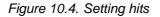

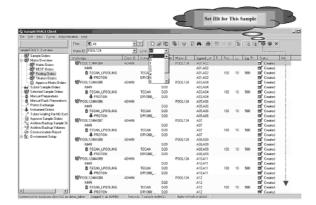

Press **Select All** or **CTRL A** and submit all the samples. The initial level is 0 with 4 substances. At this point all of the samples are ready for automation. Once all the samples have been submitted, the worksteps will be run automatically.

After the experiments are finished, the data can be viewed with the available spectra viewer and a decision can be made if the samples are hits or not. If the samples are hits, then the corresponding checkboxes can be selected for these

samples. The next step (level), with 2 substances per sample, can be executed by pressing the Execute button. Protein Imaging is finished when all the samples with hits have been measured with only one substance.

# Environment Setup

SampleTrack<sup>™</sup> Environment Setup

The SampleTrack<sup>™</sup> Environment Setup module of the SampleTrack<sup>™</sup> Overview allows existing information from the user, instruments etc., to be defined or modified. To avoid unwanted deletions or modifications to the setup parameters, access to this section should be limited to the System Administrator.

#### Inserting a New User in the Companies Setup Module 11.2

The Companies setup module allows you to insert new users in the Sample-Track<sup>™</sup> system. This process can basically be broken down into three steps:

Step One: Inserting a New Company" Inserting a New Company" on page 86

Step Two: Entering a New Department" Entering a New Department" on page <u>87</u>

Step Three: Entering a New User" entering a New User" on page 88

11.1

# Inserting a New Company

As some laboratories have users from various companies, the first step is to determine if the new user's company exists on the **Companies** list. When not, a new company entry can be added to the list:

- Highlight **Companies** in the Environment Setup.
- Press the **New** button.
- Fill out the data blocks.
- Press **Save** to enter the new information in the list.

<mark>∛Sample Track Client Tool - [Compar</mark> ♥ Eile Edit View Options Window <u>F</u>			
SampleTRACK Overview			Beload from DB
Sample Orders Verify Sample Orders Manual Preparations Protein Kachenge Frish Maruad Experiments Experiments Wating (Delayed) Tubes Wating for HD Exchange Order Appioval Companies Solvents Proten Exchange Solvents Solvent Mathods Experiment Types Experiment Substitution for In Instruments	Companies Company DEMO_COMPANY UNKNDWN	Basic Information       Company Short ID     Company Name       DEM0_COMPANY     Demo Company       Description     Demo Company       Address Info     Street       Demo I4     ZIP Code       Z1P Code     City       [2345]     Demo City       Country     Demo Land       State     Phone       Phone     Fax       [300 000 0000     [300 000 0001]	
prinected to database wgs_rmren102_orcl a	s sampletrackuser.	Logged in as ADMIN. Auto refresh enabled.	

Figure 11.1. Environment Setup

# Entering a New Department

11.4

The next step in inserting a new user is to determine if the new user's department exists on the **Departments** list. As before, if the Department doesn't exist a new department can be added to the list:

- High-light **Departments** in the Environment Setup.
- Press the **New** button.
- Fill out the data blocks.
- Press **Save** to enter the new information in the list.

Sample TRACK Overview         Image: Sample Orders         Image: Sample Order Sample Order Approval         Image: Sample Order Approval	
<ul> <li>→ DEMD_COMPANY</li> <li>→ DEMD_DEPARTMENT</li> <li>→ TEST_DEP</li> <li>→ - UNKNDWN</li> <li>→ TEST_DEP</li> <li>→ Departments</li> <li>→ Departments</li> <li>→ Departments</li> <li>→ Demo_Departments</li> <li>→ Demo_Departm</li></ul>	Demo Depatmeni     Company Short ID
Contraction Contra	Cty Cty Fax 000 000 0001 New Deete Cancel Save
Connected to database wgs: miren102 orcl as sampletreckuser. Logged in as ADMIN.	

Figure 11.2. New Department

# Entering a New User

Once you have selected the new users company and department, enter the new user:

- High-light **Users** in the Environment Setup.
- Press the **New** button.
- Fill out the data blocks.
- Press Save to enter the new information in the list.



Note: In each of the cases mentioned above you can speed the entry process by finding a company, department or user with similar data and using Copy to create a new entry, whereas you must only edit the data. To insert a new entry DO NOT simply select a present user and edit the entry, as this will only edit that entry, not create a new one. Always use **New** or **Copy** to create an entry!

#### Figure 11.3. New User


Sample Orders       Users         Gr Verity Sample Orders       ADMIN         Proton Exchange       GUEST         Besic Information       User Short ID         Guest Speriments       GUEST         Last Name       First Name         Guest User       Guest         Order Approval       Users         Companies       Guest         Companies       Guest User         Proton Exchange Solvents       Departments         Solvents       Guest Spectrometer User         Proton Exchange Solvents       Guest Spectrometer User         Proton Exchange Solvents       First Name         Proton Exchange Solvents       Guest User         Proton Exchange Solvents       Guest Spectrometer User         Proton Exchange Solvents       First Name         Proton Exchange Solvents       Guest User         Proton Exchange Solvents       First Visible         Prone/Fax       Phone         Phon	Exit
Copy     New     Delete     Cancel     Save       Set to Visible     Set to Invisible     Set to Invisible     Set to Invisible       Connected to database was immenia     Connected to database was immenia     Auto refresh enabled.	

# Deleting a User, Department, or Company

A company can only be deleted when all of the departments have been deleted. In turn, a department can only be deleted when all users within the department have been deleted. In any case discretion should be used in who has access to this section. Generally, only the system administrative should be assigned this privilege.

To delete an entry, highlight the entry with your mouse and press **Delete**. You will prompted for a confirmation, press Yes if you are sure you want to delete the entry.

Figure 11.4. Message: Delete?

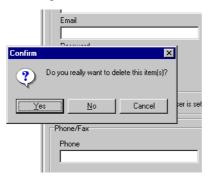

## Inserting a New Instrument

11.7

11.8

11.6

Important: The insertion of new instruments should normally be carried out only by trained Bruker personnel. An improper insertion or modification of an instrument can result in a communication failure.

A new instrument is added to the **Instruments** list by pressing the New button and completing the necessary entries for **Common Settings**, **Communication Paths & Files** and **Spectrometer Settings**. The new instrument is saved to the list by pressing the **New** button.

Again a word of warning, incorrect parameters will result in improper or no communication!

# Inserting a New Method Type

The **Method Type** serves the purpose of providing a short description of the various type of preparation methods that are available. As new methods are supported, they can also be described here.

#### Inserting a New Solvent

In most cases solvents and experiments are maintained within the spectrometer software, e.g. ICON-NMR<sup>™</sup> or external LIMS. They are imported automatically into the SampleTrack<sup>™</sup> system.

The **Solvents** block under NMR provides information on the instruments and their solvents that are currently defined. To insert a new solvent:

- Highlight Solvents.
- Select the appropriate instrument (e.g. DPX300) from the **Instruments & Solvents** block.
- Press **New** and fill in the required data.
- Press **Save** to add the new solvent entry to the instrument that you have selected.

#### Solvent is Visible

When entering a new solvent you will notice two checkboxes, **Solvent is Visible** and **Solvent is Valid**. At times it may be desirable to make a solvent temporarily unavailable to a user. Deselecting the **Solvent is Visible** box will hide this option from the user, the user will not be able to select the solvent. Checking the box will return the solvent to the list of options.

#### Solvent is Valid

When a solvent will no longer be used for experiments you can not simply delete it, as previous experiments may still have a reference link to the solvent. Instead the **Solvent is Valid** box should be deselected, which will delete the solvent from the list of solvents that are available. The solvent can be reinstated at a later time by rechecking the box.

Image: Second Stress	
SampleTRACK Overview Sample Orders Verify Sample Orders Manual Preparations Proton Exchange Finish Manual Experiments Delay Experiments Experiments Waiting For HD Exchange Order Approval Experiments Setup Companies Delay Experiments Delay Experim	I Exit Reload from DB trument Short ID X300

Figure 11.5. New Solvent

# Inserting a Proton Exchange Solvent

The solvents listed in this module represent all of the proton exchange solvents that have been defined for manual preparations (only).

To insert a new solvent entry:

Highlight Proton Exchange Solvents.

Press **New** and enter the requested data.

Press Save to add the new solvent to the list.

<b>W Sample Track Client Tool - [Proton Exchan</b> <b>B</b> rile Edit View Options Window Help	ange Solvents]			_ 🗆 ×
				<u>I</u> Exit
SampleTRACK Overview	5 1 01 1			Reload from DB
Verify Sample Orders	Description Descr	ange Solvent Short ID In:	ANUPREP	
Connected to database wgs_nmren102_orcl as sample	letrackuser. Logged in as ADMIN.	Aut	to refresh enabled.	li.

Figure 11.6. New Proton Exchange Solvent

11.10

# Inserting a New Preparation Method

This module allows you to pre-define new preparation methods. To configure a new preparation:

#### Highlight Preparations.

Select the relevant instrument.

Press the New button.

Fill out the requested data.

Press Save to add the preparation method to the list.

#### Preparation is Visible, Preparation is Valid

As was true with the insertion of a solvent module, the **Preparation**, **Solvent Methods**, and **Experiments** modules also have **Visible and Valid** checkboxes.

The same rules apply for those checkboxes as those for the solvent module.

Sample Track Client Tool - [Preparation File Edit View Options Window Help		×
SampleTRACK Overview		<u>R</u> eload from DB
Verify Sample Orders     Manual Preparations     Proton Exchange     Frish Marual Experiments     Delay Experiments     Delay Experiments Waiting (Delayed)     Tubes Waiting For HD Exchange     Order Approval     Companies     Operatments     Delay Experiment Substitution for In     Preparations     Experiment Substitution for In     Experiment Substitution for In     Instruments	Autopreparations	Basic Information         Preparation Short ID         ACETON_M         Instrument Short ID         MANUPREP         Description         Manual Aceton preparation         Preparation Type         MANUPREPTYPE         Preparation Information         Sample Amount         Solvent Short ID         Solvent Short Information         Sample Amount         Min. Temp.         Max. Temp         Max         Temp         Manual Aceton preparation         Preparation Information         Sample Amount         Min. Temp.         Max. Temp         Max         Solvent Short ID         Solvent Amount         Solvent Short ID         Solvent Amount         Solvent Mix Temp.         Solvent Mix Temp
Connected to database wgs_nmren102_orcl as s	sampletrackuser.	Logged in as ADMIN. Auto refresh enabled.

Figure 11.7. New Preparation Method

# Inserting a Solvent Mixing Method

### 11.12

The **Solvent Methods** section under **Preparations** allows you to configure solvent mixing methods for a preparation. Again this is accomplished by highlighting **Solvent Methods**, selecting the appropriate instrument, pressing New and filling out the request data blocks. The entry is added to the list by pressing the Save button.

To do this:

- Highlight Solvent Methods.
- Select the appropriate instrument.
- Press **New** and fill in the required data.
- Press **Save** to add the new solvent mixing method.

Sample Track Client Tool - [Solvent Elle Edit View Options Window H					
					<u>Exit</u>
SampleTRACK Overview					<u>R</u> eload from DB
Sample Orders Verify Sample Orders Proton Exchange Finish Manual Experiments Delay Experiments Experiments Waiting (Delayed) Tubes Waiting For HD Exchange Order Approval Companies Companies Companies Companies Solvents NMR Solvents Proton Exchange Solvents Proton Exchange Solvents Experiment Types Experiment Substitution for Im Instruments	Instruments & Solvent Mixing Methods → AUTOPREP → UNKNOWN → USONIC → MANUPREP	Basic Information Preparation Short USONIC Description	D for automatic preparation	Instrument Short ID AUTOPREP	BYE
Connected to database wgs_nmren102_orcl a	s sampletrackuser. Log	ged in as ADMIN.		Auto refresh enabled.	

Figure 11.8. New Solvent Mixing Method

# Inserting New Experiments

In most cases solvents and experiments are maintained within the spectrometer software, e.g. ICON-NMR<sup>™</sup> or external LIMS. They are imported automatically into the SampleTrack<sup>™</sup> system.

When the spectrometer AU program is started, data on the experiments that are available are exported to SampleTrack<sup>TM</sup>. The **Experiments** module allows you to define new experiments for spectrometers (only). This step requires that the new experiment is recognized by the spectrometer, i.e. it must be an experiment predefined for that spectrometer.

To insert a new experiment:

Highlight Experiments.

Select the required instrument (a spectrometer in any case).

Press New and enter the appropriate parameters.

Press Save to add the entry to the list of experiments.

Sample Track Client Tool - [Experim					
🧩 Eile Edit ⊻iew Options Window Ł	1eib				
SampleTRACK Overview Sample Orders Sample Orders Shanual Preparations Proton Exchange Finish Manual Experiments Leaveriments Waiting (Delayed) Experiments Waiting For HD Exchange Order Approval Environment Setup Companies Departments Users Users NMR Leavers Solvents	Instruments & Experiments	Basic Information Experiment Short ID [C13CPD Experiment Type [13C Description [C13 exp. comp. pulse d Plot Title Description Start Status	■c. 1024 scans [14]	Instrument Short ID DPX300	<u>I</u> <u>Exit</u> <u>Reload from DB</u> ▼
Proton Exchange Solvents     Preparations     Solvent Methods     Experiment Types     Experiments     Experiments     Instruments		C Start Time 18:00	_	00:00:00	
Connected to database wgs nmren102 orcl a			ew <u>D</u> elete Visible Set to Invisib	Cancel Save	aid

Figure 11.9. New Experiment

# **Experiment Substitution for Imported Files**

11.14

In some cases it may be necessary or desirable to use a substitute name for an experiment. For example, with a C13 experiment, the problem exists that the LIMS lists the experiment as C13, while the spectrometer recognizes the experiment as C13CPD. This naming problem can be eliminated by assigning an alias through this module. This module creates, in effect, an Experiment Translation Table.

Furthermore, this module allows you to assemble a group of experiments together under one experiment name. This allows you to perform a variety of experiments under one experiment title. For example, the figure below illustrates an experiment "Test". Grouped under this experiment are two experiments, a **PROTON** and a **C13CPD**. When you create a new grouping (highlight **Experiment Substitution for Imported Files**, select an Instrument, press New) the order in which the experiments will be carried out is determined as you add new experiments (in the **Order** box). Upon saving this experiment group you must only select the experiment "Test" to carry out both the PROTON and C13CPD experiments.

#### Assembling a Group of Experiments Under One Title

Sample Track Client Tool - [Experim           W Elle         Edit         View         Options         Window         Edit		es]			× 8×
					<u> </u>
SampleTRACK Overview	Instruments & Experiment Substitution				<u>R</u> eload from DB
Sample Urders     Verify Sample Orders     Manual Preparations     Proton Exchange     Finish Manual Experiments     Delay Experiments     Delay Experiments     Verify Sample Orders     Delay Experiments     Delay Experiment Setup     Order Approval     Solvents     Departments     Departments     Departments     Solvents     Preparations     Solvent Methods     Experiment Substitution for Inr     Instruments	Instruments & Experiment Substitution	Basic Informati Substitution N NMR Order Ir 2 [ Description		Experiment Short ID C13CPD 32	Save
Connected to database wgs_nmren102_orcl a:	s sampletrackuser.	ged in as ADMIN.	A	Auto refresh enabled.	

Figure 11.10. Experiment Substitution Table

# **Environment Setup**

# *The Object Manager*

# 12

# Introduction to the Object Manager

When SampleTrack<sup>™</sup> is first installed, a set of default user groups is installed in the Object Manager. The administrator must then adjust these group client profiles (or add new groups) to meet the needs of their individual laboratory. The Ob-

files (or add new groups) to meet the needs of their individual laboratory. The Object Manager allows the administrator to quickly and easily administer the privileges and work areas of groups of users. Normally each user belongs to at least the most basic group, the **All User Group**.

This group is assigned a basic set of user privileges which allow them to perform the most common tasks in SampleTrack<sup>TM</sup>. Additional privileges can be granted to a user by assigning them to another existing group that has these privileges, by changing the rights of the entire group that they are currently in, or by creating a new group that contains the privileges that they require.

When modifying an existing group, or adding a new group, there are four main configuration areas that must be addressed. These include the Client Profile, Table Defaults, Easydialog and Filter Modules.

The Client Profile is basically a set of rights and privileges that are assigned to a group of users in regards to their use of SampleTrack<sup>™</sup> and the laboratory automation. The rights that are assigned a group are based on the needs of the group of users and the standard policies of the individual laboratory.

The Table Default settings are used to create default starting values for various field entries within the tables in SampleTrack<sup>™</sup>.

The EasyDialog configuration is the pre-defined EasyDialog configuration that the group of users will be using (see the chapter on configuring EasyDialog <u>"Opening</u> <u>the EasyDialog Configuration Window" on page 123</u>).

The Filter Modules are a selection of pre-defined or user-defined filters that are assigned to the group.

# **Opening the Object Manager**

To open the object manager you must be logged-in to the SampleTrack<sup>™</sup> Client Tool as administrator. Select Options - Object Manager from the main pull-down menu as shown in the figure below:

Figure 12.1. Menu Object Manager



The Object Manager window shown below will open. If you are opening the Object Manager for the first time, e.g. right after the installation of SampleTrack<sup>™</sup>, the window will contain default entries in the Group Tree and Content frames. These entries will need to be adjusted based on your client needs.

Figure 12.2. Object Manager Main Window

File Edit View Help         Group Tree         Content         Spectrometer User         Guests         Easy Dialog User         Recycler	🞯 Object Manager	×
Administrators     Object       Spectrometer User     Image: Constraint of the second	<u>File E</u> dit <u>V</u> iew <u>H</u> elp	
Spectrometer User     Generation User     Generation User     Generation User     Generation User     Generation     Gene	Group Tree	Content
Administrators Group, Userlevel: 99	<ul> <li>Preparation User</li> <li>Preparat</li></ul>	Image: Client Profile

## The Object Manager Main Menu

The Object Manager main menu contains three main sections: File, Edit, and View; and a Help section. The File menu section contains a Microsoft Windows® standard Printer Setup, a Print selection that allows you to print the present screen, and a Close selection which closes the Object Manager.

#### Figure 12.3. Maintaining Groups

	Edit View Help Insert Group Delete Group	
<u>File</u> <u>E</u> dit <u>V</u> iew <u>H</u> e	<u>P</u> aste	<u>V</u> iew <u>H</u> elp
Printer Setup Print Close	Insert Table <u>D</u> efaults Insert <u>C</u> lient Profile	<u>F</u> ilter Information <u>U</u> pdate View F5
	Effective <u>R</u> ights <u>M</u> anage Users	Show <u>E</u> nvironment Variables
	Group Properties	

The main menu option Edit contains a group of selections that relate directly to the administration facilities of the Object Manager. The following list is a quick description of each selection.

Insert Group:	Used to create a new user group.
Delete Group:	Used to delete a work group.
Paste:	This is a Windows® standard paste routine that can be used to paste data from one group or individual to another, or from an external source (e.g. Clipboard).
Insert Table Defaults:	Used to create default starting values for various field entries within the tables in SampleTrack™.
Insert Client Profile:	Used for establishing a new client profile.
Effective Rights:	Used to provide a quick overview of the rights that are currently assigned to an individual or group.
Manage Users:	Used to add or delete users from a group.
Group Properties:	Used to assign properties to a group (e.g. group name, description and user level).
The main menu option View	Contains selections that relate to the overview of the Object Manager window.
Filter Information:	Used to select which information will be displayed in the Object Manager window.
Update View:	Refreshes the information in the Object Manager window.
Show Environment:	Displays the current attributes that are assigned to the groups
Variables:	The present environment.

# The Group Tree and Content Sub-windows

The Group Tree and Content sub-windows function similarly to the File Manager of Microsoft Windows<sup>®</sup>. The groups are displayed in the Group Tree, and the details (or contents) of the groups are displayed in the Content window.

When SampleTrack<sup>™</sup> is first installed a number of default groups are created. As the following figure illustrates, these groups contain various sub-categories including Client Profiles, Module Sample Orders (filter information), Table Defaults and EasyDialog configurations.


 Group Tree
 Content

 Image: Administrators
 Image: Object

 Image: Object
 Im

Figure 12.4. Content Sub-Window


Figure 12.5. Default Groups

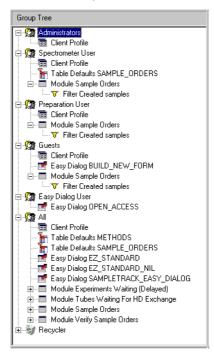

# Viewing Group Details in the Content Window

12.5

When you select one of the groups or group sub-category in the Group Tree, the contents will appear in the Content window. For example, if I select the Client Profile for group **All** the following contents will be displayed in the Content window:

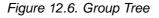

💞 Object Manager		×
<u>Eile E</u> dit <u>V</u> iew <u>H</u> elp		
Group Tree	Content	
Administrators     Spectrometer User     Preparation User     Sector Secto	Communication Report	pup Rights
•		Cancel <u>S</u> ave
		/ii


# Client Profile Example


When viewing the Client Profile, four separate window views containing various rights are available in the Content window, as shown in the figure below:

By selecting/deselecting the various rights, the administrator can quickly establish a SampleTrack<sup>™</sup> Client Profile. You will notice that each group has a client profile, where the administrator can assign specific rights for that specific group. The concept of user classification and client profiles will be discussed in further detail in the next section.

When you click on another sub-category of the Group Tree a different set of data will be displayed in the Content window. Each sub-category has its own unique details, thus the Content window will change as you move from one tree node to another.

This should be familiar for anyone who has worked with the Microsoft® Windows® File Manager.

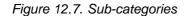

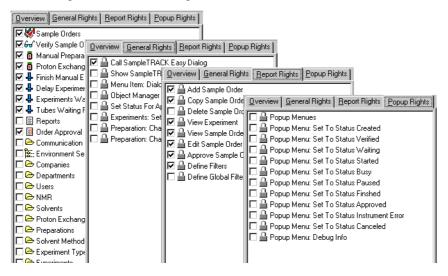

**User Classification and Client Profiles** 

12.7

Each user group has its own Client Profile, which consists of a selection of access privileges to various elements of SampleTrack<sup>™</sup>. These client access privileges are granted to all the users within a group. In most cases, a basic access configuration is established that applies to every user, then additional privileges are assigned to specific workgroups as needed. This allows the administrator to control the flow of samples to the various laboratory automation (e.g. Spectrometers) by assigning users to various groups based on their access requirements.

Each client profile can be assigned a user level. The lowest user level **00**, is the **AII** profile. The highest user level, **99** is generally reserved for the system administrator. A new group client profile is automatically assigned the next available user level number when the profile is created. This level can be adjusted later on by the administrator in the Group Properties window.

The user level privileges are inheritable, whereby a user who is part of a higher level user group (parent) inherits the privileges from any lower level group (child) that they are also a member of. For example, if I have a user "Joe Tester" that is in a group called "Spectrometer User" which has user level 04, and also is a member of the group "Preparation User" which has a user level 03, the user "Joe Tester" would have all of the privileges of the Spectrometer User group, the Preparation User group and the basic group: **All**. This will become more apparent as we discuss the details of the object manager.

### Adding a User to an Existing Group

12.8

When SampleTrack<sup>™</sup> is first installed a default configuration of pre-defined groups is loaded in the Object Manager.

A user will automatically be added to the **All** group when the user is created in the Environment Setup of the SampleTrack<sup>™</sup> client. If a user needs additional privileges, the first step is to determine if the access rights that they need are contained in an existing group. For example, if the new user needs access to reports that the Preparation User group uses, then they probably need to be added as a member to that group. If an existing group does not contain the necessary rights that a user needs, and you do not wish to add the rights that are lacking in an existing group, then you can create a new group for the user. There is no minimum requirement for the number of users that need to be in a group.

To add a user to a group, the user account must first be created in the **Environment Setup**. The user can then be inserted into a group by selecting the group with your mouse, and then selecting Edit - Manage Users, from the Object Manager main menu. This will open the group management tool for the group that you have selected. For example, if you click on the Preparation User group and open the management group tool the window shown below will appear.

Manage Grou	p: Preparation User				>
Members			Available users		
Short ID	Name		Short ID	Name	-
🕵 TESTO		<u>R</u> emove Users	🕵 TEST92		
🕵 TEST10		I	🕵 TEST93		
		<u>A</u> dd User	🕵 TEST94		
			🕵 TEST95		
			🕵 TEST96	,	
			🕵 TEST97	,	
			🕵 TEST98	,	
			🕵 TEST99	,	
			🕵 Test_stan	Tester, Joe	
			😥 UNKNOWN	Company, Unknown	_
			1		2
			OK	Cancel	

Figure 12.8. Adding a New User (Or Group of Users)

To add a new user (or group of users) to the group you must only select the individual(s) from the available users list and press the **Add User** button. This will copy the user you have selected into the members side of the window. Likewise, to delete a user, select the user from the Members window and press the **Remove Users** button.

# The Object Manager

# Example: Create a User Group

# Creating a New Group

If you do not find a user in the list of available users, then the user has probably not been created in the Environment Setup.

If the available groups do not fit the requirements of a user you have the choice of modifying an existing group configuration, or adding a new group. The choice which is made really depends on your individual laboratory requirements. For example, it may not always be desirable to give an entire group access to a specific filter if only one person or a small number of the group require the filter. In this case the individual or small group can be granted access to the filter by creating a new group with this privilege. By also keeping them as members of the original group they maintain their current rights, but also have the new privilege.

#### Adding a Client Profile

👹 Object Manager

The next step in adding a new group is to create a client profile. Select the new group that you have created. For an example I have created a group called OM\_Test as shown in figure <u>"Create a Client Profile" on page 105</u>. Select Edit - Insert Client Profile from the Object Manager main menu. A new Client Profile block will appear.

Image: Sector meter User         Image: Se		Group Tree	Content
🔚 🗁 Proton Exchange Solvents	trators meter User tion User alog User	Administrators     Administrators     Sectore and a sector and a	Qverview       General Rights       Report Rights       Popup Rights         Image: Sample Orders       Image: Sample Orders         Image: Sample Order Approval       Image: Companies         Image: Sample Orders       Image: Sample Orders         Image: Sample Order Orders       Image: Sample Order Orders         Image: Sample Order Order Orders       Image: Sample Order Orders         Image: Sample Order Order Order Order Order Order       Image: Sample Order Order Order Order Order Order Order         Image: Sample Order Order Order Order Or
Proton Exchange Solvents     Deparations     Deparations     Deparations			Preparations

Figure 13.1. Create a Client Profile

13.2

13.1

# **Example: Create a User Group**

The check boxes will all be shaded-gray when you first insert the new profile. You will need to go through each section of the Content window (Overview, General Rights, Report Rights and Popup Rights) and place a check mark in the box of the rights that you wish to assign to the new group. Likewise, if you want to disallow a right to a group, you will need to check the box so that the gray-shade disappears and a white box is displayed.


If a right has been assigned to the members of the group in a lower level (e.g. the **All** group) then you can leave the boxes gray-shaded.

When you finished making the appropriate selections, click anywhere in the group tree. You will be prompted as to whether you want to save the current client configuration, press **Yes** to save the new client profile (or **No** to exit without saving).

If you have saved the new client profile, the following window will be displayed, indicating that the changes (new client profile) will take place the next time you login to SampleTrack<sup>™</sup>:

You can continue to the next step of adding a new group without logging in. You should now see a new client profile under the new group that you have created in the **Group Tree** window.

Figure 13.2. Program Message

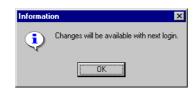

# Adding Table Defaults to the New Group

The next step in adding a new client group is to add any required table default values that apply to the group. This tool is useful if the group repeatedly uses the same values in the Sample Orders, Methods, Preparations, or Experiment windows. By creating table defaults you can place the most often used values as the default value for a given field. For example, if the new group you have created uses the same project number in the Sample Order Details table (window) you can assign a default value to the project number.

To create a new table default value, select the group you want to add the table default to (click on the group with your mouse), and select Edit - Insert Table Defaults from the Object Manager main menu. A pop-up Edit Table Defaults window will appear as shown in the figure below:

13.3

Select a table from the pull-down menu **Table** that you wish to work with. As an example select the SAMPLE\_ORDERS table, as shown in the following figure.

💥 Ubject Manager				
<u>F</u> ile <u>E</u> dit <u>V</u> iew <u>H</u> elp				
Group Tree		Content		
Administrators     Administrators     Administrators     DM_Test     Spectrometer User     Guests     Guests     All     All	Edit Table Defaults Table SAMPLE_ORDERS Column PROJECT_NUMBER Default Value OM_TEST 001	Object Client Profile	OK Cancel	

Figure 13.3. Adding Table Default Values for a Group

Select the **PROJECT\_NUMBER** field from the **Column** pull-down menu. Now enter the default value OM\_TEST 001 in the **Default Value** input window. When you press the OK button the new table default value will be listed under your new group. To view this press the **+ Box** next to your group in the Group Tree, then click once on the new **Table Defaults SAMPLE\_ORDERS** entry listed under the Client Profile. You should now see the OM\_TEST 001 value listed under the Value column of the Content window as shown in the following figure:

Figure 13.4. Defaults in the Content Window

💞 Object Manager		X
<u>F</u> ile <u>E</u> dit <u>V</u> iew <u>H</u> elp		
Group Tree	Content	
Administrators     Administrators     Administrators     Client Profile     Table Defaults SAMPLE_ORDERS     Spectrometer User     Preparation User     All     SAI     Recycler	Field Name	Value OM_TEST 001
		l li

The next time that a user logs into this group, and makes an entry in the Sample Order Details window, the new default value will for example, appear in the Project Number field as shown below:

Figure 13.5. Example: Default Project Number

💞 Sample Ord	ler Details				
	Pre <u>v</u> ious	<u>N</u> ext	Search in DB	<u>R</u> eload from DB	<u>C</u> ancel
Header Sample ID	Sample Title		Project	No. Group	
99033051				EST 001	
Client ID	Client Name		Client Department	Client Com	

In the same manner the Project No. default value will appear in the group EasyDialog Window (if the Project No. field is used in EasyDialog).

The number of default values that can be used is unlimited, however it is only advisable to use these default table values when the value will normally be repetitive.

Inserting a New Group EasyDialog	g Configuration	13.4
----------------------------------	-----------------	------

There are two methods available to assign an EasyDialog configuration to a user group. If an EasyDialog configuration already exists, such as the EZ\_STANDARD shown in the following figure, you only need to drag & drop the configuration into your new group as illustrated below (or copy and paste if you prefer).


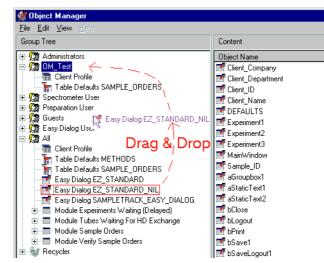


Figure 13.6. Drag and Drop EasyDialog Configuration to a Group

If an existing EasyDialog configuration does not fit the needs of the new user group that you have created, you can create a new EasyDialog configuration in the EasyDialog configuration tool and save it directly into the new group as illustrated in the next figure.

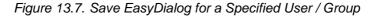

ample Order	
ave Configuration	×
Configuration	User/Group
EZ_STANDARD	All
EZ_STANDARD_NIL	All
SAMPLETRACK_EASY_DIALOG	All
C OPEN ACCESS	Easy Dialog User
BUILD_NEW_FORM	Guests
Config Name EZ_STANDARD_NIL	Save
User / Group	Cancel
2. Experiment m Easy Dialog User	<u> </u>


In either case a member of the group can then use this configuration when they log into EasyDialog.

#### Inserting a New Filter Method

13.5

In SampleTrack<sup>™</sup>, user specific filters can be stored under a user defined name and are available after each new log-in. In order to assign a new filter to a group, you must first select a module(s) for a group. The filter module is a group of related filters that are used for a specific activity, such as a sample order module, a module for experiments waiting etc.

If you want to apply the standard filters for Sample Orders, for example, you can drag & drop the Module Sample Order from the **All** group to your new group (see the figure below).

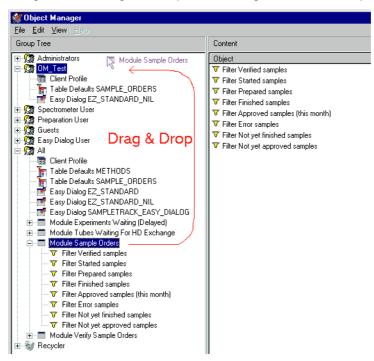



The group can then use the filters that accompany this group. You can add or delete existing filters to this module as needed. To create a new filter you must utilize the **Customize Filter Settings** in the SampleTrack<sup>™</sup> Client Tool. Refer to the section **Filter & Search Tools** for details.

#### Inserting a New Group Name

13.6

To create a new group select Edit - Insert Group from the Object Manager main menu. An input window will appear asking for a new Group Name and Description, as shown in the figure below. By default the next available user level will be displayed in the User Level field. You can either use the default value, or select a different user level from the pull-down list. If you anticipate needing intermediate user levels in the future you should select a higher user level, e.g. 10. In any case you can change this user level at a later time as needed.

When you press OK, the new group will be added to the Group Tree.

Figure 13.9. User Level in the Insert New Group Window

Insert new group	×
Group Name OM_Test Description For Object Manager Test Only	OK Cancel
User Level 05	

## Handling the Tube

The Tube ID and TubeID.txt File

The Tube ID is the temporary identifier used to identify the sample after the preparation during the NMR measurement. This Tube ID is part of the NMR method table of each sample order and is entered either manually by the preparer or during the communication, e.g. with the NMR B-ACS system.

The TubeID.txt file contains a list of all Tube ID's in use.

#### Automatic Entering of the Tube ID

During the sample preparation performed by the NMR B-ACS, the barcode collar with the associated Tube ID is taken at random from the collar dispensing station. If the NMR B-ACS receives a collar with a barcode already in use, the collar is rejected and a new one will be taken from the dispensing station. The Tube ID number is then recorded by NMR B-ACS in the preparation report file.

During the file import by the SampleTrack<sup>™</sup> software import module, the sample Tube ID is entered in the method table of the associated data set in the Oracle<sup>®</sup> database, and the TubeID.txt file is updated.

During the manual preparation of an order, the preparer must manually enter the Tube ID number.

#### Updating the Jobs in the Tube ID File

The Tube ID file is updated through the triggering of the SampleTrack<sup>™</sup> communication server by the database, which results in the scanning of all the orders in the database. The Tube ID's of all orders with header status equal or greater than **Started** and below **Approved** are written to the file.

Through this process a Tube ID of a deleted or approved order is removed from the file and can be reused.

14.4

14.2

14

### Handling the Tube ID

## System Parameters

# 15

#### SampleTrack<sup>™</sup> Client Parameters

/MultiInstance:	Allows more than one instance from SampleTrack <sup>™</sup> Cli- ent to be opened.	
/Username: <usr>:</usr>	The login user name: e.g. GUEST	
/Password: <passwrd>: Password for a corresponding user.</passwrd>		
Example 1:	C:\Programs\Bruker\SampleTrack\bin\STClient.exe / UserName:JOE /Password:test	
Example 2:	C:\Programme\Bruker\SampleTrack\bin\STClient.exe / Multiinstance /Username:admin /Password:mypass- word	

#### EasyDialog Parameters

/Username:<usr>: The login user name: e.g. GUEST /Password:<Passwrd>: Password for a corresponding user. /Admin: The program will be opened in EasyDialog configuration mode (the Toolbars appear). /Config: EasyDialog is opened with the Load Configuration window to allow EasyDialog configuration selection. /Config:<Name>: EasyDialog is opened with the <name> as a parameter. When this name does not exist the program will be halted. /SysButtons: Displays the system Windows button and system menu when EasyDialog is opened.

15.2

### System Parameters

## Shortcuts in SampleTrack™

# 16

The following shortcuts are used in most SampleTRACK windows. Some of them are not yet implemented.

- Ctrl-A Select all
- Ctrl-C Copy content
- Ctrl-D Delete method
- Ctrl-F Search
- Ctrl-I Version Info
- Ctrl-O Start Object Manager
- Ctrl-P Print screenshot
- Ctrl-R Start DB Registry Editor
- Ctrl-V Paste content
- Ctrl-X Cut content
- F1 Help
- F3 Continue search
- F5 Refresh
- F6 New record
- F7 Copy record
- F8 Save to database
- F9 Submit
- F12 Superuser Mode on / off

### Shortcuts in SampleTrack™

## SampleTrack™ Tables

#### Introduction to SampleTrack Base Tables

17.1.1

17.1

The base tables are used internally by both EasyDialog and the Object Manager. When using database version V2.033 there are four base tables used to setup a sample order:

SAMPLE\_ORDERS (header information)

METHODS (method specific header information, e.g. NMR, MS, IR ...)

NMR\_PREPARATION\_WS (NMR preparation information)

NMR\_EXPERIMENT\_WS (NMR experiment information)

The base tables are linked via primary and foreign keys as illustrated below:

SAMPLE\_ORDERS

| -----> METHOD (NMR)

| -----> NMR\_PREPARATION\_WS

|----->NMR\_EXPERIMENT\_WS

The Fields Overview section of the EasyDialog configuration window contains these very same tables and fields:

#### The EasyDialog Fields Overview

17.1.2

Refer to the chapter on EasyDialog specific help for details on the selection and use of the individual fields. see also: <u>"The Fields Overview Window" on page 127</u>

The tables that follow describe the meaning of each field found in the base tables.

The **Field** column in each table represents the name of the field as it appears in SampleTrack<sup>™</sup> (e.g. in the EasyDialog Fields Manager).

Likewise the **Description** column is a short explanation of what the field is.

#### Figure 17.1. Field Overview

Fields Overview							×
Field Overview	Fields						
	Fieldname	Value	Required by DB	Default Value	Used	Set From DB	
• METHODS	SAMPLE_FRACTION		No		No		
MMR PREPARATION WS	SAMPLE_ID	99031603	Yes		Yes		
MR EXPERIMENT WS	SAMPLE_NATURE		No		No		
	SAMPLE_ORDER_NUMBER		Yes		No	Yes	
	SAMPLE_STRUCTURE		No		No		
	SAMPLE_TITLE		No		No		
	STATUS	30	Yes	30	No	Yes	
	STATUS_COMMENT		No		No		
	URGENT SAMPLE ORDER		No		No		
1						▶	11.

#### The SAMPLE\_ORDER Table

#### 17.1.3

Field	Description
Sample_Order_Number_Key	The unique, internal number used as the primary key for the SampleTrack™ SAMPLE_ORDERS table.
Sample_ID	The unique sample ID used for the sample in SampleTrack™.
Sample_Title	Free entry (could be used for the plot title in older versions).
Sample_Nature	Free entry.
Sample_Charge	Free entry.
Sample_Fraction	Free entry.
Project_Number	Free entry.
Date_Entry	Date when the record was created.
Date_Start	Date when the sample preparation was started.
Date_End	Date when the last experiment was finished.
Status	The status of the sample. The following status are available:
	10 Created
	20 Verified
	30 Started
	40 Working
	50 Pause
	60 Finished
	70 Approved

80	Archived	
----	----------	--

95 Instrument error

The SampleTrack<sup>™</sup> Client ID abbreviation (a required field).

#### The METHODS Table

17.1.4

Field	Description
Sample_Order_Number_Key	The foreign key used for the sample orders table.
Method_Type_Key	The method type (currently only NMR).
Plot_Title	The title for the spectra plot.
Journal_Number	The journal number field.
Sample_Changer_Holder	The last position of a tube in the sample changer holder.

#### The NMR\_PREPARATION\_WS Table

User\_Name\_Key

17.1.5

Field	Description
NMR_Preparation_WS_Key	The unique, internal number used as the primary key for the SampleTrack™ NMR_PREPARATION_WS table.
Sample_Order_Number_Key	The foreign key used for the methods table, used in conjunction with the Method_Type_Key.
Method_Type_Key	The foreign key used for the methods table, used in conjunction with the Sample_Order_Number_Key.
Sample_Amount	The quantity of the sample being used.
Solvent_Amount	The quantity of the solvent being used.
Solvent_Name_Key	The solvent name used in the SampleTrack™ preparation (a required field).
Instrument_Name_Key	The instrument name used in the SampleTrack™ preparation (a required field).
Preparation_Name_Key	The preparation name used in the Sample- Track™ preparation (a required field).
User_Name_Key	The user who has prepared the sample.
Date_End	The date when the preparation was finished.
Status	The status of the preparation. Refer to the avail- able status in the SAMPLE_ORDERS table

The Experiment Table

Field	Description
NMR_Experiment_WS_Key	The unique, internal number used as the primary key for the SampleTrack™ NMR_EXPERIMENT_WS table.
NMR_Preparation_WS_Key	The foreign key used for the NMR_PREPARATION table.
Experiment_Name_Key	The unique name of the SampleTrack <sup>™</sup> experi- ment (also composite).
Composit_Experiment	The name of the sub-experiment.
Experiment_Type_Name_Key	The type of experiment (1H, C12).
Experiment_ID	The ID used for each experiment.
Instrument_Name_Key	The instrument name used in the SampleTrack™ experiment (a required field).
Result_Data_Path	The local path for the instrument data results.
User_Name_Key	The user who has performed the experiment.
Date_Start	The date the experiment was started.
Date_End	The date when the experiment was finished.
Status	The status of the experiment Refer to the avail- able status in the SAMPLE_ORDERS table

#### VCU Request Tables

Introduction to VCU REQUEST Tables	17.2.1
	17.2.1

The VCU\_REQUEST\_ALL view is utilized when you desire to compile statistical or report data to view with an external program such as Crystal Reports<sup>™</sup> or Microsoft® Excel®. For example, you can compile a report in Microsoft® Excel® that shows all of the Proton experiments that have been completed.

A basic understanding of ODBC, SQL and tables is assumed when working with external programs. Refer to your external program handbook for assistance.

The fields that are listed in the following tables will assist you in configuring a link from your external program (using e.g. ODBC or SQL) to various data in Sample-Track<sup>™</sup>.

#### The VCU Sample Order Table

Field	Description
User_Name_Key	The client short ID, must exist in SampleTrack™.
Last_Name	Last name of the client.

17.2

17.2.2

#### **VCU Request Tables**

First_Name	First name of the client.
Department_Name_Key	Department name, must exist in SampleTrack™.
Sample_ID	The unique sample ID used for the sample in Sample- Track™.
Sample_Title	Free entry (could be used for the plot title in older versions).
Sample_Nature	Free entry.
Sample_Charge	Free entry.
Sample_Fraction	Free entry.
Project_Number	Free entry.
Date_Entry	Date when the record was created.
Date_Start	Date when the sample preparation was started.
Date_End	Date when the last experiment was finished.
Status	The status of the sample. The following status are available:
	10 Created
	20 Verified
	30 Started
	40 Working
	50 Pause
	60 Finished
	70 Approved
	80 Archived
	95 Instrument error

#### The VCU Preparation Table

Field	Description
Method_Type_Key	Method Type (only NMR in the moment).
Plot_Title	Title for the plot of the spectra.
Sample_Amount	Amount of the sample used.
Solvent_Name_Key	Solvent name for the preparation, must exist in Sample- Track™.
Instrument_Name_Key	The instrument name used for the preparation, must ex- ist in SampleTrack™.
Preparation_Name_Key	The name of the preparation, must exist in Sample- Track™.
User_Name_Key_1	User who started the preparation.
Date_Start_1	Date when the preparation was started.
Date_End_1	Date when the preparation was finished.

### SampleTrack<sup>™</sup> Tables

Status\_1

Status of the preparation. The status could be the same as one of the above.

#### The Experiment Table (VCU\_REQUEST\_ALL)

17.2.4

Field	Description
FIEIU	Description
Composit_Experiment	The name of the sub-experiment.
Experiment_Name_Key	The unique name of the SampleTrack™ experiment (al- so composite).
Experiment_Type_Name	e_KeyThe type of experiment (1H, C12).
Experiment_ID	The ID used for each experiment.
Instrument_Name_Key_	1The instrument name used in the SampleTrack <sup>™</sup> experiment (a required field).
Result_Data_Path	The local path for the instrument data results.
User_Name_Key_2	The user who has performed the experiment.
Date_Start_2	The date the experiment was started.
Date_End_2	The date when the preparation was finished.
Status_2	The status of the experiment. The status could be the same as one of the above.


## EasyDialog Configuration

# 18

#### **Opening the EasyDialog Configuration Window**


18.1

You will need to have administrator privileges to perform the EasyDialog configuration. To open the EasyDialog designer open the Administration pull-down menu and select: *Start EasyDialog Designer* 

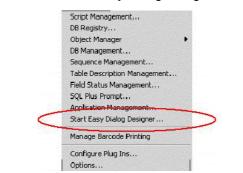





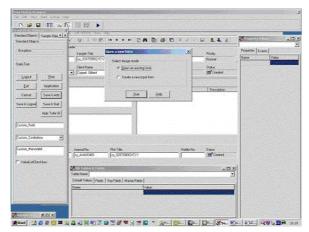

#### Opening a Form / Creating a New Form

18.2

The configuration of EasyDialog is accomplished through a object-oriented, drag, drop & modify methodology. To prepare a user input or request form you must only select an object from a list, drag it onto a design screen, and edit the specific

parameters for that object. By selecting various pre-programmed object modules in this manner you can adjust the form to fit the individual needs of your user or group of users.

There are currently two modes in the EasyDialog configuration window: Open an Existing Form or Create a New Input Form.

Upon entering the configuration window you will be presented with the following window where you can select the design mode that you wish to use:

Figure 18.3. Open / Create a New Form

en a new form	
Select design mode	
Open an existing form	
C Create a new input form	
Start	

If you wish to open an existing form to revise or to use as a template select Open an Existing Form. You will then be able to modify the form, and to save the changes that you make to the form as the original form name, or under a new name.

If you wish to create an entirely new input form, select Create a New Input Form.



Note: An input form is a form (or input screen) where a user supplies information, for example, for a preparation or an experiment that they want to perform.

#### The Configuration Window

18.3

There are five main panels which make up the EasyDialog Configuration Tool:

- The Administrator Toolbar
- The Drag & Drop Objects
- The Property Editor
- The Fields Overview
- The new form itself

#### Drag & Drop Objects


18.4

The Drag & Drop Objects window consists of a collection of pre-defined objects. These objects can then be used to build new input forms.

Figure 18.4. Drag & Drop Objects

andard Obje Sa	nple Objec Metho	d Object Prep Ob	ects Exp Objects
Standard Objects			
Groupbox			
StaticText			
Print	Print HTML		
Logout	Application		
Exit	Cancel		
	Save & Buffer		
Save & Add	Save & Quit		
Save & Logou	Appr Tube ID		
Custom_Field			
Custom_Combo	bax 👻		
Custom_Memofi	eld		
C Custom_Radi			
Custom_Cheo	kbox		
ValueListChe	ckbax		
C Printer_Selec	tion		
Active_Printer	Charles and Charles and		

The objects are seperated systematically within Tab windows, grouping standard objects like 'printer selection', 'Save' buttons or custom fields, custom combo boxes, check boxes or radio buttons, that are not defined in the standard database fields.

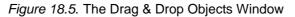

Standard Obje Sample Object Method Object Prep Objects	Exp Objects
Sample Order	- 1
Sample_ID	
Contraction of the second s	
Sample_Title	
[Sample_Time	
Client_ID •	
Client_Name	
Tonen Trane	
ClientDepartment	
ClientCompany	
Urgent Sample	
Tip_0t_Day	
Land the second s	
ABCDEFGH	
12000000012	
"00000000"	
******	
**************************************	
**************************************	
*0000000e	
NOOOOOOO0	
**************************************	
**************************************	
NOCOCOCOC	

The fields in the sample order, method ,preparation and experiment tabs are typical fields used in an adapted EasyDialog form. Custom fields are saved in the database in the same manner as the standard fields.

#### How to Drag & Drop Objects

This object-oriented programming approach to form building, allows a non-programmer to quickly and easily build new EasyDialog forms. To use these objects in a new form you must only select one, click once on the object with your left mouse button and hold the button down while you drag the object onto your new form. The figure below illustrates this simple drag and drop process.

Figure 18.6. Drag - Drop Objects

	Aannii	nistrator i colbar - New_Lotti (Alij	
l	<u>F</u> ile	⊻iew <u>O</u> ptions <u>H</u> elp	
	Ľ	🛩 🖬 🔳 🖬 🔛	
Ī	_ [	Drag&Drop Objects 🛛 🛛 🛛	
I	11	Standard Objects Input Objects	SampleTRACK Easy Dialog
I	17	Sample Order	
I	14	Sample_ID	
I	LI	Sample_Title	Sample ID
I	18		99031504
I			
I	19	Client_Name Click	
I		ClientDepartm Select	Drog 4
I			Drag Drop
I	20	ClientCompany	ыор
I		General	
I	21	Groupbox	
I			
1		StaticText	

#### Grids

18.6

In order to simplify the alignment of the objects in the new form, a grid system has been implemented. By enabling the grid (Options - Grid - Enable Grid) you can align the new objects on the X and Y coordinates of the grid lines.

The distance between the grid lines is adjustable. To change the grid distance select Options - Grid - Change X or Change Y Grid Distance. An input window will appear which will allow you to enter the new grid distance:

Figure 18.7. Changing the Grid Distance

Grid Distance	X
X Distance:	
20	
OK	Cancel

#### The Property Editor

18.7

Once an object has been dropped into a new form you can configure the object field or title using the Property Editor. To configure a new object, click on the object in the new form with your mouse. The Property Editor will display the current attributes of the object. You can then make any changes to these attributes as needed. When you have several objects on your form you must only click on the corresponding object to change its attributes.

SampleTRACK Easy Dialog	💶 🗖 🗙 🛛 Field Ed	Field Editor		
	Sample	Sample_ID : TDBEdit		
	Caption	n Sampl	e_ID	
Sample ID	Datase	t qSamp	pleOrder	
•	Fieldna	ime SAMP	'LE_ID	
	Enable	ed True		
	Height	21		
	Left	99		
	Name	Sampl	e_ID	
	Тор	83		
	TabOr	der 0		
	Tag	0		
	Visible	True		
	Width	169		
	Requir	ed False		
	SaveT	oBuffer False		

Figure 18.8. Configuring a New Object for EasyDialog

#### The Fields Overview Window

18.8

The Fields Overview window is a second source of objects for your new form. The fields listed here are the actual fields that are found in the SampleTrack<sup>™</sup> database. Normally these should only be used if an object is not found in the Drag & Drop Objects, or if a special definition is required.

Figure 18.9.	The Field	Overview	Window
--------------	-----------	----------	--------

Fields Overview							×
Field Overview	Fields						
	Fieldname	Value	Required by DB	Default Value	Used	Set From DB	
■	SAMPLE_FRACTION		No		No		
■ NMR PREPARATION WS	SAMPLE_ID	99031603	Yes		Yes		
• NMR_EXPERIMENT_WS	SAMPLE_NATURE		No		No		
	SAMPLE_ORDER_NUMBER		Yes		No	Yes	
	SAMPLE_STRUCTURE		No		No		
	SAMPLE_TITLE		No		No		
	STATUS	30	Yes	30	No	Yes	
	STATUS_COMMENT		No		No		
	URGENT SAMPLE ORDER		No		No		_ <b>_</b> _
	•						• //

The Fields Overview window in the example is broken down into four main sections: SAMPLE\_ORDERS, METHODS, NMR\_PREPARATION\_WS, and NMR\_EXPERIMENTS\_WS. Each section has several corresponding fields, as shown in the Fields box in the figure above, which can be dragged & dropped onto a form.

The SAMPLE\_ORDERS fields represent the main information that will be included in EasyDialog. The METHODS section contains the fields that pertain specifically to methods. Likewise, NMR\_PREPARATION\_WS and NMR\_EXPERIMENTS\_WS contain fields that pertain specifically to worksteps for preparations and experiments, respectively.

#### Fields Overview Required Values

Certain fields are required by the database, and are identified with a **Yes** in the **Required by DB** column. When a field is required by the database, then a value must be provided. This value can be derived from three sources. The first is that the value is set by the database automatically. In this case the **Set From DB** column will contain a **Yes** value. The second method is to define a default value for the field (see <u>"Default Values in the Fields Overview Window" on page 128</u>). If the value is not derived from the first two methods, then the field must be present in the EasyDialog form so that the user can enter this value.

#### Default Values in the Fields Overview Window

18.10

When you double-click with your mouse on one of the fields, a window will appear that will allow you to enter a default value to the corresponding field:

Figure 18.10.Default Field Value in the Fields Overview Window

SAMPLE_ID	×
Default Value	
Test001	
OK Cancel	

The default value is then displayed in the **Default Value** column of the Fields Overview window.

Figure 18.11.Default Value in the Fields Overview Window

Fields				
Fieldname	Value	Required by DB	Default Value	Used
LIMS_IMPORT		No		No
MODIFIED_BY		No		No
PROJECT_NUMBER		No		No
SAMPLE_CHARGE		No		No
SAMPLE_FRACTION		No		No
SAMPLE_ID		Yes	Test001	No
SAMPLE_NATURE		No		No
SAMPLE_ORDER_NUMBER		Yes		No
SAMPLE STRUCTURE		No		No
•				

This default value will then be displayed when you start EasyDialog.

🜃 SampleTRACK Easy Dialog	
Sample ID	
Test001	

#### The STATUS Field

There are several pre-defined status entries used by SampleTrack<sup>™</sup>. The table below lists the current status codes and their meaning:

Status Code	<b>Description</b>
10	Created
20	Verified
25	Waiting
30	Started
60	Finished
70	Approved

In the Fields Overview window, the SAMPLE\_ORDERS, the NMR\_PREPARATION\_WS, and the NMR\_EXPERIMENTS\_WS categories each contain a STATUS field. By using this STATUS field you can assign a default status code value to each category from the table above.

The point at which a Sample Order, Preparation or Experiment is assigned one of these status's is completely definable by the administrator in the EasyDialog configuration. For example, in one laboratory setup, the user might bring the sample to the laboratory where further preparations will be carried out, and the responsible laboratory personnel will place the sample on the sample changer. In this system the user's EasyDialog should set the Sample Order to status "Created" or "Verified" when the user presses OK. The laboratory personnel would then use the Verify Sample Orders or Manual Preparations windows in the Client Tool to upgrade the status's when the preparation has been completed.

In an open access system, for example, where the user has access to a spectrometer and can place the samples that they have prepared on the sample changer themselves, EasyDialog must be configured so that the user can manually enter a Tube ID and Sample ID. In this case, when the user presses the OK button in EasyDialog, the Sample Order status should be set to "Started", the Preparation status to "Finished", and the Experiment status to "Verified".

There are of course many different versions of these scenarios that could be used in a laboratory. EasyDialog can be configured for your individual laboratory situation.

In any case the method used to configure which status will be achieved at what point, is completely programmable by the administrator in the EasyDialog configuration.

#### Configuring a Status

18.12

To configure the status of the Sample Order, Preparation or Experiment, doubleclick with your mouse on the "Status" field located in the Fields Overview panel under SAMPLE\_ORDER, NMR\_PREPARATION\_WS, or NMR\_EXPERIMENT\_WS, respectively.

A default value window will appear (see the figure below) where you can place one of the status codes. For example, to set the Sample Order to **10** where it will be placed on status **Created**, when the user clicks the OK button in EasyDialog. Double click with your mouse on the **Status** field under SAMPLE\_ORDER in the

**BRUKER BIOSPIN** 

Fields Overview window. When the default value window appears enter the status code for **Created**, which is **10** in the input box, as shown in the figure below.

Figure 18.13. Entering a Default Status Value

STATUS			X
Default Value			
10			
	ОК	Abbrechen	]


Press the **OK** button in the default value window to accept this new value. You should now see your new default value listed under the default value for SAMPLE\_ORDER - Status in the Fields Overview window.


You can likewise set a default value listed in the Sample Status Code table to the preparation status and experiment status in this manner.

#### The Administrator Toolbar

#### 18.13

The Administrator Toolbar consists of a pull-down menu and several buttons which allow you to select various routine (e.g. Open, Save) and specialized tasks or options (e.g. environment variables, grid settings). You can also choose which panels are displayed, by (de)selecting the corresponding button from the button bar.

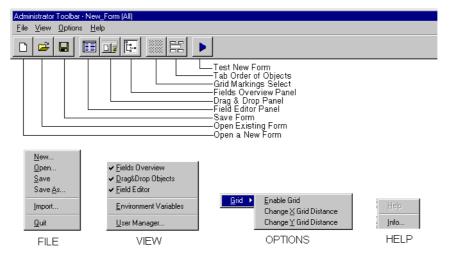

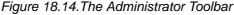

#### Example: Creating a New Input Form

18.14

The following form example provides a detailed step-by-step description of how to build a new input form. If you have not already done so, read the help section <u>"Opening a Form / Creating a New Form" on page 123</u> to select a design mode before proceeding with this example.

#### Creating a New Form: Opening the Form

To create a new form select Create a New Input Form from the opening prompt, or File - New from the EasyDialog menu, then Create a New Input Form.

A new form will appear, similar to the one in the following figure:


Figure 18.15.A New Input Form

Administration of the streag Part (AI)												×
For Sines Types media												
🗅 🛎 🖬 🖬 🗊 🕼	- 13	831	•									
Darffing Lade	1	Sample I RA	CE Escue Du	aleu							Felc Editor	R.
Stander: Insetuncer											Herea Talo	-
Schold Jilder	100									11		
Sunger L'	le -											
Sunge 106	I											
	le -	::::	:::	:::				:::		11		
										- 11		
Qua Sanz 🔹												
ClexDecalment												
										11		
ClientCompany												
Lierere	il											
a finage as	1.1											
	1 · · ·											
csitate cet												
· · · · · · · ·										11		
_opol Diri										11		
	1.1											
:		::::	:::	:::								
								111		111		
L	1							• • •				
		::::		:::				111		111		
	pa .									- 11		
											1	
										11		
	1235	<pre>kLevey</pre>										a.
		Overdes.		-	ılıc	_		_		_		
		SOMELS: OF	1.FBS	- I - 2	c drame		Volue		Reputed by .	њ ре	la.#Value Lood	Section DB -
		VETHERS			.CIC:0001	LE9			Чc		\c	
		We			CHIVE RULE				1c		<u>\c</u>	
	÷.	willout	) MERTAR	- IS	HEN DAL HEN NGCIER THE	-MILAH			16 36		5	
				12	ти млангала и	>			i.		No.	
					ik tet di Ne		DCHO_C	X0HPM	°s:		\c	
					SHC _1 HI				Чc		NC	. E

#### Creating a New Form: Establishing Group Boxes

#### 18.14.2

In creating the outline of a new form it is normally a good design methodology to first establish certain input areas which group related fields. This is accomplished most easily by using a "Group Box". In our example here I will use Sample Order, Preparation and Experiments. Drag three group boxes from the Drag & Drop Objects window onto you new form. Fill-in the title for each group box by clicking once on the box to select it, and by filling in the Caption block of the Field Editor. When you have completed these steps your form should look something like the following figure.

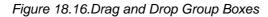

12	SampleTRACK Easy Dialog	_ 🗆 ×
	r Sample Order	
:		
:		
:		
:		
÷	Preparation	
:		
2		
:	L	
:	Experiments	

#### Creating a New Form: Adding the First Object

18.14.3

Drag & drop a Sample\_ID object from the Drag & Drop Objects onto the Sample Order group box.

Figure 18.17.Drag and Drop Field Sample ID

SampleTRACK Easy Dialog		
Sample Order		
. 99031710		
· · · · · · · · · · · · · · · · · · ·	· · · · · · · · · · · · ·	<u>.</u>

Don't be alarmed if the object doesn't fit properly, you can adjust the size of the Sample Order group box by clicking once on the group box to highlight it, then by pulling on one of the highlighted corners to adjust the size; or by changing the respective values (Left, Top, Width, and Height) in the Fields Object window.

#### Creating a New Form: Adjusting the Group Box

18.14.4

Figure	18.18.Adjusting	the Size	of an	Object

📽 SampleTRACK Easy Dialog	 	
Sample Order		
Sample ID		•
99031710		:
		÷
		:
Preparation		· · · · · · ·

You can also adjust the size of an object in the same manner as described above.

To continue with our example, add the following objects to your Sample Order group box: Client\_ID, Client\_Name, Client\_Department and Client\_Company. Adjust the size of the objects and group box until they similar to those shown in the figure below. If you desire a Client Department and Client Company title, drag & drop two StaticText objects to the Sample Order group box and fill in the appropriate Caption in the Field Editor.

#### Creating a New Form: The Completed Group Box

18.14.5

Figure 18.19. The complete Group Box

🜃 SampleTRACK Eas	sy Dialog		_ 🗆 ×
Sample Order			
. Sample ID			
99031712			
Client ID	Client Name	Client Department	Client Company
	Administrator, User 💌	DEMO_DEPARTMEN	DEMO_COMPANY

You will notice that the values (ADMIN, Administrator, User, DEMO\_DEPARTMENT and DEMO\_COMPANY) that are shown in the figure above are not in the field objects that you have entered. These will appear the first time you run this new form. The values represent the internal field values from the SampleTrack<sup>™</sup> database.

Creating a	New Form:	Adding the	e Other	Group	Boxes
------------	-----------	------------	---------	-------	-------

18.14.6

Now fill in the fields as described previously until the Preparation and Experiment group boxes are completed as shown in the following figure:

🜃 SampleTRACK Easy Dialog						_ 🗆 ×
. Sample Order						
. Sample ID						
99040162						· · · · · · · · · · · · · · · · · · ·
Client ID Client Name		Client Depart	ment	Client Co	mpany	
ADMIN Administrator,	User 💌	DEMO_DEP	ARTMEN	DEMO_0	OMPANY	
C Chloroform		Tube ID				
C DMSO						· · · · ·
Experiments						
1.Experiment		•		:::::		
		-				
2. Experiment mit D20						
. 3. Experiment						
:						
				::::		

Figure 18.20.Adding Group Boxes

#### Creating a New Form: Adding Buttons

#### 18.14.7

The final configuration features that needs to be added to complete our example are the task buttons. The Static Text buttons are pre-defined and configurable, you need only to drag and drop the objects onto your new EasyDialog form and adjust the parameters in the Field Editor. The exact buttons that you use depend on the tasks that you want the user to perform. In the following example I have added some of the normal buttons that are used in EasyDialog.

Sample Order Sample ID	
99040162	
·	
Client ID Client Name	Client Department Client Company
ADMIN Administrator, User	DEMO_DEPARTMEN     DEMO_COMPANY
Preparation	
Chloroform	Tube ID
C DMSO	
Experiments	
1.Experiment	:: <u>Print</u> : <u>Logou</u>
<b>_</b>	<u>Exit</u>
2. Experiment mit D20	Save & Add
·	
	Save & Logout
3. Experiment	
▼	

Figure 18.21.Adding Buttons in a Form

Saving an EasyDialog Configuration

18.14.8

To complete the EasyDialog form select File - Save As from the EasyDialog main menu. A "Save As" window will appear, as shown in the following figure, that will prompt you for a Configuration Name and User/Group under which to save the ile.

Figure	18.22.Save	EasvD	Dialog	confiaura	ation

Configuration	User/	Group
EZ_STANDARD	All	
🕼 EZ STANDARD NIL	All	
Config Name EZ_STANDARD_0M_Test	_	Save

Enter a configuration name such as EZ\_STANDARD\_<new name>, where <new name> is a name that you give the file. If you want to save the configuration under an existing name (e.g. EZ\_STANDARD), select the name from the configuration list - this will overwrite the existing file.

For the User/Group select a group or user from the pull-down menu. If you want to make the new EasyDialog configuration available for all users, select "All", otherwise select the specific user or group that you want the configuration assigned to.

This completes the steps for creating a new form.

/Username:<usr>: The login user name: e.g. GUEST

/Password:<Passwrd>: Password for a corresponding user.

/Admin: The program will be opened in EasyDialog configuration mode (the Toolbars appear).

/Config: EasyDialog is opened with the Load Configuration window to allow Easy-Dialog configuration selection.

/Config:<Name>: EasyDialog is opened with the <name> as a parameter. When this name does not exist the program will be stopped.

/SysButtons: Displays the system Windows button and system menu when Easy-Dialog is opened.

### EasyDialog Configuration

## E-Mail Support

# 19

#### E-Mail Support

An additional feature of SampleTrack is the ability to notify users of finished samples or experiments per E-mail. In order to notify users per E-mail, the E-mail addresses must first be added to the E-mail listing in the Environment Setup – Users – Basic Information window. New users can either be inserted manually or through file import.

The generation of E-mails is executed by the database and the SampleTrack Communication Server. The Communication Server handles E-mail support like an instrument, no user interaction is necessary. To activate E-mail notification, check **Served By Communication Server** in the **Envoronment Setup / Instruments / E-mail** window.


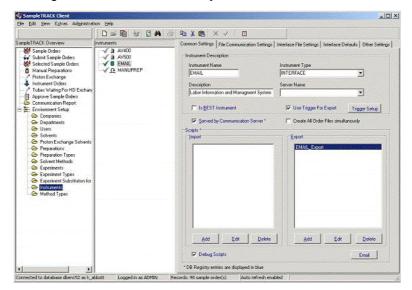


Figure 19.1. E-mail Served by the Communication Server

#### Installation of E-mail support

Start the installation in the Administration menu, DB Management / Load Components.

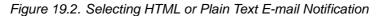

Summer and the second s	Components		
	Name	Current	New 🔺
	BEST Spectrometer Samp Print Sample Order Details HTML E mail Notification S Plain Text E mail Notificatio Web Configuration for San	Sample Configuration ample Configuration on Sample Configuration apleTRACK	

#### Installation steps

- Definition of the Interface in the Instrument Setup for E-mail support.
- Definition of the trigger which exports data to the E-mail program.
- Definition of parameters for E-mail support.

#### System Requirements

SampleTrack V2.35 ScriptLib V3.3 DB V2.039

#### E-mail Internal Processes

The following internal processes are performed when using the E-mail notification in SampleTrack. First a database trigger generates a dataset in the table INTERF\_TRIG\_TAB, which provides a list of samples, for which an E-mail notification needs to be performed. This trigger is activated in the database each time a sample or an experiment is set to the **Finished** status. The SampleTrack Communication Server scans the INTERF\_TRIG\_TAB table for available datasets. If a dataset is found, the Communication Server runs the predefined script which executes the E-mail notification. The script contains commands, that are integrated with the Communication Server, which prepare data for E-mail distribution. After finishing this process the current dataset is deleted.

19.2.2

Automatic E-mail notification is supported by the SampleTrack Communication Server. As the Communication Server only supports instruments, an instrument interface must be defined before automatic E-mail notification can be used. The following steps must be performed in order to define an instrument interface:

Select Environment Setup - Instruments from the tree, and press the New button to define a new instrument.

Assign the name EMAIL to the new instrument (all instruments should have capital letters).

Select the instrument type INTERFACE.

Select the script EMAIL\_Export from ScriptLib (ScriptLib V3.3 is required).

Press the Save button to save all the data to database.

#### E-mail Trigger Setup

19.2.4

The button **Trigger Setup**, in the Common Settings window, is used to define the event which determines when the E-mail should be sent. Possible events include: experiments, preparations, methods and sample orders. E-mail notification is activated following the completion of one of these events. The event option can be selected from the pull-down box **Fire Trigger When**. After selecting the event option, you must also select one of the checkbox options under **Has Following Status**, which indicates the status of when the trigger should be fired. Standard options for these two parameters are **Status Finished** for the status, and **Sample Order** for the table. On the right side of the Trigger Setup window is the internal database trigger, displayed as SQL text.

To activate the trigger, the checkbox **Use Trigger For Export** must be checked in the Common Settings window.

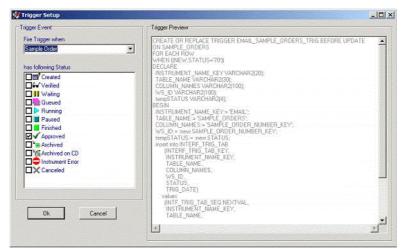


Figure 19.3. The E-mail Trigger Setup Window

#### E-mail Parameters

To edit the E-mail parameters press the button **E-mail**, which will open an E-mail parameter dialog window.

Figure 19.4. E-mail Parameters

💞 Email Setup			\$2.45 A		
SMTP Server					
Host	F	ort	Charset		
mailserver.bruker.de		25	us-ascii		
Mail					
From Address	To Address	MaHD		Ok.	
samtrack@bruker.de	[USERS.E_MAIL]	[SAMPLE_ORDERS	SAMPL	Cancel	
Subject		Mail Ts	0e		
Report for ISAMPLE_01	RDERS.SAMPLE_ID]	Plain	•	Import	
				Export	
Body Attachments				C. C	
NMR_EXPER<br /NMR_EXPER<br Your SampleTF	RIMENT_US>	MR_EXPERIMENT_WS	EXPERIME	NT_NAME_KEY]	(INNR_EXPERIMENT_VS.EXPERIM
SampleTRACK S	iystem, [ProductV	ersion]			

E-mails can be configured as HTML or plain text format. The following information must be modified:

Field	Value	Description
Host	<your host="" server="" smtp=""></your>	Host where the E-mails will be sent.
Port	25	Port for SMTP server.
Charset	US-ASCII	Different locales have differ- ent character sets. Default is US-ASCII.
User	Samtrack	Name of E-mail sender
Mail ID	SAMPLE_ORDERS.SAMPLE_ID]	Unique identifier, to identify E-mail.
Subject	Sample '[SAMPLE_ORDERS.SAMPLE_ID]' finished	Title for the E-mail.
Attachments	[SAMPLE_ORDERS.DOCU_FILES]	Attachments that should be sent with the E-mail.
	The following fields are contained in the mail con body entries (dependant on the number of experiments)	-
	Haadaa	

Header

The following experiments in the sample order [SAMPLE\_ORDERS.SAMPLE\_ID] are finished:

#### Body

[NMR\_EXPERIMENT\_WS.EXPERIMENT\_NAME\_KEY] on [NMR\_EXPERIMENT\_WS.INSTRUMENT\_NAME\_KEY]

Text for bodies. Mail can consist of several entries.

#### Footer

Please remember to pick up you plots.

Thank you for choosing SampleTrack.

Text for footer information.



Note: Once you have finished the configuration, you must restart the Communication Server in order for the changes to take effect.

#### E-mail Interface Trigger Table

19.2.6

#### Description for INTERF\_TRIG\_TAB

The following fields have been predefined for the INTERF\_TRIG\_TAB table:

Field Name	Туре	Description
INTERF_TRIG_PK	Number (22)	Unique internal number.
INSTRUMENT_NAME_KE	YVarchar (20)	Interface (instrument) name; dependant on the trigger instrument (will be set for E-mail support instrument: E-mail).
TABLE_NAME	Varchar (30)	The name of the table that has fired the trigger.
COLUMN_NAMES	Varchar (100)	Reference fields that identify the dataset (SAMPLE_ORDER_NUMBER_KEY, METHOD_TYPE_KEY, NMR_PREPARATION_WS_KEY, NMR_EXPERIMENT_WS_KEY).
WS_ID	Varchar (100)	Reference values that identify the dataset (1234,NMR, 1235,1236).
STATUS	Varchar (4)	The status that has fired the trigger.
TRIG_DATE	Date	The date and time when the trigger was fired.

#### E-mail Address Registering

The E-mail address is registered in the environment user setup

SampleTRACK Client		
Eile Edit View Extras Administration Help		
] D 🛩 🛍 😽 🖬 🖂	新 🖻 🕹 🛍 🛛 🗙 🧏 🖗	
SampleTRACK Overview SampleTRACK Overview Matrix Order BEST Order Pooling Order Pooling Order Approve Ma CSI CC	Basic Information         User Short ID         GDH         Last Name         Hiss         Description         gdh         Email         gdh@bruker.de         Password         Date Password Changed         [19.01.2001         User Status         © Invalid user         Phone/Fax	
Connected to database dben152 as demo_labor Logged in as AD	MIN Records: 25 sample order(	s) Auto refresh enabled //

Figure 19.5. E-mail Address in the Environment User Setup

## Data Exchange & Directory Structure

# 20

#### Introduction to Data Exchange

SampleTrack<sup>™</sup> contains several predefined table structures that are used for the organization and display of key data both internally and in external programs. This chapter provides a description of the individual fields used in these table structures.

There are several fields found in the four base tables used internally by Sample-Track<sup>™</sup>. Refer to Introduction to SampleTrack Base Tables for details.

There are also several fields that are used by the view VCU\_REQUEST\_ALL. This view is used when viewing data from an external program such as Crystal Reports<sup>™</sup> or Microsoft® Excel®. Refer to the Description of the VCU\_REQUEST\_ALL Tables for details.

#### SampleTrack™ bstsdata Directory Structure

When using the SampleTrack<sup>™</sup> software it is essential that the following directory structure is created as described below:

All instrument directories must be located in a directory labeled bstsdata.

For each instrument there must be a separate subdirectory with the specific instrument name. These instrument subdirectories must consist of the necessary in and out directories which are used for file communication.

The **In directory** is where instrument input files are to be located.

Instrument	File Type
LIMS	SampleTrack <sup>™</sup> report files for the company LIMS
PREP	Preparation order files from SampleTrack™
SPECT	Spectrometer order files from SampleTrack™

The **Out directory** is where the instrument reports are to be located.

Instrument	File Туре
LIMS	LIMS work lists for SampleTrack <sup>™</sup> import
PREP	Preparation report files from SampleTrack™
SPECT	Spectrometer report files from SampleTrack™

**BRUKER BIOSPIN** 


20.2

20.1

Each In or Out directory must contain a save and an error sub-directory.

When the files have been accessed successfully by the SampleTrack<sup>™</sup> system, or any other instrument, they are moved to the associated save directory.

When the access was unsuccessful (e.g. a syntax error) the files are moved to the error directory.

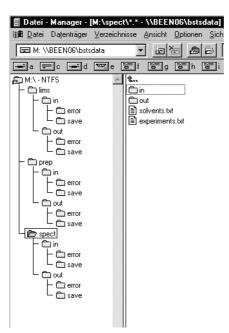

# JCAMP Definition File

# 21

21.1

# JCAMP Definition File Introduction

The communication files serve as the exchange between SampleTrack<sup>™</sup> and the served instruments and systems like spectrometers, preparation robots or LIMS.

Each instrument has a separate In and Out directory defined for its information exchange.

#### In Directory:

Order-File created by SampleTrack

Information:SampleTrack --> Instrument

#### **Out Directory:**

Report-File created by Instrument

The **Report File** contains the status or result of a measurement or preparation order.

Several files can be generated for one order.

#### Information:Instrument --> SampleTrack

The control software for each instrument scans the In directory for current orders, reads the incoming files and reports the results into the out directory. The Out file structure is the same as the incoming file structure completed with the result and order status information.

#### Workflow Example:

SampleTrack generates an order file for the preparation robot for each preparation .

File name: Sample-ID.[\*]p(valid file nomenclature: abc123.p - abc123.10p - acb123.11p - abc123.102p)

After the preparation the preparation robot writes the report file:

File name:Sample-ID.[\*]r

SampleTrack generates for each experiment / measurement an order file for the spectrometer.

File name: Tube-ID.10r; Tube-ID.11r

The spectrometer generates for each experiment / measurement 2 report files.

**BRUKER BIOSPIN** 

File name: Tube-ID.11s (**Status**-Info 1.Exp); Tube-ID.12s (**Result**-Info 1. Exp.) Tube-ID.13s (**Status**-Info 2.Exp); Tube-ID.14s (**Result**-Info 2. Exp.)

A communication file includes two parts: Header and variable part.

The header rows are preceded by two hash marks (##). The header information is defined by the universal JCAMP format.

The variable part is preceded by the hash and dollar symbols (##\$). Variables preceded by ##\$ are user defined and in alphabetcal order.

#### Header:

##TITLE= Setup File, SampleTrack Software Version x.x

(x.x The current version number)

##JCAMPDX= 5.0

(JACMP Version)

##DATATYPE= Parameter Values

(Data or parameter file)

##ORIGIN= BRUKER BIOSPIN GMBH

(Responsible company name)

##OWNER= Client company, that uses the data

#### Variable Part:

##\$ACAMREQUEST= NOT USED

(Table: NMR\_EXPERIMENT\_WS . AC\_AM\_REQUEST)

Special requests for AC or AM instruments.

'YES' or empty

Format: string (max. 40 digits).

--> Generated by SampleTrack.

--> Interpreted by spectrometer.

#### ##\$ACCOUNT= NOT USED

(Table: DEPARTMENTS.COST\_CENTER)

The client cost center

Format: string (max. 40 digits).

--> Generated by SampleTrack.

--> Interpreted by spectrometer --> JCAMP data file.

##\$ANALYST=

(Table: NMR\_EXPERIMENT\_WS . USER\_NAME\_KEY)

Spectrometer operators ID (Not necessarily the logged in user).

**BRUKER BIOSPIN** 

Format: string (10 digits).

--> Generated by SampleTrack.

--> Interpreted by spectrometer --> JCAMP data file.

##\$BESTMETHOD=

(Table: SAMPLE\_ORDERS . USER\_NAME\_KEY)

Client user ID

Format: string (10 digits).

--> Generated by SampleTrack.

--> Interpreted by spectrometer --> JCAMP data file.

##\$CLIENTCODE=

(Table: SAMPLE\_ORDERS . USER\_NAME\_KEY)

Client ID

Format: string (10 digits).

--> Generated by SampleTrack.

--> Interpreted by spectrometer --> JCAMP data file.

##\$CLIENTDEPARTMENT=

(Table: USERS . DEPARTMENT\_NAME\_KEY)

Client department ID

Format: string (max. 40 digits).

--> Generated by SampleTrack.

--> Interpreted by spectrometer --> JCAMP data file.

##\$CLIENTEXPERIMENT= NOT USED

(Table: NMR\_EXPERIMENT\_WS.CLIENT\_EXPERIMENT)

Format: string (max. 40 digits).

--> Realized with parameter ##\$EXPERIMENT.

##\$CLIENTNAME=

(Table: COMPANIES . COMPANY\_NAME)

Company name

Format: string (max. 40 digits).

--> Generated by SampleTrack.

--> Interpreted by spectrometer--> JCAMP data file.

##\$CLIENTSOLVENT= NOT USED

(Table: NMR\_PREPARATION\_WS.CLIENT\_SOLVENT)

Client solvent: H2O,THF,DMSO,D2O,CDCL3,...

Format: string (max. 40 digits).

--> Realized with parameter ##\$SOLVENT.

#### ##\$COMPOSIT\_BASE\_EXPNO =

(Table: NMR\_EXPERIMENT\_WS . COMPOSIT\_BASE\_EXPNO)

Virtual basis experiment number of a composit experiment. No spectra.

--> Generated by spectrometer.

--> Interpreted by SampleTrack.

When the composit experiment is restarted, this field is transfered to ICON-NMR<sup>™</sup>:

- --> Generated by SampleTrack
- --> Interpreted by spectrometer.

#### ##\$COMPOSITEXPERIMENT=

#### (Table: NMR\_EXPERIMENT\_WS . COMPOSIT\_EXPERIMENT)

Composit experiment name as defined in the spectrometer software.

This experiment is not necessarily defined within the SampleTrack experi

ment tables.

Format: string (max. 256 digits)

- --> Generated by spectrometer.
- --> Interpreted by SampleTrack.

#### ##\$COMPOSITMULTIORDERFILES

(Table: BSTS\_REGISTRY -- DEFAULT\_JCAMP\_VALUES - Value)

Switches between the possibility of decomposing a composite experiment within SampleTrack (1) or the spectrometer software (0). When the decomposition takes place in the spectrometer software (0), SampleTrack generates the order with one experiment.

--> Generated by SampleTrack.

--> Interpreted by spectrometer.

#### ##\$COMPOSITSEQNUMBER=

#### (Table: NMR\_EXPERIMENT\_WS . COMPOSIT\_SEQ\_NUMBER)

The composite experiment sequence number followed by comma (,) and the number of experiments belonging to the composite experiment.

Format: n,m

--> Generated by spectrometer.

--> Interpreted by SampleTrack.

#### ##\$DANGER= NOT USED

(Table: SAMPLE\_ORDERS . DANGER\_HINT)

Indicates dangerous attributes of a sample like

toxic, biohazard, radioactive, irritant, flammable, cancer hazard, ...

Format: string (max. 40 digits).

--> Possibly generated by SampleTrack.

#### ##\$DATASETNAME= NOT USED

Directory path name of the result file

Format: string (max. 40 digits).

--> Generated by SampleTrack.

--> Interpreted by spectrometer.

#### ##\$DISKUNIT=

(Table: NMR\_EXPERIMENT\_WS . DISKUNIT)

Target directory (default: /u)

Format: ASCII string (max. 50 digits).

--> Generated by SampleTrack .

--> Interpreted by spectrometer (ICON-NMR™).

##\$DISPOSE= NOT USED

Format: ASCII string (max. 40 digits).

--> Generated by SampleTrack.

--> Interpreted by AutoPREP.

##\$EXPERIMENT\_WS\_ID=

(Table: NMR\_EXPERIMENT\_WS . NMR\_EXPERIMENT\_WS\_KEY)

Unique experiment ID

Format: ASCII string.

--> Generated by SampleTrack

--> Interpreted by SampleTrack.

#### ##\$EXPERIMENT=

(Table: NMR\_EXPERIMENT\_WS . EXPERIMENT\_NAME\_KEY)

Experiment name: 13C,31P,1H,1R,GC-IR,GC-MS,MS,...

Format: ASCII string (max. 256 digits).

--> Generated by SampleTrack.

--> Interpreted by spectrometer.

##\$EXPERIMENTEND= <Stunde : Minute Monat / Tag / Jahr>

(Table: NMR\_EXPERIMENT\_WS . DATE\_END)

Format: ASCII string (max. 40 digits)

--> Generated by Spectrometer.

--> Interpreted by SampleTrack.

##\$EXPERIMENTNUMBER=

(Table: NMR\_EXPERIMENT\_WS . EXPERIMENT\_NUMBER)

ExpNo

Part of RESULTDATAPATH.

--> Generated by Spectrometer.

--> Interpreted by SampleTrack.

**Restart Experiment:** 

--> Generated by SampleTrack.

--> Interpreted by spectrometer.

##\$EXPERIMENTSTART= <Hour : Minute Month / Day / YEARr>

(Table: NMR\_EXPERIMENT\_WS . DATE\_START)

Format: ASCII string (max. 40 digits).

--> Generated by Spectrometer.

--> Interpreted by SampleTrack.

#### ##\$EXPERSTATUS=

(Table: NMR\_EXPERIMENT\_WS . STATUS)

Order status SampleTrack --> Spectrometer.

Possible entries: NEWPROBE or ADDITIONAL.

(SampleTrack --> Spectrometer).

measurement status Spectrometer --> SampleTrack.

Possible entries: ENDOK, ENDERR, WORKING.

(Spectrometer --> SampleTrack).

The error message can be attached in case of an error (ENDERR),

separated by a space.

Format: ASCII string (max. 40 digits).

--> Generated by SampleTrack or Spectrometer.

--> Interpreted by SampleTrack or Spectrometer.

#### #\$FILTER=

(Table: NMR\_PREPARATION\_WS . FILTER)

Indicates the use of a paper filter.

Possible entries: 0, 1.

--> Generated by SampleTrack.

--> Interpreted by AutoPREP.

##\$INSTRUMENT=

(Table: NMR\_EXPERIMENT\_WS . INSTRUMENT\_NAME\_KEY)

Indicates the analysis instrument for an experiment e.g. DPX300 Format: ASCII string (20 digits).

--> Generated by SampleTrack.

--> Interpreted by spectrometer.

#### ##\$JCAMPRESULTFILE=

field exists for compatibility reasons.

Use JCAMPRESULTFILE\_REAL instead.

(Table: NMR\_EXPERIMENT\_WS . JCAMP\_RESULT\_FILE)

Path and file name of processed data (REAL) in JCAMP format.

(source: [1D: 1r] [2D: still open]).

Format: ASCII string (max. 40 digits).

--> Generated by Spectrometer.

--> Interpreted by SampleTrack.

Example JCAMPRESULTFILE.

##\$JCAMPRESULTFILE\_RAW=

(Table: NMR\_EXPERIMENT\_WS . JCAMP\_RESULT\_FILE)

Path and file name of raw data (FID, SER) in JCAMP format.

(source: [1D: fid] [2D: still open]).

Format: ASCII string (max. 40 digits).

--> Generated by Spectrometer.

--> Interpreted by SampleTrack.

See example JCAMPRESULTFILE.

#### ##\$JCAMPRESULTFILE\_REAL=

(Table: NMR\_EXPERIMENT\_WS . JCAMP\_RESULT\_FILE)

Path and file name of processed data (REAL) in JCAMP format.

(source: [1D: 1r] [2D: still open]).

Format: ASCII string (max. 40 digits).

--> Generated by Spectrometer.

--> Interpreted by SampleTrack.

See example JCAMPRESULTFILE

#### ##\$JCAMPRESULTFILE\_REAL\_IMAG=

(Table: NMR\_EXPERIMENT\_WS . JCAMP\_RESULT\_FILE)

Path and file name of processed data (REAL and IMAG) in JCAMP format.

(source: [1D: 1r und 1i] [2D: still open]).

Format: ASCII string (max. 40 digits).

--> Generated by Spectrometer.

--> Interpreted by SampleTrack.

See example JCAMPRESULTFILE

#### ##\$JOURNAL=

(Table: METHODS . JOURNAL\_NUMBER)

Indicates the labor journal number.

Format: ASCII string (max. 40 digits).

--> Generated by SampleTrack.

--> Interpreted by spectrometer.

#### ##\$MAXTEMP=

(Table: NMR\_PREPARATION\_WS . MAX\_TEMP)

Maximum sample or solvent temperature.

Format: ASCII string (max. 40 digits).

--> Generated by SampleTrack.

--> Interpreted by the preparation robot.

#### ##\$MINTEMP=

(Table: NMR\_PREPARATION\_WS . MIN\_TEMP)

Minimum sample or solvent temperature.

Format: ASCII string (max. 40 digits).

--> Generated by SampleTrack.

--> Interpreted by the preparation robot.

#### ##\$NMRPARAMS=

(Table: NMR\_EXPERIMENT\_WS . PARAMS)

Experiment parameters seperated by semicolons.

Parameter name and its value are seperated by a blank.

Example:ns 20;sw 10.

Format: ASCII string (max. 40 digits).

--> Generated by SampleTrack.

--> Interpreted by the preparation robot.

#### ##\$ORDERNUMBER=

(Table: SAMPLE\_ORDERS . SAMPLE\_ORDER\_NUMBER\_KEY)

Internal unique order number.

Format: ASCII string (20digits).

--> Generated by SampleTrack.

-->Interpreted by SampleTrack.

#### ##\$PHDEMAND= NOT USED

(Table: NMR\_PREPARATION\_WS . )

Required PH value.

Format: ASCII string (max. 40 digits).

- --> Generated by SampleTrack .
- --> Interpreted by the preparation robot.

#### ##\$PREPARATION\_WS\_ID=

(Table: NMR\_PREPARATION\_WS . NMR\_PREPARATION\_WS\_KEY)

Unique key of the preparation record.

Format: ASCII string.

--> Generated by SampleTrack.

--> Interpreted by SampleTrack.

##\$PREPARATIONDATE= <Hour : Minute Montht / Day / Year>

(Table: NMR\_PREPARATION\_WS . DATE\_END)

End of preparation.

Format: ASCII string (max. 40 digits).

--> Generated by the preparation robot.

--> Interpreted by SampleTrack.

#### ##\$PREPARATIONOPERATOR=

(Table:NMR\_PEPARATION\_WS.USER\_NAME\_KEY)

(or NMR\_PEPARATION\_WS.PREPARATION\_OPERATOR\_NAME)

ID of the chemist doing the preparation.

Format: ASCII string (20 digits).

--> Generated by SampleTrack.

#### ##\$PREPSTATUS=

(Table: NMR\_PREPARATION\_WS . STATUS)

Order status preparation robot --> SampleTrack.

Possible entries: ENDOK, ENDERR, WORKING.

The error message can be attached in case of an error (ENDERR),

separated by a space.

Format: ASCII string (max. 40 digits).

--> Generated by the preparation robot.

--> Interpreted by SampleTrack.

#### ##\$PROTONEXCHANGE=

(Table: NMR\_PREPARATION\_WS . PROTON\_EX\_NAME\_KEY)

Indicates the solvent for a proton exchange.

Example: D2O, ...

Used by the spectrometer software by extending.

#### ##\$SAMPLETITLE with [##\$PROTONEXCHCHANGE]-Exchange for the plot

Format: ASCII string (20 digits).

--> Generated by SampleTrack.

--> Interpreted by spectrometer.

#### ##\$RESULTDATAPATH=

Example RESULTDATAPATH.

(Table: NMR\_EXPERIMENT\_WS . RESULT\_DATA\_PATH)

FID path name (raw data) after measurement.

A combination of several spectrometer variables like SPECTROMETER

OPERATOR, SAMPLEID and EXPERIMENTNUMBER.

Example: /u/data/guest/nmr/AB000134/10.

SPECTROMETEROPERATOR = guest.

SAMPLEID = AB000134.

EXPERIMENTNUMBER = 10.

Format: ASCII string (max. 40 digits).

--> Generated by Spectrometer.

--> Interpreted by SampleTrack.

##\$RETURNDATE= NOT USED

Format: <Monat/Tag/Jahr> ASCII string (max. 40 digits)

##\$ROUT\_POS= NOT USED

(NMR\_EXPERIMENT\_WS.WS\_POSITION)

Format: ASCII string (max. 80 digits).

--> Generated by SampleTrack.

--> Interpreted by SampleTrack.

#### ##\$SAMPLEAMOUNT=

(NMR\_PREPARATION\_WS . SAMPLE.AMOUNT)

Sample substance in mg.

Format: ASCII string (max. 40 digits).

--> Generated by SampleTrack .

--> Interpreted by the preparation robot.

##\$SAMPLECHANGERHOLDER=

(Table: NMR\_EXPERIMENT\_WS . SAMPLE\_CHANGER\_HOLDER)

Sample position within the B-ACS carousel.

Format: ASCII string (max. 40 digits).

--> Generated by Spectrometer (when reading the tubeID-Barcode).

or manually entered (MANUPREP) with urgent sample orders.

--> Interpreted by SampleTrack or Spectrometer (URGENT).

#### ##\$SAMPLEID=

(Table: SAMPLE\_ORDERS.SAMPLE\_ID)

Identification of the sample substance.

Format: ASCII string (10 digits).

--> Generated by SampleTrack.

--> Interpreted by spectrometer.

#### ##\$SAMPLENATURE= NOT USED

(Table: SAMPLE\_ORDERS.SAMPLE\_NATURE)

State of sample aggregation like solid, liquid, semifluid, crystalline.

Format: ASCII string (max. 40 digits).

--> Generated by SampleTrack .

##\$SAMPLETITLE=

(Table: SAMPLE\_ORDERS . SAMPLE\_TITLE)

Plot title.

Format: ASCII string (max. 240 digits).

--> Generated by SampleTrack.

--> Interpreted by spectrometer.

##\$SENSITIVE= NOT USED

Sample qualities like light, hygroscopic...

Format: ASCII string (max. 40 digits).

--> Generated by SampleTrack.

#### ##\$SETUPDATE= NOT USED

(Table: SAMPLE\_ORDERS.DATE\_ENTRY)

Order setup or entry date.

Format: ASCII string (max. 40 digits).

--> Generated by SampleTrack.

#### ##\$SETUP\_FILE=

ICON-NMR<sup>™</sup> uses this variable temporarily.

It holds the order file name in a report file.

Example: 1234.10r

Format: ASCII string (max. 40 digits).

--> Generated by Spectrometer.

--> Evaluation for debugging only.

#### ##\$SPECTROMETEROPERATOR=

(Table: NMR\_EXPERIMENT\_WS . INSTRUMENT\_LOGIN\_NAME)

Spectrometer User. If the variable #USER is empty, it fills up the

#RESULTDATAPATH with the UserID.

Format: ASCII string (20 digits).

--> Generated by Spectrometer.

--> Interpreted by SampleTrack.

#### ##\$SOLVAMOUNT=

(Table: NMR\_PREPARATION\_WS . SOLVENT\_AMOUNT)

Solvent amount in ml.

Format: ASCII string (max. 40 digits).

--> Generated by SampleTrack.

--> Interpreted by the preparation robot.

#### ##\$SOLVENT=

(Table: NMR\_PREPARATION\_WS . SOLVENT\_NAME\_KEY)

Used solvent ID like H2O,THF,DMSO,D2O,CDCL3,...

Format: ASCII string (max. 40 digits).

--> Generated by SampleTrack.

- --> Interpreted by preparation robot.
- --> Interpreted by spectrometer.

##\$SOLVMETHOD=

(Table: NMR\_PREPARATION\_WS .

SOLVENT\_MIX\_METHOD\_NAME\_KEY)

Used solvent method like USONIC, HEATER...

Format: ASCII string (max. 40 digits).

--> Generated by SampleTrack .

--> Interpreted by the preparation robot.

#### ##\$SOLVTEMP=

(Table: NMR\_PREPARATION\_WS . SOLVENT\_MIX\_TEMP)

Solvent temperature.

Format: ASCII string (max. 40 digits).

--> Generated by SampleTrack .

--> Interpreted by the preparation robot.

#### ##\$SOLVTIME=

(Table: NMR\_PREPARATION\_WS . SOLVENT\_MIX\_TIME)

Solvent duration.

Format: ASCII string (max. 40 digits).

--> Generated by SampleTrack .

--> Interpreted by the preparation robot.

#### ##\$TUBEHANDLING= NOT USED

Tube handling like 4mm,top melting, coding, vacuum, ...

Format: ASCII string (max. 40 digits).

--> Generated by SampleTrack.

#### ##\$TUBEID=

(Table: METHODS . TUBE\_ID)

Temporarily sample identifikation during the measurement:

It is the ID on the barcode collar tube. The tube ID is entered when the

preparation takes place, either manually, by hand scanner or with the B-BCE instrument.

Format: ASCII string (max. 40 digits).

--> Generated by the preparation robot.

--> Interpreted by SampleTrack und Spectrometer.

#### ##\$ URGENT =

(Table: SAMPLE\_ORDER . URGENT\_SAMPLE\_ORDER)

Urgent orders are measured before others.

Valid entries: 1 or 0

--> Generated by SampleTrack

--> Interpreted by spectrometer

#### ##\$ USER =

(Table: NMR\_EXPERIMENT\_WS .USER\_NAME\_KEY)

User ID - known in the spectrometer software and ICON-NMR™

Part of RESULTDATAPATH.

Format: ASCII string (20 digits).

--> Generated by SampleTrack.

--> Interpreted by spectrometer.

# 22

22.1

22.2.1

# Introducing Scripts for Automation

The package **stscript.bpl** (Module: dmScript.pas) is used for SampleTrack Client, Easy Dialog and Communication Server.

It includes routines for automation to store custom specific configurations.

Available commands are most of the commands used in Delphi 6.0


The following sections describe the common variables and functions.

Common Variables and Functions	22.2
--------------------------------	------

#### Variables with Instrument Type: Interface

vFileNameFile name for Import or Export (read/write).vMethodMethod Name, default is the content from vDefMethodName(usually 'NMR').vScriptName (RO)Name of the running scripts.vInstrumentName(RO)Name of the current instrument.vTransactionError(RO)Boolean-Variable within the program (StComSrv) that indicates a transaction error.


Figure 22.1. Instrument Type Interface

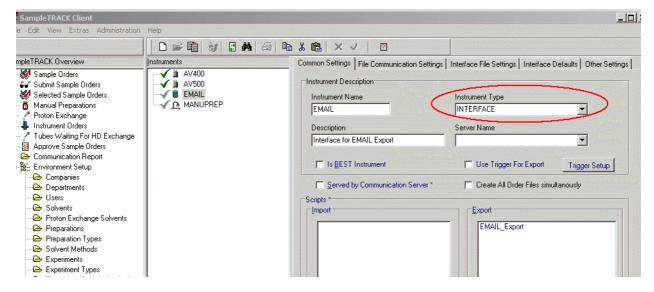

#### Columns on Tab 'Interface File Settings'

vClientIdCol (RO) vExperimentCol (RO) vHdExchangeCol (RO) vSampleIdCol (RO) vSampleTitleCol (RO) vSolventCol (RO) vPrepInstrCol (RO) vExpInstrCol (RO)

Figure 22.2. Interface File Settings

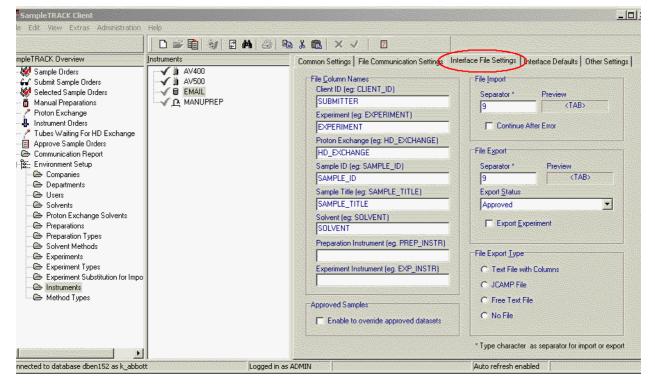

fClientId	Field content of vClientIdCol (global variable, editable)	
fExperiment	Field content of vExperimentCol (global variable, editable)	
fHdExchange	Field content of vHdExchangeCol (global variable, editable)	
fSampleId	Field content of vSampleIdCol (global variable, editable)	
fSampleTitle	Field content of vSampleTitleCol (global variable, editable)	
fSolvent	Field content of vSolventCol (global variable, editable)	
fPrepInstr	Field content of vPrepinstrCol (global variable, editable)	
fExpInstr	Field content of vExpInstrCol (global variable, editable)	
vDefSoStartStatus (RO)Start status Sample Order for the current instrument		

vDefMNmrStartStatus (RO)Start status methods (NMR), for the current instrument

vDefNpwStartStatus (RO)Start status Nmr\_Preparation\_Ws for the current instrument

vDefNewStartStatus (RO)Start status Nmr\_Experiment\_Ws for the current instrument

vDefMMsStartStatus (RO)Start status methods (MS) for the current instrument

vDefMpwStartStatus (RO)Start status Nmr\_Preparation\_Ws for the current instrument

vDefMewStartStatus (RO)Start status Nmr\_Experiment\_Ws for the current instrument

vDefCompanyName (RO)Company Name

vDefDepartmentName (RO)Department Name

vDefNmrExpInstrName (RO)Default NMR Spectrometer Instrument

vDefNmrPrepInstrName (RO)Default NMR Preparations Instrument

vDefMethodName (RO)Default Method

#### **Common Terms for Script Processing**

- SO 'SAMPLE\_ORDERS'
- M 'METHODS'

NPW 'NMR\_PREPARATION\_WS'

NEW 'NMR\_EXPERIMENT\_WS'

UNCHECKED'0'

#### CHECKED'1'

When a table is opened multiple times, the record can be identified using an alias name + '\_' + a consecutive number, rather than using table names.

Example: (always beginning with '1')

SAMPLE_ORDERS_1
SAMPLE_ORDERS_2
NMR_EXPERIMENT_WS_1
NMR_EXPERIMENT_WS_2
etc.

Functions for Script Processing

Procedure Self.Abort#

Stops the current processing.

Transaction and post\_transaction\_script will be performed.

22.3

22.3.1

Procedure Self.Exit#

Ends the current script. The script(s) that follow will be carried out.

Procedure Self.SetGlobVar#(VariableName, Value: String). Allows the definition of global variables valid within a list of scripts.

Function Self.GetGlobVar#(VariableName: String): String Delivers the value of a variable that was set previously. If not defined, the value returned is an empty String.

Function Self.ConvertToFieldName#(Text: String): String Converts a text into data base field name format and returns it as a String.

Function ShellExecute#(Operation, FileName, Parameter, Directory: String.

ShowMode: Integer): Boolean.

Executes a command.

**Operation** defines the action to be executed, e.g. **Open**. Filename is the name of the commando file or the file to be opened. **Parameter** are the parameters passed to **FileName**. **Directory** is the working directory. **Showmode** indicates the size of the new window.

1	= Minimized
2	= Maximized
No entry	= Normal

#### Functions for File Import/Export

22.3.2

Function Self.GetFileFieldCount#: Integer.
Returns the number of fields from the standard I/O file.
Function Self.GetFileFieldName# (Index: Integer): String.
Returns the column name from the I/O file.
Function Self.GetFileValue#(ColumnName: String): String.
Returns the column content from the I/O file.
Function Self.SetFileValue#(ColumnName, Value: String): Boolean.

Writes a value into the column of the I/O file.

Function Self.OpenJCAMPFileTable#(FileName, AliasTableName: String): Boolean.

Opens an optional JCAMP file. This file can be used additionally to the standard I/O file. Fields are handled as table fields.

Function Self.OpenTextFileTable#(FileName, AliasTableName: String.

HasHeadline: Boolean; Separator: Char): Boolean.

Opens a text file as a table.

This file can be used additionally to the standard I/O file.

Fields are handled as table fields.

Function Self.SetTextColumns#(Table, FieldNames: String): Boolean.

Defines, which columns to be used.

Column separator is ";".

Function Self.ReadTextLine#(Table: String; var Line: String): Boolean. Adds a line to the file.

Function Self.WriteTextLine#(Table, Line: String): Boolean. Writes a line to the file

#### **String Manipulation**

Function Self.Extract#(var OrgString: String; Sep: String): String. Extracts the next column field from a field list with column separator **Sep** 

# Functions for Tables

	Function Self.GetSequenceNumber(SequenceName: String): String.		
	Returns the next sequence number		
	Function Self.ExecStoredProc(aProc: String): String.		
	Returns the return value of a stored procedure		
	Function Self.OpenTable(Table, LookupColumnNames, LookupValues, OrderBy: String.		
ReadOnly: Boolean): Boolean.			
	Opens a table for repeated read or write.		
	The first record found with the where clause is the current record.		
	(Conditional on LookupColumnNames, LookupValues)		
	Return value: Record exists (True/False)		
	Function Self.CloseTable(Table: String): Boolean.		
	Closes the table		
	Function Self.GetTableState(Table: String): Integer.		
	Returns the table status:		
	0=EOF; 1=browse; 2=insert; 3=edit		
	Function Self.GetTableCountOf(Table: String): Integer.		
	Returns the number of tables in the list <b>aTable</b>		
	Function Self.ReadFromTable(Table, ColumnName: String.		
	var ResultString :String): Boolean.		
	Returns a value of an opened table. Additionally field contents from the ta- ble EXT_FIELDS can be returned.		
	Those fields must be predefined in the DB Registry.		
	Return value: Record exists (True/False)		
	Function Self.WriteToTable(Table, ColumnName, Value: String): Boolean.		
	Writes into an opened table.		
	Additional fields can be written into table EXT_FIELDS as defined in the DB Registry.		

Function Self.SaveTable(Table: String): Boolean. Saves the current Record into the Database

Function Self.InsertToTable(Table: String): Boolean. Saves the current Record into the Database

Function Self.Locate(Table, LookupColumnNames, LookupValues,

OrderBy: String): Boolean)

Jumps to the first record in a table with a hit on the where clause.

(Conditional on LookupColumnNames, LookupValues)

Function Self.CopyFields(SourceTable, DestinationTable: String): Boolean. Copies the content of all fields with the same column field from source table to destination table.

Destination table columns containing values will not be overwritten.

Function Self.DeleteRecord(Table: String; CloseTable: Boolean): Boolean. Deletes the current record and closes the table when CloseTable is TRUE.

Function Self.DeleteFromTable(Table, LookupColumnNames,

LookupValues: String): Boolean.

Deletes one or more records.

Function Self.SetMethodsStatus(aSampleOrderNumberKey, aStatus: String): Boolean.

Sets the status for all worksteps and methods.

Returns TRUE on success.

The Tube ID is released when the status >= 70 und status<>90.

Function Self.SetWSStatus(aWSTable, aWSKey, aStatus: String): Boolean.

Sets the status of the workstep.

Worksteps are preparations and experiments. The superior status must be set by the function RecreateStatus.

Function Self.RecreateStatus(aTable, aKey: String): Boolean.

Checkes the method and sample order status. Conditionally modifies the status when required.

Releases the tube ID when the status >= 70 and status <> 90.

Sample Order: self.Recreate(SO, '1234')

Methods: self.Recreate(M, '1234;NMR')

Function Self.GotoFirstRecord(aTable: String): Boolean.

Sets the cursor to the first record in the table aTable. The table must be opened before. Returns FALSE when the table was not found.

Function Self.GotoLastRecord(aTable: String): Boolean.

Sets the cursor to the last record in the table aTable. The table must be opened before. Returns FALSE when the table was not found.

Function Self.GotoNextRecord(aTable: String): Boolean.

Sets the cursor to the next record in the table aTable. The table must be opened before. Returns FALSE when the table was not found.

Function Self.GotoPrevRecord(aTable: String): Boolean.

Sets the cursor to the previous record in the table aTable. The table must be opened before. Returns FALSE when the table was not found.

Function Self. GotoRecordNumber (aTable: String; aLine): Boolean.

Sets the cursor to the specified record in the table aTable. The table must be opened before. Returns FALSE when the table was not found.

Function Self.BOF(aTable: String): Boolean.

Returns TRUE when when the first record is hit by the function GotoPrev-Dataset or is hit by the function GotoFirstDataset.

Returns TRUE when the file set is empty or the table was not found.

Function Self.EOF(aTable: String): Boolean.

Returns TRUE when when the last record is hit by the function GotoNext-Dataset or is hit by the function GotoLastDataset.

Returns TRUE when the file set is empty or the table was not found.

Function Self.RecordCount(aTable: String): Integer.

Returns the number of rows of a file set

Function Self.CancelTable(aTable: String): Boolean.

Cancels the changes

Function Self.RenameTable(aOldTable, aNewTable: String): Boolean. Renames the table for further processing

# Lookup Functions Database

Function Self.LookupTableValue(Table, LookupColumnNames, LookupValues,
ResultColumnName: String; var ResultString: String): Boolean.
Looks up a value in a table and closes the table
Returns: Record found (True/False)
Function Self.GetInstrumentDefaultValue(Instrument, Title: String.
var ResultString: String): Boolean.
Returns the title within the section I_ <instrument></instrument>
Returns: Record found (True/False)
Function Self.GetDefaultValue(Title: String; var ResultString: String): Boolean.
Returns the title content within the section I_ <aktuelles instrument=""></aktuelles>
Returns: Record found (True/False)
Function Self.GetGlobalDefaultValue(Method, Title: String;
var ResultString: String): Boolean.
Returns the title content from the section
GLOBALS or GLOBALS_ <method>.</method>
In case the String <b>Method</b> is empty, the section GLOBALS is inquired in the following.
Returns: Record found (True/False)
Function Self.ConvertToUNC(aInstrument, aPath: String): String.
Converts an instrument path to UNC
Function Self.GetDBTime(Dummy: Integer): String.
Returns the current database date and time
Format: '10:00 03/22/1999'
Function Self.GetDBRegistryEntry(aUser, aSection, aTitle, aDefault: String.
aField: Integer): String.
Returns a value from the DB Registry
aField indicates the kind of information: 0=Value, 1=Desc, 2=Memo1, 3=Memo2
Function Self.GetAliasName(aTableName, aFieldName: String): String.
Returns the column alias name

BRUKER BIOSPIN

# Field / Value Evaluation

	Function Self.IsValidString(aText: String; aUseUpperLetters, aUseLowerLe aUseFigures: Boolean; aCustomChars: String): Boolean; Checks for a valid String	tters,
	Function Self.IsValidFloat(aText, aRangeString: String.	
	aRangeIncrement: String): Boolean.	
	Checks for a valid floating point number	
	Function Self.IsValidInteger(aText, aRangeString: String;	
	aRangeIncrement: String): Boolean.	
	Checks for a valid number	
	Procedure Self.CheckValidFieldString(aTableName, aFieldName: String;	
	aUseUpperLetters, aUseLowerLetters, aUseFigures: Boolean;	
	aCustomChars: String).	
	Checks for a valid field String	
	Procedure Self.CheckValidFieldFloat(aTableName, aFieldName: String;	
	aRangeString: String; aRangeIncrement: String).	
	Checks for a valid field floating point number.	
	Procedure Self.CheckValidFieldInteger(aTableName, aFieldName: String;	
	aRangeString: String; aRangeIncrement: String).	
	Checks for a valid field number.	
Show Messages		22.

Procedure Self.ShowMessage(MessageString: String). Shows a message.

Procedure Self.ShowError(ErrorString: String).

Shows an error message. The script processing is aborted for the current loop.

(Internal Exception).

Procedure Self.ShowDatasetFields(aTable: String). Displays the current fields and their content.

# Function: Send Emails

Function Self.SetMailParams(aHost, aPort, aUserID, aCharSet: String): Boolean. Configures E-mail with common parameters.

Function Self.SetMailID(aSampleID: String): Boolean.

Places an identifier to send the E-mail automatically when the identifier changes.

Function Self.SetMailFrom(aFromAddress, aFromName, aReplyTo: String): Boolean.

Defines the E-mail sender

Function Self.SetMailTo(aToAddress, aToCarbonCopy, aToBlindCarbonCopy: String): Boolean.

Defines the E-mail receiver.

Function Self.SetMailSubject(aSubject: String): Boolean. Defines the E-mail title.

Function Self.SetMailHeader(aHeader: String): Boolean. Defines the E-mail header.

Function Self.SetMailFooter(aFooter: String): Boolean. Defines the E-mail footer.

Function Self.AddMailBody(aBody: String): Boolean. Defines an E-mail body block.

Function Self.ClearMailBody: Boolean. Deletes an E-mail body block.

Function Self.AddMailAttachment(aAttachment: String): Boolean. Adds an attachment to the E-mail.

Function Self.ClearMailAttachments: Boolean. Deletes the E-mail attachment list.

Function Self.SendMail: Boolean. Sends the E-mail.

Function Self.ClearMessage: Boolean. Deletes the complete E-mail content.

### Special Functions

22.9

Function Self.GetExperimentID(SO, NEW: String):String.

Returns a new experiment ID for an experiment.

The experiment type must be known, otherwise the sample ID is returned. The format is defined in the DB Registry:

Globals\Experiment\_ID\_Mask.

Function Self.CheckForDelayedExperiment(NEW: String):Boolean.

Checks if the experiment (current record) must be set to the status DE-LAYED. In that case the status and the start date is modified. The function returns TRUE.

Function Self.DoExperimentSubstitution(NEW, Experiment: String):Boolean.

Performs an experiment substitution if an appropriate value in the table Exp\_Substitions exists.

This Function may generate several experiments as defined in Exp\_Substitions.

Returns TRUE when a substitution takes place.

Function Self.DoHDExchange(NPW, HDInstrument, HDSolvent: String):Boolean.

Adds a preparation with HD Exchange (if not already existing).

Returns TRUE if already existing.

Function Self.CheckForCompositExperiments:Boolean.

Checks the internal dataset list for a composite experiment. If not, sub experiments are added.

Returns TRUE if COMPOSITMULTIORDERFILES=1

Function Self.ClearInvalidOrderFiles(aInstrument: String):Boolean.

Delete invalid order files for the given instrument. When the parameter alnstrument is empty all the order files for all the instruments of type **EXP** are checked for invalid orders.

Function Self.GetNewMethod(aSOKey, aMethodType: String): String.

Generates a new method of type aMethodType for the sample order key aSOKey.

Example: Self.GetNewMethod(1234, 'NMR')='NMR-2').

Function Self.CheckTubeIDStatus(aMethodTable: String): Boolean.

Checks, if the Tube ID is already used in the method table and inserts it where applicable.

## Function: Archiving

22.10

Function Self.GetCurrentDir(aDir: String; aLevel: Integer = 0): String. Returns the last directory name of a directory path name. aLevel > 0 indicates the number of parent directories up to the root. Example: C:\a\b\c\d\e getcurrentDir(Dir,0) returns e getcurrentDir(Dir,1) returns d\e

Function Self.GetParentDir(aDir: String; aLevel: Integer = -1): String.

Returns the complete path of the parent directory.

Example: Example: C:\a\b\c\d\e

GetParentDir(dir,1) returns C:\a\b\c\d

GetParentDir(dir,2) returns C:\a\b\c

Function Self.ArchiveData(aTable, aFieldName, aDestPath, aFileMask: String;

aDescend: Boolean): Boolean.

Archives all files in aFieldName. aFieldName can be a path or a list of file names. aDestPath indicates the destination path. aFileMask indicates the filter for the files to be moved. aDescend indicates, if underlying directories must be archived.

Caution: All the data is deleted in the source directories!

Function Self.CopyData(aTable, aFieldName, aDestPath, aFileMask: String;

aDescend: Boolean): Boolean.

Copies all files in aFieldName. aFieldName can be a path or a list of file names. aDestPath indicates the destination path. aFileMask indicates the filter for the files to be copied. aDescend indicates, if underlying directories must be copied. Copy Data can be used as backup routine.

Function Self.DeleteDataFiles(aTable, aFieldName, aFileMask: String;

aDescend: Boolean): Boolean.

Deletes all files in aFieldName. aFieldName can be a path or a list of file names. aFileMask indicates the filter for the files to be deleted. aDescend indicates, if underlying directories must be deleted.

#### Caution: Data are deleted!

Function Self.DeleteDataDirs(aTable, aFieldName, aFileMask: String.

aDescend: Boolean; aParentDir: Integer = 0): Boolean.

Deletes all directories when empty. aFileMask indicates the filter for directories to be deleted. aDescend indicates, if underlying directories must be deleted. aParentDir indicates if parent directories should be deleted. aParentDir is the first parent directory.

#### Caution: Data are deleted!

Function Self.ExecuteArchivingForSO(SOKey, Config: String): Boolean.

Archives SampleOrder with the key SOKey with archiving Method Config.

#### Caution: This function works only within Import or Export Scripts.

## **Obsolete Functions**

22.11

Function Self.DeleteDataset(Table: String): Boolean. Deletes the current record and closes the table. Replaced by: Self.DeleteRecord verwenden.

Function Self.GotoFirstDataset(aTable: String): Boolean.Sets the cursor to the first record in table aTable.The table must be opened before. Returns FALSE if the table is not found.Replaced by: Self.GotoFirstRecord

Function Self.GotoLastDataset(aTable: String): Boolean.Sets the cursor to the first record in table aTable.The table must be opened before. Returns FALSE if the table is not found.Replaced by: Self.GotoLastRecord

Function Self.GotoNextDataset(aTable: String): Boolean.Sets the cursor to next record in table aTable.The table must be opened before. Returns FALSE if the table is not found.

Replaced by: Self.GotoNextRecord

Function Self.GotoPrevDataset(aTable: String): Boolean.

Sets the cursor to the previous record in table aTable.

The table must be opened before. Returns FALSE if the table is not found. Replaced by: Self.GotoPrevRecord

Function Self.WriteFileHeader(Header: String): Boolean.

Writes the file title.

Use Self.OpenTextFileTable and Self.WriteTextLine for Exports except for JCAMP and Table Files.

Function Self.WriteFileLine(Line: String): Boolean.

Adds a line at EOF.

Use Self.OpenTextFileTable and Self.WriteTextLine for Exports except for JCAMP and Table Files.

## Alphabetical List of Function and Procedures

22.12

Abort AddMailAttachment AddMailBody ArchiveData BOF CancelTable CheckForCompositExperiments CheckForDelayedExperiment CheckTubeIDStatus CheckValidFieldFloat CheckValidFieldInteger

**BRUKER BIOSPIN** 

ClearInvalidOrderFiles ClearMailAttachments ClearMailBody

ClearMessage CloseTable ConvertToFieldName ConvertToUNC CopyData CopyFields DeleteDataDirs DeleteDataFiles DeleteFromTable DeleteRecord DoExperimentSubstitution DoHDExchange EOF ExecStoredProc ExecuteArchivingForSO

Exit

Extract GetAliasName GetCurrentDir GetDBRegistryEntry GetDBTime GetDefaultValue GetExperimentID GetFileFieldCount GetFileFieldName GetFileValue GetGlobalDefaultValue GetGlobVar GetInstrumentDefaultValue

GetParentDir

GetSequenceNumber

GetTableCountOf

GetTableState

GotoFirstRecord

GotoLastRecord

GotoNextRecord GotoPrevRecord

GotoRecordNumber

InsertToTable

IsValidFloat

IsValidInteger

Locate

LookupTableValue

OpenJCAMPFileTable

OpenTable

OpenTextFileTable

ReadFromTable

ReadTextLine

RecordCount

RecreateStatus

RenameTable

SaveTable SendMail

SetFileValue

SetGlobVar

SetMailFooter

SetMailFrom

SetMailHeader

SetMailID

SetMailParams

SetMailSubject SetMailTo SetMethodsStatus SetTextColumns SetWSStatus

ShellExecute ShowDatasetFields ShowError ShowMessage WriteTextLine WriteToTable

# System Requirements

# 23

**General Hint** 



For server, PC and spectrometer software requirements see also the BRUKER SampleTrack<sup>™</sup> homepage for Server and Client PC requirements: <u>http://</u>www.bruker-biospin.de/NMR/automat/samtrac/samtrac5.html

### Server Requirements

23.2

23.1

At least one server is required exclusively for SampleTrack<sup>™</sup> based on the following minimum requirements:

Pentium IV, 2 GHz (or faster), 1 GB RAM or better.

60 GB harddisk or larger. As an option an additional harddisk for mirroring. This does not include storage for spectra data.

100/1000 Mbit/s Ethernet connection.

A Windows® 2000 compatible graphics adapter with minimum 1024x768 pixel resolution.

A 17" (or larger) monitor with at least 1024x768 pixel resolution and 75 Hz vertical frequency.

A Windows® 2000 compatible mouse and keyboard.

Microsoft Windows® 2000 / 2003 Server.

ORACLE® Standard Edition 8.1.7 or higher (can be delivered by BRUKER).

# **Client PC Operation System Requirements**

SampleTrack<sup>™</sup> can be installed on existing spectrometer PC's or other compatible systems based on the requirements below:

Pentium IV, 1GB RAM or better.

2 GB free space on harddisk (minimum)

100/1000 Mbit/s Ethernet connection.

A Windows® 2000 compatible graphics adapter with minimum 1024x768 pixel resolution.

A 17" (or larger) monitor with at least 1024x768 pixel resolution and 75 Hz vertical frequency.

A Windows® 2000 compatible mouse and keyboard.

Microsoft Windows® 2000 / 2002 Professional.

ORACLE® Client 8.1.7 or higher (can be delivered by BRUKER).

## Spectrometer Software Requirements

- \* XWIN-NMR<sup>™</sup> 3.1 patch level 11 or higher or TOPSPIN® 1.2 or higher
- \* ICON-NMR<sup>™</sup> 3.1 or higher

#### Directory Sharing with UNIX Instrument Computers

In order for data to be exchanged between the instruments in your laboratory and the SampleTrack<sup>™</sup> data management system, several defined directories have to be shared. This definition may be handled by various NFS software products or using the SAMBA daemon on your UNIX computer. For installation and support please contact your network administrator.

23.4

23.5

# Oracle® 9 Server and Client Installation

24

24.1

# **ORACLE® 9i Database Server Installation**

Refer to your Oracle® documentation and provider information for the database installation. The following section may help to prevent common errors that may occur during installation.

For the Oracle® 9i client installation see <u>"ORACLE® 9i Client Installation"</u> on page 186.

The following installation screenshots are dedicated for the database server PC.

Insert the  $\ensuremath{\mathsf{Oracle}}\xspace{\mathbbmath{\mathbb{R}}}$  CD into CDROM drive - the following Welcome dialog will appear



Figure 24.1. The Oracle® Welcome Window

Accept the name OraHome92 and the path C\oracle\ora92



Figure 24.2. The Oracle® File Locations Window

Choose Oracle® 9i database

Figure 24.3. The Oracle® Available Products Window



Select Standard Edition



Figure 24.4. The Oracle® Installation Type Window

#### Select General Purpose

Figure 24.5. The Oracle® Database Configuration Window



Select standard port number 2030

Figure 24.6. Microsoft® Transaction Server Window



#### Select Gobal Database Name dbstrack and SID dbstrack

Figure 24.7.	The Oracle®	Database	Identification	Window
--------------	-------------	----------	----------------	--------

Oracle Unive	ersal Installer: D	atabase Identification	-	<u> </u>
Database	ldentificati	on		
	abase is uniquely id I Database Name fo	entified by a Global Databas r this database.	e Name, typically of the forr	n "name.domain".
Global Databas	e Name: dbstrack			
instance on this		one Oracle9i instance whic cle System Identifier (SID). prefer.		
SID:	dbstrack			
Exit	Help	Installed Products	Previous	Next


Select database file directory C:\oracle\oradata



Figure 24.8. The Oracle® Database File Location Window

Use the default character set

Figure 24.9. The Oracle® Database Character Set Window

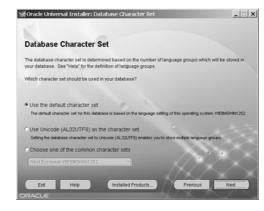

Before the installation a summary of the selected items is shown

Figure 24.10. The Oracle® Summary Window



Installation is running



Figure 24.11. The Oracle® Installation Status Window

The configuration assistant lets you modify the sys and system passwords (recommended)

Figure 24.12. The Oracle® Database Configuration Assistant .

🛱 Database Configuration Assistant	×
Database creation complete. Check the logfiles at C:\oracle\admin\dbstrack\create for details.	
Database Information: Global Database Name: dbstrack System Identifier(SID): dbstrack Server Parameters Filename: C:\oracle\ora92\database\spfiledbstrack.ora	
Change Passwords	
For security reasons, you must specify a password for the SYS and SYSTEM accounts in the new database.	
SYS Password:	
Confirm SYS Password:	
SYSTEM Password: ++++++++	
Confirm SYSTEM Password: +++++++	
Note: All database accounts except SYS, SYSTEM, DBSNMP, and SCOTT are locked. Select the Password Management button to view a complete list of locked accounts or to manage the database accounts. From the Password Management window, unlock only the accounts you will use. Oracle Corporation strongly recommends changing the default passwords immediately after unlocking the account. Password Management)	
(OK)	

Database installation has finished



Figure 24.13. The Oracle® End of Installation Window

Find the SampleTrack database within the Enterprise Manager Console


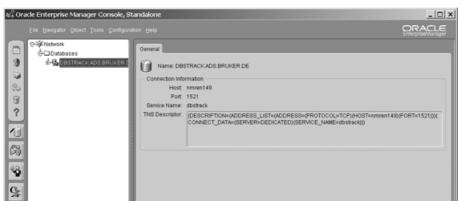



Figure 24.14. Enterprise Manager Console, Standalone Window

The Path in the Environment System

24.1.1

The path "C:\oracle\ora92\bin" must be in the environment system variable, see advanced system properties

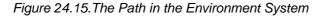

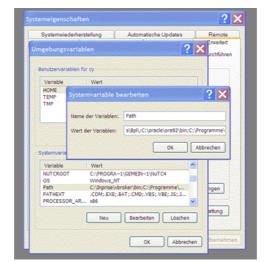

## **ORACLE® 9i Client Installation**

Installing Oracle® 9i client on a Windows® PC

Refer to your Oracle® documentation and provider information for the database installation. The following section may help to prevent common errors that may occur during installation.

Version 8 and SampleTrack installation:

For the installation of Oracle® 8 client installation see the file <u>Getting Start-</u> <u>ed.pdf</u>.

The following installation screenshots are dedicated for each PC, where the Oracle® 9i client together with SampleTrack client is needed.

(For the installation of SampleTrack client installation see the file <u>Getting</u> <u>Started.pdf</u>.

Insert the Oracle® CD into CDROM drive - the following Welcome dialog will appear



Figure 24.16.The Oracle® Welcome Window

Accept the name OraHome92 and the path C\oracle\ora92

Figure 24.17.The Oracle® File Locations Window



#### Select the Oracle® client installation

Figure 24.18. The Oracle® Available Products Window



Select the installation type Administrator

Figure 24.19.The Oracle® Installation Type Window



Before the installation a summary of the selected items is shown



Figure 24.20. The Oracle® Summary Window

# Oracle® 8 Server and Client Installation

25

Oracle® 8.1.7. Database Server on a Windows PC

Refer to your Oracle® documentation and provider information for the database installation. The following section may help to prevent common errors that may occur during installation.

# Installation Method

25.2

25.1

Insert the Oracle® CD into CDROM drive - the following dialog will appear



Figure 25.1. The Oracle® Installation Window:


**Choose: Install Products** 

🔭 Oracle Un	versal Installer	
	File Locations	
	Source Enter the full path of the file representing the product(s) you want to install:	
	Path: E:\stage\products.jar Browse	
	Destination Enter or select an Oracle Home name and its full path:	
	Name: OraHome81	
	Path: D:toracletora81	
	About Oracle Universal Installer	)
Exit	Help Installed Products Previous Next	


Figure 25.2. The File Locations Window

Press: Next

Figure 25.3. The Loading Product Information Window

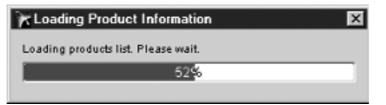

Wait...

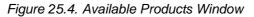



Select: Oracle® Enterprise Edition, then press Next



Figure 25.5. Installation Types Window

Choose: Typical Press: Next

🔭 Oracle Un	iversal Installer		
	Database Identification		
	An Oracle8i database is uniquely identified "name.domain". Enter the Global Databas		lly of the form
	Global Database Name: dbstrack	-	
	<ul> <li>A database is referenced by at least one O other instance on this computer by an Orac</li> </ul>		
	been entered which you can accept or char		led of D flas
	SID: dbstrack		
Exit	Help Installed Produ	cts Previous	Next

Figure 25.6. Database Identification Window

For Global Database Name and Instance Name enter: dbstrack

Figure 25.7. Summary Window



A summary of the products to be installed is shown.

A summary of the products to be installed is displayed. Press Install to continue.

🔭 Oracle Universal Installer \_ 🗆 X Install 0 talling Oracle Enterprise nager Common Files 2.2.0.0.0 Onaste WeisbB to: op web b acting vt/update\_res\_r ethods\_time estamp htm 35% y and access v (Cancel) id ma ou can find a log of this install session at: C:Program FilesOracleUnvertoryVogsVinstallActions.log you install another product, the above log will be copied to C:Program FilesOracleUnvertoryVogsVinstallActions200 10: 102-06-25\_10-05-26-AMJo

Figure 25.8. Install Progress Window


please wait ...

Wait...

Database creation completed will appear.



Figure 25.9. Oracle® Database Configuration Assistant Alert Window

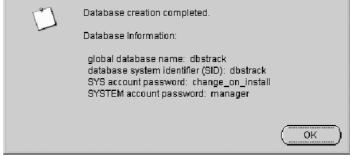

Press OK to end installation

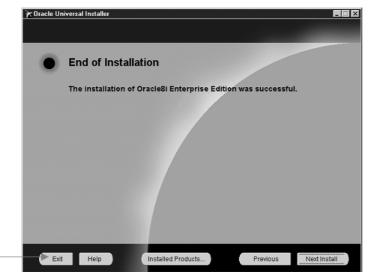


Figure 25.10.End of Installation Window

Press exit to end installer

Press Exit to close the installer.

# Installing Oracle 8.1.7. Client on a Windows PC

25.3



Figure 25.12.Welcome Window

🔭 Oracle Uni	iversal Installer	
	Welcome	
	The Oracle Universal Installer guides you through the Installation and configuration of your Oracle products.	
	Click "Installed Products" to see all installed products.	
	Deinstall Products	
	About Oracle Universal Install	er)
Exit	Help Installed Products Previous Next	)


Press Next

🔭 Oracle Univ	versal In	staller	_ 🗆 🗙
	File	Locations	
	Source Enter th	e he full path of the file representing the product(s) you want to install:	
	Path:	D:\oracle817\stage\products.jar Browse	
		nation	
	Name:	r select an Oracle Home name and its full path: OraHome81	
		OraHome81	
	Path:	D:\oracle\ora81 Browse	)
		About Oracle Universal Instal	ler
Exit		Help Installed Products Previous Next	

Figure 25.13. File Location Window

Select the source and destination, then press Next



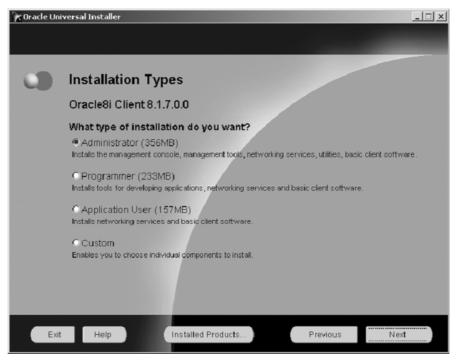

Select either "Administrator" or "Application User", press Next

Figure 25.15.Summary Window



A summary of the programes to be installed is shown.

Press Install to start the installation.

Press OK to end installation.

...

# *Client/Server Network Connection*

# Net8 Assistant

26

26.1


There are several possibilities in creating client/server network connections

Figure 26.1. Client Server Network Connection

	à.	Open Office Document		HP DLA	۲			
4	77			HP RecordNow	۲.			
110	l	WinZip	G.	Oracle - OraHome81	۱G	Application Development	F	
	•	-		Oracle Installation Products	<b>ا</b> ل	Database Administration	F	
8 <b>-</b>	141	Programs •	۲	SampleTRACK	<u>ا</u>	Enterprise Manager	F	
<b>.</b> 2 3	-	Documents +			G	Extended Administration	F	
ž.	_				C.	Network Administration	Þ	* Enterprise Login Assistant
1 H L	5	Settings •			Œ	] Oracle Internet Directory	•	Microsoft ODBC Administrator
٦ đ	Ē.	e			Œ	Quick Tours	•	g 🐔 Net8 Assistant
8	2	Search •			Q.	Release Notes	•	🥳 Net8 Configuration Assistant
	9	Help			ø	Installation Guide and Release Notes		Oracle ODBC Help

The best one is the Net8 Assistant.

Figure 26.2. Service Naming

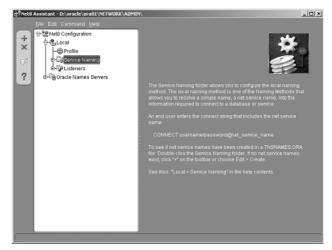


Mark "Local/Service Naming" and click the plus symbol.

Figure 26.3. Net Service Name Window




Enter the netservice name "dbstrack".

Figure 26.4. Protocol Window

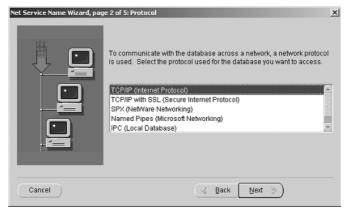


Select the network protocol TCP/IP.

#### Figure 26.5. Host Name Window

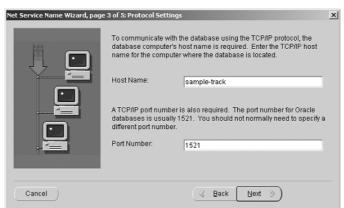

Enter the host name of the computer where the Oracle Database has been installed.

Figure 26.6. Service Name Window

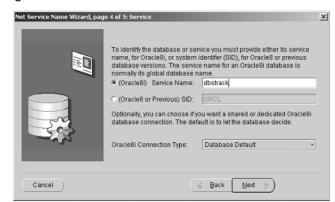

For Oracle 8.1.7 or higher use the Oracle 8i Service Name: "dbstrack"

Figure 26.7. Test Window

Net Service Name Wizard, page	5 of 5: Test	×
	Press Test if you would like to verify that the database can be reached using the information provided. When you are finished, or if you want to skip testing, press Finish to create the net service name or, if enabled, Next to continue. Test	
Cancel	( Back Next > Finish	5

Press Finish to finish the installation.

Figure 26.8. Result Window



After pressing **Finish** the results should look like those above.

# SampleTrack™ System and Client Installation

# 27

## SampleTrack<sup>™</sup> System vs. SampleTrack<sup>™</sup> Client

27.1

There are only a few differences between the SampleTrack<sup>TM</sup> System and the SampleTrack<sup>TM</sup> Client Installation. With the system installation the SampleTrack Communicator moves to the Autostart folder and starts automatically with each PC start. The Communicator often runs on a PC where the database was installed. The PC running the Communicator must have network connection to the database and to all served spectrometer PC's. ST Config is an additional program for the administrator that controls basic connectivity parameters <u>"ST Config - A Program for the Administrator" on page 214</u>.



Figure 27.1. The SampleTrack Installation CD

# SampleTrack<sup>™</sup> System and Client Installation

# Installation of the SampleTrack<sup>™</sup> System

#### Figure 27.2. The SampleTrack Installation Wizard

Run the installation program for SampleTrack called "stinstXXX\_XX.exe" which is placed on the SampleTrack-CD. (The actual setup file is named: "stinst242\_18.exe")

₿Bruker SampleTRACK V2.42	2.18 - InstallShield Wizard
	Welcome to the InstallShield Wizard for Bruker SampleTRACK V2.42.18
	The InstalShield(R) Wizard will instal Bruker SampleTRACK Y2.42.18 on your computer. To continue, click Next.
	WARNING: This program is protected by copyright law and international treaties.
	<back cancel<="" th=""></back>

Run the SampleTrack installation program "stinstXXX\_XX.exe" which is located on the SampleTrack CD.

Press Next to continue.

Enter the customer user name and organization as shown in the figure below:

Bruker S						
Custome	r Information					Citing of the
Please e	nter your informa	tion.				1000
User Na	ne:				_	
Bruker I	Instruments					
Organiza						
Bruker	Instruments					
Instal th	is application for:					
	Anyone w	ho uses this c	omputer (all use	rs)		
	C Only for g	je (Bruker Ins	truments)			
				_		
	ampleTRACK	/2.42.18 - 1	< <u>B</u> ack		Next >	Cancel
Destinati	ampleTFIACK \ on Folder It to install to this Instal Bruker Sa	folder, or dic	nstallShield V k Change to ins /2.42.18 to:	/izard		ar.
Destinati	ampleTRACK \ on Folder at to install to this	folder, or dic	nstallShield V k Change to ins /2.42.18 to:	/izard		×
Destinati	ampleTFIACK \ on Folder It to install to this Instal Bruker Sa	folder, or dic	nstallShield V k Change to ins /2.42.18 to:	/izard		ar.
Destinati	ampleTFIACK \ on Folder It to install to this Instal Bruker Sa	folder, or dic	nstallShield V k Change to ins /2.42.18 to:	/izard		ar.
Destinati	ampleTFIACK \ on Folder It to install to this Instal Bruker Sa	folder, or dic	nstallShield V k Change to ins /2.42.18 to:	/izard		ar.
Destinati	ampleTFIACK \ on Folder It to install to this Instal Bruker Sa	folder, or dic	nstallShield V k Change to ins /2.42.18 to:	/izard		ar.
Destinati	ampleTFIACK \ on Folder It to install to this Instal Bruker Sa	folder, or dic	nstallShield V k Change to ins /2.42.18 to:	/izard		ar.
Destinati	ampleTFIACK \ on Folder It to install to this Instal Bruker Sa	folder, or dic	nstallShield V k Change to ins /2.42.18 to:	/izard		ar.
Destinati	ampleTFIACK \ on Folder It to install to this Instal Bruker Sa	folder, or dic	nstallShield V k Change to ins /2.42.18 to:	/izard		ar.
Destinati	ampleTRACK V on Folder et to install to this Install Bruker SI C: (Program File	folder, or dic	nstallShield V k Change to ins /2.42.18 to:	/izard		ar.

#### Figure 27.3. Customer Information Window


Enter user name and organization. Select installation option. Press Next.

Change destination folder if required.

Press Next.


Choose Install SampleTrack System and press Next


Figure 27.4. Setup Type Window

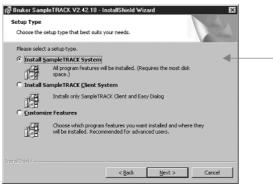

Choose Install SampleTRACK System and click next

Ready to install SampleTrack

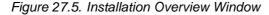

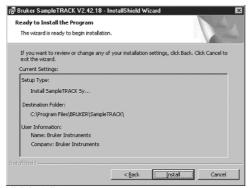

click install

Press Install.

Wait....

#### Finish the InstallShield Wizard

18 Bruker SampleTRAUK V2	.42.18 - InstallShield Wizard	×
	InstallShield Wizard Completed	
A	The InstallBheld Witzard has successfully installed Bruker SampleTRACK V2.42.18. Click Finish to exit the witzard.	
	<back cancel<="" finish="" th=""><th>1</th></back>	1

Figure 27.6. InstallShield Wizard Completion Window

please wait..

Press Finish.

#### SampleTrack<sup>™</sup> Patches

27.3

Installing the newest SampleTrack Patch.

SampleTrack patches (stbinxxx\_xx.exe) contain all the binary files of a complete SampleTrack installation.

They are usually placed in the directory "C:\Program Files\bruker\Sample-Track\bin\" .

To install the patch copy the file (stbinXXX\_XX.exe) to the SampleTrack directory

"C:\Program Files\bruker\SampleTrack\".

The start the executable file.

The new files will be copied to the ./bin directory. It is allowed to remove, rename or delete the old bin-directory before installing the patch to get rid of any invalid files.

🚔 C:\Program Files\BRUKER	\S 💶 🗙
<u>File E</u> dit <u>V</u> iew <u>H</u> elp	
SampleTRACK	- E ä
Din 242_18 data dbstruct Help Manual stbin242_24.exe	
1 object(s) selected 5.82	мв //

Figure 27.7. Installing the SampleTrack Patch

27.4

The actual SampleTrack patch is also available on the Bruker FTP-server: ftp://ftp.bruker.de/pub/nmr/SAMTRACK/INSTALL/V2.X

The SampleTrack Client checks the availability and the installation state of the SampleTrack database when it is first run.

In this description the behavior of the SampleTrack client is shown when running the first time on an "empty" Oracle® database.

Such a system is present if Oracle® was installed as described in the first chapter.

If running the SampleTrack client for the first time the following message should appear:

Figure 27.8. Information Window

Informati	on 🗵
<b>(</b>	Program not configured
	ОК

Press OK.

Figure 27.9. Database Export Window

)atabase Expert	X
_	Welcome to the SampleTRACK System. Please follow the next steps to prepare SampleTRACK for automation features. Important: ORACLE must be installed to proceed with the following steps:
	Please select one of the following options:
9	C greate a new database for this SampleTRACK system
	connect to an giving SampleTRACK system database
Create New Database	
View Summary	Berk Concel

select "create a new database for SampleTrack System" and click next

Select Create a new database for SampleTrack System and press Next.

Enter the Standard Oracle® Host String - dbstrack

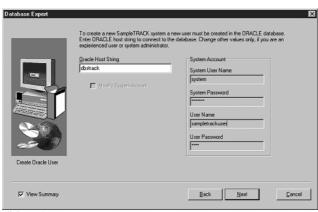


Figure 27.10.Standard Oracle® Host String

click next

#### Press Next.

A SQL script loads the default values into the database.

#### Figure 27.11.SQL Script Window



enter "exit"

Enter Exit.

When the storage of the standard values was completed successfully the following window will appear:

Figure 27.12.Confirmation Window

	Creation of the new user was successful. Now tables for SampleTRADK will be created. Please press button (Nexb) to start creation of tables.
Ceate Database Tables	
View Summary	Back Cancel

click next

Press Next.

Figure 27.13.Command Prompt Window



enter exit

Enter Exit.

Load SampleTrack components manually

Figure 27.14.Load Components Window

atabase Expert		×
_	SampleTRACK system tables are available. Load a SampleTRACK configuration to st within SampleTRACK.	art automation
	Please select one of the following options:	
-	C load sample configuration	
	C import detabase from file	
	C load SampleTRACK components manually	-
	C start with gmpty database	
Start SampleTRACK		
View Summary	Back New	Çancel

select "load SampleTrack components manually" and click next

Select the items that you want installed from the Component Selection Window.

Database Expert	Components below can be loaded into SampleTF existing changed data can lead to unpredictable		eep in m	ind that li	oading	items in
	Components					
	Name	Currer	nt Ne	w 🔺		
	Required data     DB Start Data	V2.02	4 V2.	024		
	Script Library		V4.	1		
And the second s	Default Filter Configuration					
	AutoPREP Sample Configuration					
	Delete All Tables Before Importing Data					
Load SampleTRACK	Name	0	Lurrent	New		
Components	AutoPREP Sample Configuration					
	Spectrometer Sample Configuration					
	BEST Spectrometer Sample Configura	noite			-	
	Print Sample Order Details Sample Cor	nfiguration				Cancel
View Summary						
View Summary	HTML Email Notification Sample Confi					Gancer
View Summary	HTML Email Notification Sample Confi Plain Text Email Notification Sample C					Zanco

Figure 27.15.Component Selection Window

select the items as shown

The items selected depend on the user's instrument environment.

The SQL statements output will be displayed:

Database Expert x Results ٠ 7 lines updated in 'ESTS\_RE mporting tables finished. mporting tables ..... elected items will be imported. In morting data 'Easy Dialog Sample 'BSTS\_REGISTRY Importing Easy Dialog pdated in 'H tables finis itens will be inported. data 'Default Filter Co ele Importing onfiguratio rtir be be be Ī 1 Next

Figure 27.16. Output of SQL Statements

finally the output of the used SQL-statements is shown (for checking only) click next

Press Next.

This completes the database installation.


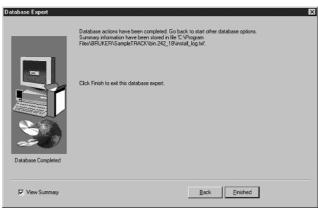


Figure 27.17. Program Completion Window

click finished

Press Finished.

The changes will be available with the next login.

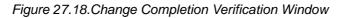

I he following windows come i	up	_ <i>8</i>   ×
SampleTRACK Client  SampleTRACK Client  SampleTRACK Overview	Sample IACK Client	
	Auto refrech disabled	
Page 13 Sec 1 13/13 At 4.3" Ln 5 0	tol 1 REC TRX EXT OVR LCK	

click ok

Press OK.

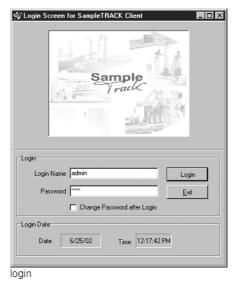


Figure 27.19. The Login Window

This completes the procedure.

🞯 SampleTRACK Client \_ 🗆 🗡 <u>E</u>dit ⊻iew E<u>s</u>tras Filter R Not yet finished samples 💌 🗅 🛩 🛍 🖐 🖤 🖬 🗛 😂 🖬 📼 ampiet RACK Overview

Senset Doctor

Ampiet Construction

Ampiet

Ampi Experi... Sample Hold Client ID Instrument Solvent Tube ID Worksteps • se Logged in as ADMIN Auto refresh enabled o database dbstrack as sample done

Figure 27.20. The Sample Track Client Window


# SampleTrack<sup>™</sup> Program Icons

27.5

After the system restart the following new entries will be found:

Figure 27.21.SampleTrack Program Icons



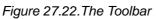


al 1914: 1914: 1914: 1914	Windows Update Windows Update New Office Document Open Office Document WinZp	대 HP RecordNow 대 Oracle - OraHome81 대 Oracle Installation Products	> > >	
Windows 2000 Professional @ 1월: 🔊 🔊 🖓 🕞 街 🕮 🗗 🖽	<ul> <li>Minutial Space</li> <li>New Office Document</li> <li>Open Office Document</li> <li>WinZp</li> <li>Brograms</li> <li>Rocuments</li> <li>Settings</li> <li>Seargh</li> <li>Heb</li> <li>Bun</li> <li>Shut Down</li> </ul>	WinZip  Acrobat Reader 5.0  HP DLA  HP RecordNow  Gr Oracle - OraHome81	• • •	
Sta	6 0 6 5 1			
		/		
Start	1000000 🖬 🔎 🛛			0.440.000 P

It is recommended that you place the icons ST Client and ST EasyDialog for each user into the toolbar by pressing theControl key and moving the icon.



The program ST Configuration tool is meant for use by the administrator only.

# ST Config - A Program for the Administrator

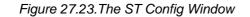

SampleTRACK Options		
Connectivity		
Database Network Configuration         C SQLNet 1.x (Oracle 7.1.x)         C SQLNet 2.x (Oracle 7.3.x)         Net8 (Oracle 8 and higher)         Override BDE Settings         Oracle Service Name         dben152	Login Settings Oracle Database Username  demo_labor Oracle Password 	•
Info	OK Cancel	Apply

Important connectivity settings are set in this stand-alone program. It is identical to the tab of the SampleTrack options found within the client menu - Administration - SampleTrack Options - Tab Connectivity.

The entries here play a role when a database connection error occurs after the installation.



Figure 27.24. The SampleTrack Administration Menu Option

27.7

## Database Connection Error After New Installation

This is a common problem with the BORLAND® BDE (Borland® Database Engine). The easiest way to handle the connectivity settings is by clicking the button **Override BDE Settings** see the figure <u>"The ST Config Window" on page 214</u> within the SampleTrack Connectivity window - either within the ST Configuration program or under SampleTrack client - Administration - Options.

You can also change the BDE settings in the control panel

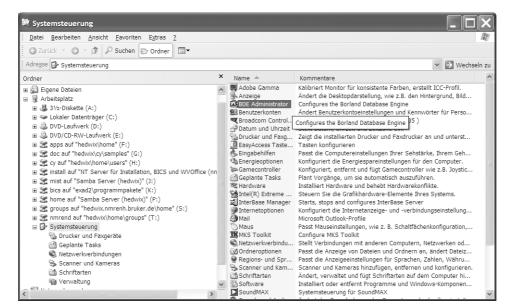


Figure 27.25.BDE Administrator Window

On the tab Configuration - Drivers - Native - ORACLE change the entry in the OR-ACLE definition window DLL32 to "SQLORA8.DLL" and the VENDOR INIT to "OCI.DLL".

bject <u>E</u> dit <u>V</u> iew O <u>p</u> tions <u>H</u> elp			
plet fut view officius Helb			
≥ X ∽ a			
Drivers and System	Definition of ORACLE		
Databases Configuration	Definition		
∃- 🔁 Configuration			
≓-On Drivers		4.0	
- On Native		SERVER	
@ PABADOX		SQLORA8.DLL	
@ DB2		OCIDLL	
O DBASE	VENDORMAN	OGLEEL	
OFOXPRO	TRACE MODE	0	
- O INFORMIX	I I I VIGE MODE		
- 🚱 INTRBASE	BATCH COUNT	200	
G MSACCESS		32	
MSSQL	11	64	
ORACLE		FALSE	
O SYBASE		FALSE	
🗈 🔁 ODBC	ENABLE SCHEMA CACHE		
🗄 💻 System	LANGDRIVER		
		NONE	
		-1	
		TNS	
		FALSE	
		READ/WRITE	
		20	
	SCHEMA CACHE DIR		
		8	
		-1	
		ORA SERVER	
		SHARED AUTOCOMMIT	
		SERVER	
		MYNAME	

Figure 27.26.BDE Administrator Window: DLL32 and VENDOR INIT

Apply your changes and login to the SampleTrack Client.

## Installation of Instruments

This section describes how to install the different types of instruments. SampleTrack supported instruments include:

- NMR spectrometer using BACS with barcoded tubes
- NMR spectrometer with BEST and barcoded Gilson racks
- Bruker AutoPrep (ASP)
- Gilson Preparation Robots controlled by BRUKER software
- Other Preparation Robots using the provided interface
- Bruker SampleRail<sup>™</sup> System
- Interfaces to other data management systems

27.8

The Environment Setup Window shows the installed instruments.

SampleTRACK Client Fie Edt View Extras Administration H amplet FRACK Overview

amplet FRACK Overview

Sample Driders

Sample Driders

Sample Driders

Sample Driders

Manual Proparations

Manual Rack Preparations

Proton Exchange

Communication Report

Comparies

Departments

Departments

Departments

Departments

Departments

Departments

Department Steps

Submits

Preparation Types

Submits

Submits

Department Syste

Submits

Department Vector

Department

Department Vector

De D≌∎ ∛ 54 8 8 8 × √ 8 Common Settings | File Communication Settings | Other Settings | MANUPREP Instrument Descrip Instrument Type PREP Instrument Name MANUPREP ¥ scription anual Preparation • Is BEST Instrument Use Trigger For Export Trigger Setup Served by Communication Server.\* Create Al Order Files simultanously Script Exp ent Substitution for Imported Fi struments od Type Add Edit Delete Add Edit Delete Debug Scripts Email " DB Registry entries are displayed in blue ted to database dbstrack as sar Auto refresh enabled

Figure 27.27. Environment Setup Window


The only instrument installed here is called MANUPREP which symbolizes the manual preparation work center.

Installation of BRUKER NMR Spectrometer Using B-ACS	27.9
---	------

```
1. Installation of a Dummy Instrument from a Template.27.9.1
```

- Open the SampleTrack client
- Select menu point Administration DB Management

Figure 27.28. The Load SampleTrack Components Window
--

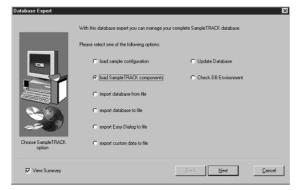

Choose Load SampleTrack Components and press Next.

 Obtahese Expect
 Components below can be loaded into SampleTRADS. Heave keep in mind that loading keem in sching blanged data can lead in urpredictable reads.

 Filter Strange Configuration
 Components below can be loaded into SampleTRADS. Heave keep in mind that loading keem in Sching blanged data can lead in urpredictable reads.

 Filter Sample Configuration
 Components below can be loaded into Sample Transport

 Filter Sample Configuration
 Spectnometrs Sample Configuration


 Filter Sample Configuration
 Spect configuration

 </tbo

Figure 27.29. Spectrometer Configuration Window

Select Spectrometer Sample Configuration and Press Next.

Figure 27.30.SQL Output Window

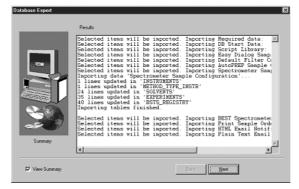

Press Next, Finished and finally login to SampleTrack client.

#### 2. Make a Copy of the Dummy Instrument "SPECT"


27.9.2

Name it in accordance to the spectrometer type (e.g. AV400, AV400\_BEST, DPX300\_1, DPX300\_2, ...)

SampleTRACK Client		동안 지수는 것을 것을 것 같은 것을 것 같은 것	
<u>File Edit View Extras Administration Help</u>			
	D ≫ 🛍 🦭 🖬 é	5 Ba & Ba × V   D	
SampleTRACK Overview	Instruments	Common Settings   File Communication Settings	Other Settings
Some Difference Sources     Sources Sources Sources     Sources Sources Sources     Manual Preparations     Manual Preparations     Manual Preparations     Manual Preparations     Proton Exchange     Instrument Orders     Active Rackup Sample Order Data     Comparise     Active Rackup Sample Order Data     Comparise     Comparise     Sources     S	SPECT	Communication Settings [ Trie Communication Settings ] Instrument Name SPECT Description Bruker Spectrometer  Fills gEST Instrument  Fills gEST Instrument  AVANCE_Jmport  Add Edt Detete  Debug Scripts  DB Registry entries are displayed in blue	Instrument Type EXP Server Name Use Trigger For Export Create AI Order Files simultanously Create AI Order Files simultanously Export AVANCE_Export AVANCE_Export Email
Connected to database dbstrack as sampletrackuse	Looped in as ADMIN	Auto refresh enabled	
uonnected to database dostrack as sampletrackuse	Loggeo in as ADMIN	Auto resean enabled	

Figure 27.31. The Instrument Setup Window

Figure 27.32. The New Instrument Name Window

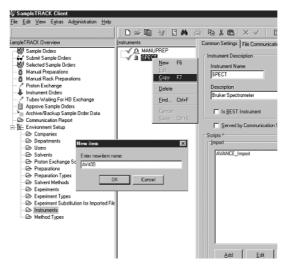

Enter the name of the new intrument (use capital letters) and click OK.Now click on the green check mark to store. The instrument will now appear in the list. Enter the import and export scripts by clicking the **Add** buttons and select the corresponding script from the list.

SampleTRACK Client			_ 🗆 🗙
<u>File Edit View Extras Administration Help</u>			
	D ≈ B   3/ E A   6	៑ Pa i @ × ✓   □	
SampleTRACK Overview	Instruments	Common Settings File Communication Settings	Other Settings
Somele Drees     Somele Drees     Somele Drees     Somele Drees     Manual Preparations     Manual Preparations     Manual Reparations     Manual Reparations     Manual Reparations     Manual Preparations     Mathematication Report     Mathematication Report     Solvents     Mathematication Report     Solvents     Solvents     Solvents     Solvent Methods     Expointert Types     Mathad Types	MANUPPEP	Instrument Description Instrument Name Av400 Description Buker Spectrometer  Geved by Communication Server * Google * Import Av4ACE_Import Av4ACE_Import Av4ACE_Import C Debug Scripts * DB Registry entries are displayed in blue	Instrument Type EXP Server Name Use Trigger For Export Trigger Setue Create AI Order Files simultanousty Export AVANCE_Export AVANCE_Export Ex
Connected to database dbstrack as sampletrackuse	Looped in as ADMIN	Auto refresh enabled	

Figure 27.33. The Import and Export Scripts Window

This completes this procedure.

27.10

Install the program stinstxxx\_xx.exe from the SampleTrack installer CD.

Figure 27.34. The Explorer Window

Ele     Ede     Yew     Favorites     Tools     Help       ↓ + Back • → • •     •     •     •     •     •       Address     •     •     •     •     •	
Address aspletrack	
	1 4 4
	• @@
Image: Stop     Image: Stop       Image: Stop <th></th>	
Type: Application Size: 15.8 MB 📃 My Computer	r //

The InstallShield Wizard will lead you through the installation.

Bruker SampleTRACK V2.42.18 - InstallShield Wizard

 Welcome to the InstallShield Wizard for
Bruker SampleTRACK V2.42.18

 The InstallShield(R) Wizard will install Bruker SampleTRACK
V2.42.18 on your computer. To continue, click Next.

 WARNING: This program is protected by copyright law and
international treaties.

 Velocity install reactions.

Figure 27.35.InstallShield Wizard Welcome Window

Press Next.


Enter your user and company name

謬Bruker SampleTRACK ¥2.42.18 - InstallShield Wizard Customer Information Please enter your information.	X
User Name:	
administrator	
Organization:	
BRUKER BIOSPIN	
Install this application for :	
<ul> <li>Anyone who uses this computer (all users)</li> </ul>	
Only for <u>m</u> e (administrator)	
InstallShield	
< <u>B</u> ack <u>N</u> ext	> Cancel

Figure 27.36.InstallShield Wizard Window

Commit or change the default destination folder.

Figure 27.37. The Destination Folder Window

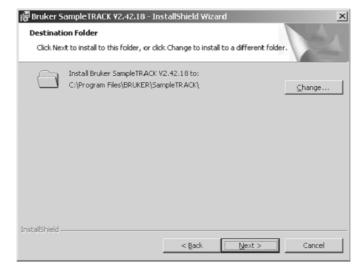

Choose "Install Sampletrack Client System".

Figure 27.38. The Setup Type Window

📳 Bruker SampleTRACK ¥2.42.18 - Ins	tallShield Wiza	rd	×
Setup Type Choose the setup type that best suits y	our needs.		
Please select a setup type.			
O Install SampleTRACK System			
All program features will space.)		quires the most disk	¢
Install SampleTRACK Client System	stem		
Installs only SampleTRA	ICK Client and Eas	sy Dialog	
C <u>C</u> ustomize Features			
Choose which program will be installed. Recom			re they
InstellShield			
	< <u>B</u> ack	<u>N</u> ext >	Cancel

Press Install to start the installation.

Figure 27.39. The Install Window

🖟 Bruker SampleTRACK V2.42.18 - InstallShield Wizard	×
Ready to Install the Program The wizard is ready to begin installation.	
If you want to review or change any of your installation settings, click Back. Click Cancel to exit the wizard.	
Current Settings:	
Setup Type:	
Install SampleTRACK Cli	
Destination Folder:	
C:\Program Files\BRUKER\SampleTRACK\	
User Information:	
Name: administrator	
Company: BRUKER BIOSPIN	
InstallShield	1

SampleTrack client installation is completed.

Figure 27.40.InstallShield Wizard Completed Window



## SampleTrack™ Web Client Installation

#### Introduction / System Requirements

The SampleTrack Web Client is a web service which offers easy access to the SampleTrack database without installing any client software on the customer's PC. The web service module is only available on

Windows® NT systems

(Windows® 2000, Windows® XP, Windows® 2003)

as a Microsoft<sup>®</sup> Internet Information Service Module or Apache Module. The Apache Webserver is automatically installed during the installation of Sample-Track. The IIS has to be installed and configured manually.

Depending on the web service usage, the server should be at least a Pentium IV with minimum 512 MB RAM.

Since the configuration must be done directly on the Web server, the PC should be available directly or via a terminal session.

Installation		28.2

Installation of Program Files


Install the SampleTrack software package on the server, that will be used as a **Web Server** later on. Be sure to select **Customize Features** and activate the item **Web Interface**.

Custom Setup	Sec. 1
Select the program features you want installed.	
lick on an icon in the list below to change how a feat	ure is installed.
D Optional	Feature Description
Applications	SampleTRACK Communicator
Archiving	
X V Instrument Simulator	
Label Designer	This feature requires 12KB on
	your hard drive.
Titration	
Web Interface	

Figure 28.1. InstallShield Wizard Custom Setup

28.2.1

After the SampleTrack installation the Web Interface module is still not activated. Login to the SampleTrack client as administrator. In the menu « Administration » choose "Configure Plug-Ins".

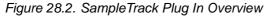


ver Plugins ampleTRACK Archiving Package	< <tr>         Vp         Down</tr>	Available Plugme SampleTRACK Atchiving Package SampleTRACK Barcode Pinting Package SampleTRACK Bart Package SampleTRACK Easy Dialog Package SampleTRACK Excel Report Package SampleTRACK Katk Support Package
nt Plugin: ampleTRACK Matrix Support Package ampleTRACK Easy Dialog Package ampleTRACK Barcode Printing Package ampleTRACK Archiving Package ampleTRACK BEST Package	<	SampleTRACK Pooling Package SampleTRACK Titation Package GempleTRACK Web bonilouration Reskone
y Dialog Plugins		
y Dialog Hugins	<	
y Dialog Plugins	< >	

Select the Web Configuration Package and add it to the "Client Plugins".

Close and restart the SampleTrack client. In the menu "Administration", select the new option "Configure Web Interface...".

The new configuration program will be displayed and shows a group tree on the left with different configuration sections. When a configuration section is selected, the current configuration is shown on the right side. Use the button "Change Data" to change the view.

Figure 28.3. Group Tree Window

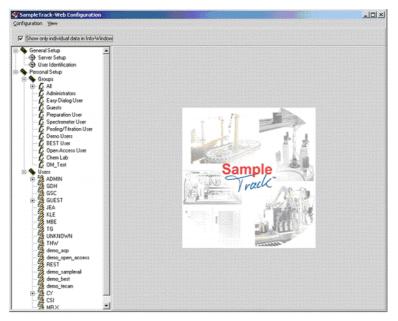

To install and start the Web Server choose "General Setup" and "Server Setup".

In the setup you must start the service as an Apache module or IIS module.

When selecting the IIS module be sure that IIS has been installed correctly before.

Click "Install" to install the service.

From this configuration window you can conveniently start and stop the Web service.

After successful installation a status window displays "Service started". Clicking on "Browser" opens the browser window with the SampleTrack Web client start page.

Secure	Socket	Layer	(SSL-HTTP)
--------	--------	-------	------------

It is possible to use the SampleTrack Web service with a secure connection (SSL).

This is independent from the SampleTrack configuration and involves the Apache or IIS server configuration. Please contact your IT department for more information.

#### Configuration

The SampleTrack Web configuration menu offers more options to customize your automation and individual interface. Apart from the "Custom-Page-Wizard" each SampleTrack client also provides individual adaption possibilities.

#### **User Identification**

The "User Identification" and "General Setup" define the kind of authentification.

With anonymous login (Autologin) a login is not required. To use the Autologin enter the user name which is used for logging into the system.

Session management: Here the session time, length a session should stay open, and the maximun number of concurrent sessions can be defined.



Hint for changing the password: When the SSL is not used, passwords are transferred unencrypted. This is also true, when passwords are being changed, even when "password encryption" was activated before.

28.3

28.3.1

28.2.3

# Login-Configurations Encrypt Password during Login (Unix-DES) Autologin Enable Autologin Username Save Default Values Session Management Session-Timeout Max Sessions [20] [min] Image: Save Changes exit

#### Figure 28.4. Login Configurations

#### Group or Individual User Settings

All further modifications to the HTML pages are group or user specific.

Select the user or group profile from the group tree on the left side. For changes concerning all users select the group "All".

When a group has no sub-details, this indicates that no specific modifications have been made for the group.

The configuration can be modified with a right click on the group and by selecting **Edit this item**.

#### **Navigation Board**

The configuration menu "Navigation" has options effecting the look and feel of the naviagtion board.

Here you can activate or deactivate the different views, or add additional links, e.g. to your own Intranet homepage. If you do not use a matrix system, you may deactivate this point.

Edit Navigation-Elem	ents				
Main Navigation					
Lounge and New	r				
✓ Sample Orders ✓ Instrument Orders					
✓ Matrix Orders					
Custom Pager					
Name		Enabled			
Custom Links					
Name	Link				
Bruker Homepage	http://www.bruke			Add	
<ul> <li>ST S0AP Interface</li> </ul>	nap://nmren166.	nmrenh bruker. de/st		Delete	

Figure 28.5. Navigation Elements Window

#### Layout

If you have modified the HTML template pages (see the section HTML packages) to adapt them to your corporate design, the layout section allows you to customize the internal design components. You can also define the color to highlight the sample orders.

#### Custom-page Wizard

The custom page wizard helps you to integrate your own HTML pages and views into the system.

Principally a HTML file is needed. It may have pictures inside. The program is optimized for HTML pages generated with Microsoft® Word ("Save as Web page").

To add a new HTML page enter a page title (it will be displayed within the navigation board) and select the HTML page.

The "Advanced Settings" options define the way the new page should be integrated:

Use **Open in new window** to open the page in a new window without the SampleTrack heading and navigation.

When using more complex HTML pages, it may happen, that the view is not the same as the original design. Try to solve the problems using the checkboxes **De-lete VML-Tags** and **Only Body&Head**.

Be certain that SampleTrack Web Module path is set correctly, e.g. when using an Apache server: "C:\Programme\BRUKER\SampleTRACK\web\apache\modules".

Select **Import** to import the HTML page.

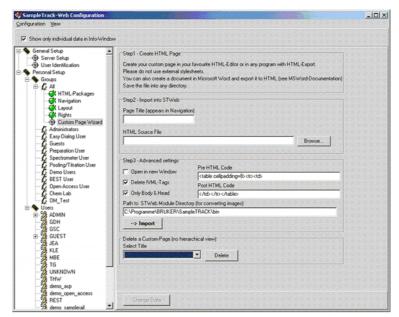


Figure 28.6. SampleTrack Web Configuration Window

#### HTML Packages

The SampleTrack Web Service are delivered as packages. While the package "Standard" consists of the necessary basics, more custom specific packages can be added. To import a package select the ".inf" file in the package main directory. Always install the standard package first, and afterwards any additional packages, so they override standard package files.

With a double click on a file you can view and edit the HTML code. After modification of a file the package status changes to "Dev" (development) and the original version is stored, so that a redo of the last modification is always possible.

To uninstall a package, right click on the file name and select "delete package" from the context menu.

Use the selection "Export Files" to export the files and edit them with a HTML editor.

Use the selection "Import file" to import new files into the system.



Remember, that a lot of HTML files include control commands for the visualization of database information. This information should not be changed when editing the file.

#### Rights

When editing the rights for the SampleTrack Web Client you can release Easy-Web Dialogs for a user, activate quick menus / Workstep Tool Boxes, mask and unmask sub menus.

In the Web configuration window all available EasyWeb Dialogs are listed. Select one and click edit.

In the window that appears you can modify the title, the authorization and the executable scripts.

In the window "Custom Rights" authorization can be changed by checking or uncheck the rights. At this development level there is only one authorization to activate or deactivate: "Applications". It allows or forbids access to view experiment results.

Easy-Dialog name	Enabled	Caption	Scripts	
ed_dialog_standard ed_dialog_standard_d2o	Yes Yes	Standard Standard with D20	-None- -None-	
Edititem ustom Rights Name of Right	Nr Status	Quickme		
<ul> <li>Applications</li> </ul>	013 Enabled	Me	mple Orders thods iparations beriments	
		Cor	fig Quickmenus	

#### SampleTrack<sup>™</sup> Web Client Installation

# Figures

1 Copyr	ight Notices	9
2 Warnii	ngs and Notes	11
3 Introd	uction to Bruker LIMS	13
Figure 3.1.	Laboratory Data Management - LIMS	13
Figure 3.2.	SampleTrack <sup>™</sup> - Worksteps	14
Figure 3.3.	The SampleTrack <sup>™</sup> Interfaces	
Figure 3.4.	SampleTrack <sup>™</sup> as Data Reference	
Figure 3.5.	Daily Routine Laboratory Functions	
Figure 3.6.	Communication Overview	
Figure 3.7.	SampleTrack <sup>™</sup> Workflow Architecture	
Figure 3.8.	A Look into the Future	20
4 Introd	uction to the SampleTrack™ Client?	21
Figure 4.1.	Login Window	21
Figure 4.2.	Tree Overview	22
Figure 4.3.	The Main Window	23
Figure 4.4.	Double Click to See the Sample Order Details	24
Figure 4.5.	Preparation Details	24
Figure 4.6.	Experiment Details	25
Figure 4.7.	Add Preparation / Experiments / Proton Exchange Buttons	
Figure 4.8.	Sample Order Window Buttons	
Figure 4.9.	Items Contracted	
-	Items Expanded	
•	Status Information	
	List of all Status Symbols	
	Filter & View Assistant Button	
-	Filter and view management	
-	The Object Manager	
-	Filter Management	
Figure 4.17.	Filter Management	32
5 Functi	ons of the SampleTrack™ Client Module	33
Figure 5.1.	Scenario 1	33
Figure 5.2.	Scenario 2	33
Figure 5.3.	Scenario 3	34
Figure 5.4.	Status Overview	34
Figure 5.5.	The Submit Button	34
Figure 5.6.	Select All	
Figure 5.7.	Tree - Manual Preparation Area	
Figure 5.8.	The Manual Preparation Window	35

Figure 5.9.	B-BCE for Easy Tube ID Barcode Reading	
Figure 5.10.	The Proton Exchange Window	
Figure 5.11.	The Finished Button	
Figure 5.12.	Example of a Manual Implemented Finish	
Figure 5.13.	Status Icon of Waiting / Delayed Experiments t	
Figure 5.14.	Delay Experiments	
Figure 5.15.	The Start Delayed Experiments Button	39
Figure 5.16.	Highlight and Start Delayed Experiments	39
Figure 5.17.	Tubes Waiting for HD Exchange	40
Figure 5.18.	The HD Exchange Window	40
Figure 5.19.	The Tubes Waiting for Exchange Window	
Figure 5.20.	The Approve Window	
	B-BCE	
Figure 5.22.	The Environment Setup Area	43
6 Workin	a with the SampleTrack™ Client	45

#### 6 Working with the SampleTrack<sup>™</sup> Client

#### 45

Figure 6.1.	SampleTrack Order	45
Figure 6.2.	Entering the Client ID	
Figure 6.3.	Defining a New Method Using Right Mouse Button	46
Figure 6.4.	Entering the Journal Number	46
Figure 6.5.	The Edit Sample Order Details Window	47
Figure 6.6.	Select a Preparation	47
Figure 6.7.	Order Details	48
Figure 6.8.	Add Experiment	48
Figure 6.9.	Add D2O for a Proton Exchange Experiment	49
Figure 6.10.	Copying an Existing Experiment	50
Figure 6.11.	A Sample Order Created Using the Copy Function	50
Figure 6.12.	The Submit Window	51
Figure 6.13.	Rerun an Experiment	52
Figure 6.14.	Adding an Additional Experiment	53
Figure 6.15.	Example of a Filter Definition	54
Figure 6.16.	Search in View - Client Tool Window	55
Figure 6.17.	Search in DB - Sample Order Details Window	55

#### 7 Interaction with Analytical Applications

Figure 7.1.	The Paperclip Icon	57
Figure 7.2.	Interaction with TOPSPIN®	57
Figure 7.3.	The Results Data Path	58
Figure 7.4.	Interaction with AMIX <sup>™</sup>	58
Figure 7.5.	MDL® ISIS/Draw <sup>™</sup> Template	59
Figure 7.6.	The More Button	59
Figure 7.7.	Creating a New UNC Formatted Link to a Document File	60
Figure 7.8.	Creating a New Report	60
Figure 7.9.	An Example of a Report	61
Figure 7.10.	Administration - Application Management	62
Figure 7.11.	Defining the Application Type	62
Figure 7.12.	Example 1: Select Application	63
Figure 7.13.	Example 2: Select Application	63

#### 8 EasyDialog

57

Figure 8.1.	EasyDialog Icon	.65
•	EasyDialog Login	
Figure 8.3.	EasyDialog Example	.66
Figure 8.4.	Entering Data with EasyDialog	.67
Figure 8.5.	EasyDialog Confirmation Message	.67
Figure 8.6.	Automatic Reading of the Barcode Tube ID	.67

#### 9 SampleTrack™ Web Client

69

Figure 9.1. Figure 9.2.	SampleTrack Web Client Login	
Figure 9.3.	Sample Orders - View	71
Figure 9.4.	Complete Order	71
Figure 9.5.	Web - Structure and Spectra Viewer	72
Figure 9.6.	Web - Structure and Spectra Viewer, Zooming In	73
Figure 9.7.	Web - Structure and Spectra Viewer, Reset Image	73
Figure 9.8.	Search	74
Figure 9.9.	Extended Search	74
Figure 9.10.	Instrument Orders	75
Figure 9.11.	Matrix Orders	75
Figure 9.12.	EasyWeb Dialog	76
Figure 9.13.	Applications	77
Figure 9.14.	Application Wizard	77
Figure 9.15.	The View Spectra Window	78

#### 10 Interaction with Automation Robots

# Figure 10.1. Install PlugIns80Figure 10.2. Pooling default values81Figure 10.3. Protein Screening / Module 'Pooling'83Figure 10.4. Setting hits83

#### 11 Environment Setup

Figure 11.1. Environment Setup86Figure 11.2. New Department87Figure 11.3. New User88Figure 11.4. Message: Delete?89Figure 11.5. New Solvent90Figure 11.6. New Proton Exchange Solvent91Figure 11.7. New Preparation Method92Figure 11.8. New Solvent Mixing Method93Figure 11.9. New Experiment94Figure 11.10. Experiment Substitution Table95

#### 12 The Object Manager

Figure 12.1.	Menu Object Manager	98
Figure 12.2.	Object Manager Main Window	98
Figure 12.3.	Maintaining Groups	99
Figure 12.4.	Content Sub-Window	100
Figure 12.5.	Default Groups	100
Figure 12.6.	Group Tree	101
Figure 12.7.	Sub-categories	102

79

85

97

Figure 12.8. Adding a New User (Or Group of Users)	103
13 Example: Create a User Group	105
Figure 13.1. Create a Client Profile	105
Figure 13.2. Program Message	
Figure 13.3. Adding Table Default Values for a Group	107
Figure 13.4. Defaults in the Content Window	107
Figure 13.5. Example: Default Project Number	107
Figure 13.6. Drag and Drop EasyDialog Configuration to a Group	
Figure 13.7. Save EasyDialog for a Specified User / Group	
Figure 13.8. Drag and Drop Filter Configuration to a Group	
Figure 13.9. User Level in the Insert New Group Window	110
14 Handling the Tube ID	111
15 System Parameters	113
16 Shortcuts in SampleTrack™	115
17 SampleTrack™ Tables	117
Figure 17.1. Field Overview	118
18 EasyDialog Configuration	123
Figure 18.1. The Start EasyDialog Designer Menu Option	123
Figure 18.2. The EasyDialog Configuration Overview Window	123
Figure 18.3. Open / Create a New Form	124
Figure 18.4. Drag & Drop Objects	
Figure 18.5. The Drag & Drop Objects Window	
Figure 18.6. Drag - Drop Objects	
Figure 18.7. Changing the Grid Distance	
Figure 18.8. Configuring a New Object for EasyDialog	
Figure 18.9. The Field Overview Window	
Figure 18.10. Default Field Value in the Fields Overview Window	
Figure 18.11.Default Value in the Fields Overview Window	
Figure 18.12. Default Value in EasyDialog	
Figure 18.13.Entering a Default Status Value	
Figure 18.14.The Administrator Toolbar	
Figure 18.15.A New Input Form	
Figure 18.16.Drag and Drop Group Boxes	
Figure 18.17.Drag and Drop Field Sample ID	
Figure 18.18.Adjusting the Size of an Object	
Figure 18.19.The complete Group Box Figure 18.20.Adding Group Boxes	
Figure 18.21.Adding Buttons in a Form	
Figure 18.22.Save EasyDialog configuration	
19 E-Mail Support	137

Figure 19.2. Selecting HTML or Plain Text E-mail Notification	
Figure 19.3. The E-mail Trigger Setup Window	
Figure 19.4. E-mail Parameters	
Figure 19.5. E-mail Address in the Environment User Setup	142
20 Data Exchange & Directory Structure	143
21 JCAMP Definition File	145
22 Script Library	159
Figure 22.1. Instrument Type Interface	159
Figure 22.2. Interface File Settings	160
23 System Requirements	177
24 Oracle® 9 Server and Client Installation	179
Figure 24.1. The Oracle® Welcome Window	179
Figure 24.2. The Oracle® File Locations Window	
Figure 24.3. The Oracle® Available Products Window	
Figure 24.4. The Oracle® Installation Type Window	
Figure 24.5. The Oracle® Database Configuration Window	
Figure 24.6. Microsoft® Transaction Server Window	
Figure 24.7. The Oracle® Database Identification Window	
Figure 24.8. The Oracle® Database File Location Window	
Figure 24.9. The Oracle® Database Character Set Window	
Figure 24.10.The Oracle® Summary Window Figure 24.11.The Oracle® Installation Status Window	
Figure 24.12. The Oracle® Database Configuration Assistant	
Figure 24.13. The Oracle® End of Installation Window	
Figure 24.14.Enterprise Manager Console, Standalone Window .	
Figure 24.15. The Path in the Environment System	
Figure 24.16.The Oracle® Welcome Window	
Figure 24.17.The Oracle® File Locations Window	
Figure 24.18.The Oracle® Available Products Window	187
Figure 24.19.The Oracle® Installation Type Window	187
Figure 24.20.The Oracle® Summary Window	187
25 Oracle® 8 Server and Client Installation	189
Figure 25.1. The Oracle® Installation Window:	
Figure 25.2. The File Locations Window	190
Figure 25.3. The Loading Product Information Window	
Figure 25.4. Available Products Window	
Figure 25.5. Installation Types Window	
Figure 25.6. Database Identification Window	
Figure 25.7. Summary Window	
Figure 25.8. Install Progress Window	
Figure 25.9. Oracle® Database Configuration Assistant Alert Win	
Figure 25.10.End of Installation Window Figure 25.11.The Install Window	
1 1901 C 20.11.111C 11131all WIIIUUW	190

**BRUKER BIOSPIN** 

Figure 25.12.Welcome Window	195
Figure 25.13. File Location Window	196
Figure 25.14.Installation Types Window	196
Figure 25.15.Summary Window	197

#### 26 Client/Server Network Connection

Figure 26.1.	Client Server Network Connection	199
-	Service Naming	
-	Net Service Name Window	
Figure 26.4.	Protocol Window	200
Figure 26.5.	Host Name Window	200
Figure 26.6.	Service Name Window	201
Figure 26.7.	Test Window	201
Figure 26.8.	Result Window	201

#### 27 SampleTrack<sup>™</sup> System and Client Installation 203

Figure 27.1.	The SampleTrack Installation CD	203
Figure 27.2.	The SampleTrack Installation Wizard	204
	Customer Information Window	
	Setup Type Window	
Figure 27.5.	Installation Overview Window	205
Figure 27.6.	InstallShield Wizard Completion Window	206
	Installing the SampleTrack Patch	
	Information Window	
Figure 27.9.	Database Export Window	207
Figure 27.10	.Standard Oracle® Host String	208
	.SQL Script Window	
	.Confirmation Window	
Figure 27.13	.Command Prompt Window	209
	.Load Components Window	
Figure 27.15	.Component Selection Window	210
	.Output of SQL Statements	
	.Program Completion Window	
	.Change Completion Verification Window	
Figure 27.19	.The Login Window	212
Figure 27.20	.The SampleTrack Client Window	212
Figure 27.21	.SampleTrack Program Icons	213
•	.The Toolbar	
	.The ST Config Window	
Figure 27.24	.The SampleTrack Administration Menu Option	214
Figure 27.25	.BDE Administrator Window	215
Figure 27.26	.BDE Administrator Window: DLL32 and VENDOR INIT	216
Figure 27.27	.Environment Setup Window	217
	.The Load SampleTrack Components Window	
	.Spectrometer Configuration Window	
	.SQL Output Window	
	.The Instrument Setup Window	
	.The New Instrument Name Window	
Figure 27.33	.The Import and Export Scripts Window	220
Figure 27.34	.The Explorer Window	221

225

Figure 27.35.InstallShield Wizard Welcome Window .	
Figure 27.36.InstallShield Wizard Window	
Figure 27.37. The Destination Folder Window	
Figure 27.38. The Setup Type Window	
Figure 27.39.The Install Window	
Figure 27.40.InstallShield Wizard Completed Window	224

#### 28 SampleTrack<sup>™</sup> Web Client Installation

Figure 28.1.	InstallShield Wizard Custom Setup	225
Figure 28.2.	SampleTrack Plug In Overview	226
Figure 28.3.	Group Tree Window	226
Figure 28.4.	Login Configurations	228
Figure 28.5.	Navigation Elements Window	228
Figure 28.6.	SampleTrack Web Configuration Window	229
Figure 28.7.	Edit Rights Window	231

#### Figures

A

activate the trigger 139 add a complete new order 27 add a Preparation 26 add an Experiment 26 Add D2O box 49 Add experiment 48 Add Experiment button 26 add new groups 97 add or delete existing filters 109 Add Preparation button 26, 47 Add Proton Exchange 26, 49 Add Sample Order button. 45 add table default values 106 Add User button 103 adding a new group 97 Adding an Experiment to a Finished Sample Order 53 Adding D20 for Proton Exchange 49 Adding Table Defaults to the New Group 106 Admin 113 administration facilities of the Object Manager 99 Administrator 31 administrator 97, 101 Administrator Toolbar 130 administrator-configurable option 66 AMIX 58 Application management 61 approve all finished orders 42 approve finished orders 42 Approved status 42 Archiving of all associated data 17 Assembling a Group of Experiments Under One Title 95 assign an EasyDialog configuration to a user group 108 Attachments 140 Automatic Entering of the Tube ID 111 automized workaround 34 Autostart folder 203 В **B-ACS 111** bar code reader 35 bar-code already in use 111 bar-code collar 111 barcode collar 36 base tables 117 **Basic filters 30** basic filters 31 **B-BCE 36** B-BCE for order approving 42

begin the measurement 27 **BEST 46** Body 141 **BSTS REGISTRY 83** bstsdata 143 buttons 27 С C13 experiments to be run over the night 38 Case Sensitive box 55 Charset 140 client access privileges 102 client ID 45 client needs 98 Client PC Operation System Requirements 178 Client Profile 101, 102 Client Profile from the Object Manager main menu 105 client profiles 101 collar 111 **Communication Overview 18** Communication Paths & Files 89 Communication Server 137 communication server 82, 83, 111 Companies setup module 85 Config 113 Configure Plugins window 80 configure solvent mixing methods 93 Configuring a Status 129 confirmation of a data 67 Content sub-windows 100 Content window 106 Content window. 101 **Contract Button 28** contracted 28 Control key 49 control the flow of samples to the various laboratory automation 102 Copy an Existing Experiment 49 copy an order 27 Copy Sample Order Button 27 create a new group 109 create a new table default value 106 Create a User Group 105 Create New Sample Order 27 Create New Sample Order Button 27 Creating a New Form Adding Buttons 133 Adding the First Object 132 Adding the Other Group Boxes 133 Adjusting the Group Box 132

Establishing Group Boxes 131 Opening the Form 131 The Completed Group Box 132 Creating a New Group 105 Creating a New Report Within the Sample Order Details Window 60 Creating a New UNC Formatted Link to a Document File 60 customer notification via E-Mail 43 **Customize Filter Settings 109** D D2O 49 Daily Routine Laboratory Functions 17 daily routine laboratory functions 17 Data Exchange and Directory Structure 143 database 111 Database Connection Error After New Installation 215 **Database Installation 179** Date\_End 118 Date\_Entry 118 Date\_Start 118 day-to-day routine operations 15 default user groups 97 default value 128 Default Value in the Fields Overview Window 128 default values 108 Define a Method and Enter the Journal Number 46 Define Filter box 31, 32 **Delay Experiments 38** Delay Experiments window 38 Delayed experiments 51 delayed experiments 39 Delete Group 99 Delete Sample Orders Button 27 delete the solvent from the list of solvents 90 Deleting a User, Department, or Company 89 Details Preparation 24 details experiments 25 Details of a Preparation 24 details of an order. 27 Details Window Buttons 26 Difference between the SampleTrack System and Client Installation 203 Directory Sharing with UNIX Instrument Computers 178 **Document Files 59** Doubleclick the experiment 57 Drag & Drop Objects 125 drag & drop the configuration 108

E

EasyDialog 65 EasyDialog configuration 123 EasyDialog configuration for a user group 108 EasyDialog configurations 100 EasyDialog form 125 EasyDialog Login Window 65 EasyDialog Parameters 113, 135 Edit - Insert Table Defaults from the Object Manager main menu 106 edit a workstep 27 Edit Item Button 27 Edit Table Defaults window 106 Editor 59 **Effective Rights 99** E-Mail 43 E-mail Address Registering 142 E-mail addresses 137 E-mail Interface 139 E-mail Interface Trigger Table 141 E-mail Internal Processes 138 E-mail notification 137 E-mail Parameters 140 E-Mail Support 137 E-mail Trigger Setup 139 Enable Administration box 31, 32 Enter Sample ID 45 Entering data in the EasyDialog Window 66 Entering the Sample and Client ID 45 Entering the Tube ID Manually 111 Environment Setup 43, 85 Example Creating a New Input Form 130 **Example Client Profile 101** Expand / Contract Button 28 expanded 28 Experiment copy 49 Details 25 experiment button 26 **Experiment Details 25** experiment group 95 **Experiment Substitution for Imported Files 95** experiment symbol 58 **Experiment Translation Table 95 Experiment Viewer 62** Experiments waiting for HD Exchange 40 F Field / Value Evaluation 168

Fields Overview Required Values 128 Fields Overview window 127 file imports and exports 82 Filter & view assistant button 30 filter & view assistant. 30 Filter and Search Tools 54 Filter and view management 30 Filter Definition 54 filter definition window 54 Filter Information 99 filter information 100 Filter management 32 Filter Tools 54 Filters & Views 30 Finish Manual Experiments 37 Finished Sample Order 53 Footer 141 Function Archiving 171 Send Emails 169 Functions for File Import/Export 162 Functions for Script Processing 161 Functions for Tables 164 G General Rights 106 **GLP 19** Gobal Database Name 182 Grid Distance 126 grid lines 126 group client profiles 97 Group Details in the Content Window 101 Group ID 45 group management tool 103 group of experiments 95 **Group Properties 99** group sub-category 101 Group Tree 100, 101 Group Tree and Content Sub-windows 100 Η HD Exchange 40 HD Exchange Solvent window 40 Header 140 hits 83 homepage 177 Host 140 host name 200 How to Drag & Drop Objects 126 Hydrogen/Deuterium 40

Ι IconNMR 83 ICON-NMR<sup>TM</sup> 3.1 or higher 178 Icons 213 import a library 81, 82 In and Out directory 145 In directory 143 In or Out directory 144 individual laboratories requirements 105 individual laboratory 97 individual laboratory requirements 67 individual order generation 51 Insert Client Profile 99 **Insert Group 99 Insert Table Defaults 99** Inserting a New Group Name 109 Inserting a New Preparation Method 92 Inserting a New User in the Companies Setup Module 85 Inserting a Proton Exchange Solvent 91 Inserting New Experiments 94 Installation of E-mail support 138 Installation of the SampleTrack System 204 Installation steps 138 Instrument 143 Instrument\_Name\_Key 119 instruments 85 instruments and systems like spectrometers, preparation robots or LIMS 145 Interaction with AMIX<sup>TM</sup> 58 Interaction with Structure Editors 59 Interaction with TOPSPIN<sup>TM</sup> 57 INTERF\_TRIG\_TAB 138 interface 15 interface for BRUKER spectrometers 14 Introduction to Data Exchange 143 ISIS/Draw<sup>TM</sup> 59 **ISISD**raw 62 ISO 9000 13 Items contracted 28 Items expanded 28 J JCAMP Definition File 145 journal number 46 Journal\_Number 119 L Laboratory Information and Management System 14 laboratory journal number 59 Ligand list 83 LIMS 13, 143

LIMS Import workaround 34 LIMS imported 51 Local/Service Naming 199 login 21 login date and time 21 Lookup Functions Database 167 Μ Mail ID 140 main window 23 Manage Users 99, 103 Manual Preparation 35 manual preparation 33 manual preparation area 35 Manual Submit Inside the Sample Order Details Window 51 mark all orders 34 mass spectrometers 13 matrix database 82 Matrix ID 82 Matrix organization 81 Matrix Settings 80 Matrix Type 80 MDL® ISIS/Draw<sup>TM</sup> 59 Measurement 51 **MESTRE-C 62** method 46 Method\_Type\_Key 119 METHODS 117 Microsoft Windows® 59 Microsoft Word® document 59 Microsoft® Word 60 modification of an instrument 89 modifying an existing group, 97 Module Pooling 82 More button 59 MS/SPR/NMR/XRAY 20 MultiInstance 113 Ν netservice name "dbstrack" 200 network protocol TCP/IP 200 new Client Profile block 105 New Department 87 new experiment 94 New Filter Method 109 new group 105, 107 New Group EasyDialog Configuration 108 New Instrument 89 New Method Type 89

new preparation method 92 New Solvent 90 New User 88 new users 85 night or weekend run over the night or weekend 38 **NMR 46** NMR laboratory 14, 15 NMR\_EXPERIMENT\_WS 117 NMR PREPARATION WS 117 NMR\_Preparation\_WS\_Key 119 0 **Object Manager 31** object manager 98 Object Manager main menu 103, 109 object-oriented, drag, drop & modify methodology 123 OK button 49, 53 only minimal rights 31 Open a form / Create a new form 123 Open EasyDialog from the SampleTrack<sup>™</sup> Client Tool 65 Open the Object Manager 98 **Opening EasyDialog 65** Opening the EasyDialog Configuration Window 123 Oracle 179 Oracle 8i Service Name 201 ORACLE® Client 8.1.7 178 Oracle® database 111 Oracle® Installation 179 **ORACLE®** Standard Edition 8.1.7 177 OraHome92 180 Order Approval 42 Order Approval window 37, 38 orders imported 34 Out directory 143 overview window 23 Ρ paperclip icon 57, 58 Password 113 Paste 99 path 61 path in the environment system 185 PC requirements 177 personal data pool 30 Plot Title 48 Plot\_Title 119 Plugins window 80 Pooling Module 79 Pooling Orders Window 82

Pooling Settings 81 Pooling Size 81 Pooling Volumes 81 Popup Rights 106 Port 140 port number 181 **PREP 143** Preparation 35 preparation and measurement 51 preparation button 26 Preparation details 24 Preparation is Valid 92 Preparation is Visible 92 Preparation Setup window 49 preparation system 79 Preparation Workstep 47 preparation workstep 24 Preparation\_Name\_Key 119 privileges 97 Project No. 45 Project No. default value 108 Project\_Number 118 Property Editor 126 Protein Imaging 84 **Protein Screening 82** protein screening 82 protein-ligand binding detection 79 Proton Exchange 37 Proton Exchange button 26 proton exchange buttons 26 proton exchange solvent 37 Proton Exchange Solvents 91 Q **QA/QC 13** quick overview of the rights 99 R release the Tube ID 37, 38 Reload or Refresh Button 27 Remove Users button 103 Report Editor 62 Report Rights 106 report within the Sample Order Details window. 60 **Required Values 128** Rerun an experiment 52 rerun the complete order 52 Rerunning a Faulty or Finished Experiment 52 Reset complete Sample Order 52 **Reset Sample Order Experiments 52** 

Resetting a distinguished experiment or a preparation 52 restart an experiment 27 result data path 58 reuse Tube ID bar-code collars 42 right edge of the sample order input window 26 rights and privileges 97 rights of the entire group 97 RIMS 20 routine laboratory functions 17 routine sample order 65 run over the night or weekend 38 S SAMBA daemon 178 Sample ID 24, 45 Sample Order Details window 45 Sample Order Experiment Details 25 Sample Order Status Information 29 Sample Order Window Buttons 27 Sample Orders module 23 Sample Title and/or Plot Title 48 Sample Title and/or Plot Title is automatically created 48 Sample Viewer 62 Sample\_Amount 119 Sample\_Changer\_Holder 119 Sample\_Charge 118 Sample\_Fraction 118 Sample\_ID 118 Sample\_Nature 118 SAMPLE\_ORDER Table 118 Sample\_Order\_Number\_Key 118, 119 SAMPLE\_ORDERS 117 Sample\_Title 118 SampleRail<sup>™</sup> 79 SampleTrack bstsdata Directory Structure 143 SampleTrack Client Installation 221 SampleTrack Client Parameters 113 SampleTrack Communicator 203 SampleTrack Environment 43 SampleTrack Environment Setup 85 SampleTrack Patches 206 SampleTrack Program Icons 213 SampleTrack<sup>TM</sup> as Data Reference 16 SampleTrack<sup>TM</sup> communication 18 SampleTrack<sup>TM</sup> Login 21 SampleTrack<sup>TM</sup> Overview 22 SampleTrack<sup>TM</sup> Workflow Architecture 19 save the spectrometer time 38 Saving an EasyDialog Configuration 134

scenario 33 order 33 Script Library 159 Script Processing 161 Scripts for Automation 159 search function 55 search parameter 55 Search Pattern field 55 search through the current list of data 55 search through the Oracle® database 55 Search Tools 55 Select the Preparation Workstep 47 selecting/deselecting the various rights 101 Server Requirements 177 setup parameters 85 shortcuts 115 Show Environment 99 Show Messages 168 software requirements 177 solvent 35 solvent for the proton Hydrogen/Deuterium exchange 40 Solvent is Valid 90 Solvent is Visible 90 Solvent Mixing Method 93 Solvent\_Amount 119 Solvent\_Name\_Key 119 special filters 54 specific filter 105 **SPECT 143** spectra data as attachment 43 Spectrometer Settings 89 Spectrometer Software Requirements 178 standard database fields 125 standard policies of the individual laboratory 97 Start delayed experiments button 39 Starting a Preparation and Measurement 51 Status 118 status codes and their meaning 129 status dependencies 33 status information 29 status symbols 29 stinstXXX\_XX.exe 204 Structure 59 Structure Editor 62 Structure Editor program 59 structure elucidation 57 stscript.bpl 159 Subject 140

Submit button 51 submit button 51 Submit Sample Order Button 27 Submit Sample Order button 51 Submit Sample Orders 34 submit the sample orders 51 Submit/Verify 51 submitted/verified status 35 switch to the next experiment 57 SysButtons 113 System Administrator 85 system passwords 184 System Requirements 138, 177 Т Table Defaults 100 Table Defaults SAMPLE\_ORDERS 107 table structures 143 TECAN® - Pooling Module 79 **TECAN®** Generating Samples 82 **TECAN®** Instrument Setup 82 **TECAN®** Matrix Settings 80 **TECAN®** Plugin Installation 80 **TECAN®** Requirements 80 template file 59, 60 The Administrator Toolbar 130 The EasyDialog Window 66 The environment setup area 43 The main menu option View 99 The METHODS Table 119 The NMR\_PREPARATION\_WS Table 119 The Object Manager 97 The right to define Filters 31 The SampleTrack<sup>TM</sup> Interfaces 15 The STATUS Field 129 The submit window 51 The Tube ID and TubeID.txt file 111 **TOPSPIN® 57** TOPSPIN® 1.2 or higher 178 Tree overview 22 Tube ID 36, 111 tube id 67 Tube ID and TubeID.txt file 111 Tube ID number 111 Tube ID release 37 TubeID.txt 111 TubeID.txt file 111 tubes waiting for exchange window 41 Tubes Waiting for Exchange window. 41

U **UNC 59 Universal Naming Convention 59 UNIX 178** Update View 99 Updating the jobs in the Tube ID File 111 **User** 140 user classification 101 User Classification and Client Profiles 102 user group 97 user level 110 user level privileges 102 user level. 102 user specific filters 30, 54 User\_Name\_Key 119 Username 113 users 103 V Variables 99 Variables and Functions 159 variety of experiments under one experiment title 95 VCU Preparation Table 121 VCU Request tables 120 VCU Sample Order Table 120 VCU\_REQUEST\_ALL 120, 143 verification of measurement results 57 Verify the order 34 View all spectra Button 28 View Experiment Button 28 Views 30 Visible and Valid checkbox 92 W waiting / delayed experiments 38 Waiting experiments 39 WEB Client 69, 231 Web Spectra Viewer 72 Web structure viewer 72 wellplate design 80 wellplate for pooling 82 wellplates 79 window buttons 27 Windows® Explorer 59 workflow 33 Workflow example 145 workgroups 102 workstep 25 Worksteps block 47 worksteps for automation 79

X x-ray crystallography 20 XWIN-NMR<sup>TM</sup> 3.1 patch level 11 or higher 178

### Notes: