BRUKER BIOSPIN AG

Low Loss Cryostats Superconducting Magnets phone: ++41 44 825 91 11 fax: ++41 44 825 93 60 e-mail: magnetics@bruker.ch

Bruker



Handle the N2-Levelsensor with extreme care!



- Remove the N2-levelsensor carefully from the cardboard box.
- Try to hold the sensor in an upright position when carrying it around.
- Bend the hinge only when necessary
- Do not apply force to the tubes of the N2-levelsensor
- The hinge is bendable just in one direction. Bending it in the wrong direction destroys the hinge.
- Introduce the sensor carefully in the nitrogen turret. If it hits any resistance do not apply force but rather try a different turret.

Handle the N2-Levelsensor with extreme care!

think forward

N2-Levelsensor