


NMR Thermometer

- Variable Temperature Control Using the 2H Lock System of AVANCE III HD Spectrometers

User Manual
Version 002



Copyright © by Bruker Corporation

All rights reserved. No part of this publication may be reproduced, stored in a retrieval system, or transmitted, in any form, or by any means without the prior consent of the publisher. Product names used are trademarks or registered trademarks of their respective holders.

This manual was written by
Application & Electronics Departments
© July 11, 2014 Bruker Corporation

P/N: Z33085

DWG: Z4D12031

For further technical assistance for this product, please do not hesitate to contact your nearest BRUKER dealer or contact us directly at:

Bruker Corporation
Industriestrasse 26
8117 Fällanden
Switzerland
Phone: + 41 44 825 91 11
FAX:+ 41 44 825 9696
E-mail: nmr-support@bruker.de
Internet: www.bruker.com

Contents

1	Introduction	5
2	Principle of NMR Thermometer	7
3	Requirements	9
4	Getting Started	11
4.1	The edlock Menu.....	11
4.2	Setting up the NMR System for the NMR Thermometer.....	13
4.3	Examples.....	15
4.3.1	Monitoring Mode.....	16
4.3.2	Regulation Mode.....	17
5	Advanced Operation	19
5.1	Define a New Solvent.....	19
5.2	Selection of NMR Thermometer Compounds.....	21
5.2.1	Predefined Solvents for the NMR Thermometer.....	22
5.2.2	Technical Considerations.....	23
5.2.3	Considerations for Shimming with Topshim.....	24
6	Applications	27
6.1	Identical Chemical Shifts at Different Spectrometers.....	27
6.2	Identical Chemical Shifts for Experiments with Different Heating due to Experimental Conditions (Temperature Compensation).....	28
6.2.1	RF Heating.....	28
6.2.2	Heating caused by Spinning Speed (HR-MAS).....	29
7	Frequently Asked Questions (FAQ)	31
8	Contact	33
	List of Figures	35
	Index	37

1 Introduction

This manual is planned as a user manual with limited technical detail. The main focus is on usability, whereas a few easy examples are provided showing how to get started with this new tool. Through use of this manual the user should get an idea for what the NMR Thermometer™ might be used for.

The comparability of data (chemical shifts) and results derived from NMR data (diffusion data, relaxation measurements), as well as the quality of the NMR spectra, depend on an accurate and precise temperature measurement. That is the reason why the temperature measurement should ideally take place inside the NMR tube and not outside of it. In general the temperature sensor of the probe is not reflecting the real situation inside the NMR tube.

The NMR Thermometer not only monitors the temperature, but also compensates for different heating effects (e.g. RF heating) that occur during an NMR experiment.

2 Principle of NMR Thermometer

The NMR Thermometer measures the temperature inside the sample by observing the chemical shifts of two ^2H signals using the lock channel (LTRX board) of the NMR system. The distance between the two signals is transferred into a temperature and directly used by the temperature control unit (BVT) for regulation. Thus, the NMR Thermometer acts as a temperature sensor (see figure below).

To obtain the second lock signal, a thermometer substance needs to be added. One of the signals should also be temperature dependent.

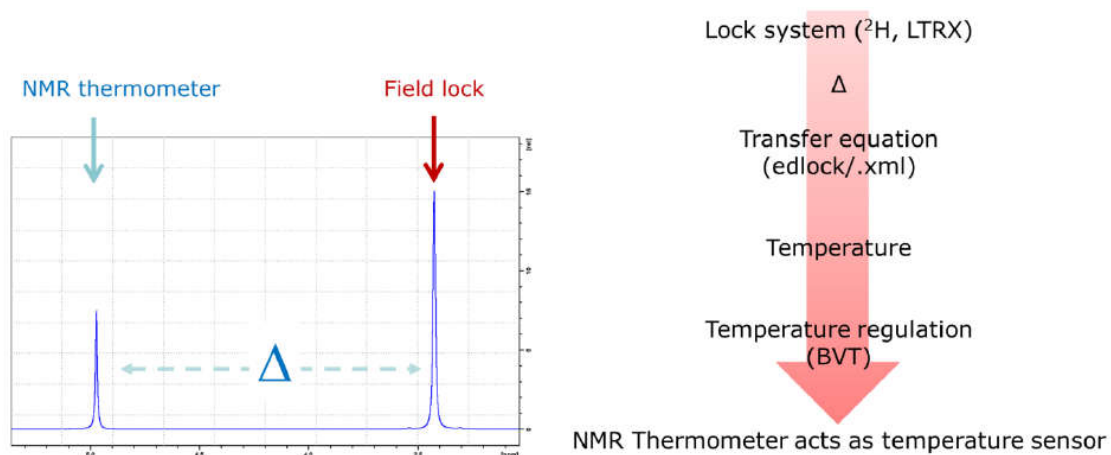


Figure 2.1: Principle of the NMR Thermometer.

3 Requirements

Hardware: Avance III HD. A hardware upgrade is required for Avance III, whereas an update to SmartVT™ and Digilock 2G is required.

Software: TopSpin version 3.2 or higher.

Firmware: Versions for Avance III HD.

4 Getting Started

This chapter provides a short and straightforward introduction on how the NMR Thermometer works and guides you through the first steps using the NMR Thermometer. As an overview, the steps include:

- Preparation of the edlock table if not yet done.
- Locking on the corresponding solvent used for the NMR Thermometer.
- Optimizing the shim and lock phase.
- Activating the NMR Thermometer in the variable temperature control.
- Performing self-tuning of the variable temperature unit. This will optimize the regulation parameters for both the VT control with the sensor and the NMR Thermometer.
- Calibration of the spectrometer for measurements of real samples after the setup of the NMR Thermometer is finished.

4.1 The edlock Menu

First we will provide a short introduction to the software implementation (*edlock*, *edte*).

Start *edlock* by typing **edlock** on the TopSpin command line. The *edlock* window opens up.

The new *edlock* (starting for TopSpin 3.0 and higher) contains both the solvent list (formerly *edsolv* in the Edit Solvent Parameters figure below) and the lock parameters for every solvent (see the The Submenu Lock figure below).

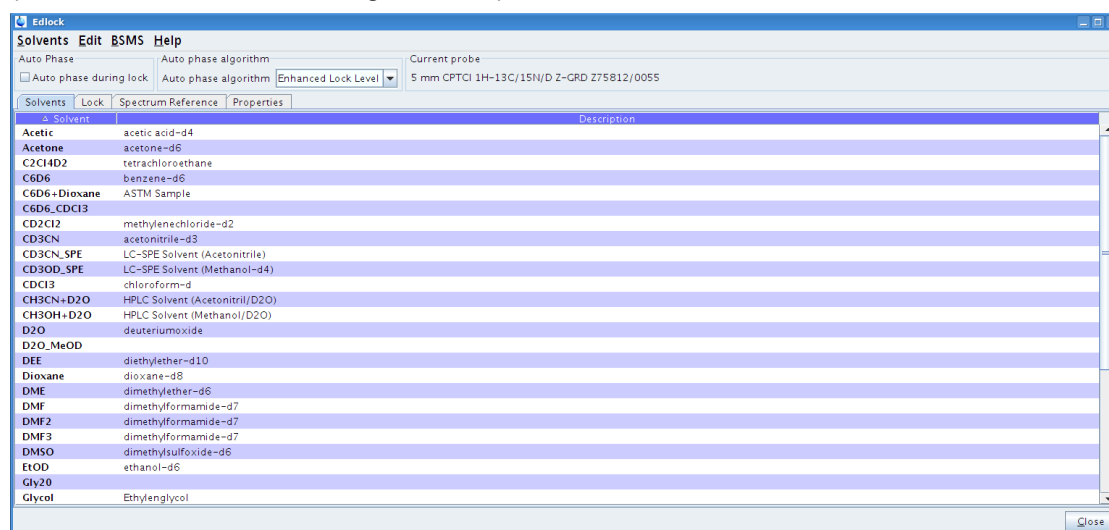


Figure 4.1: The Submenu Bar: Solvent.

The Edit Solvent Parameters window allows you to enter several parameters for the solvent, the melting point setting, and the boiling point of the solvent:

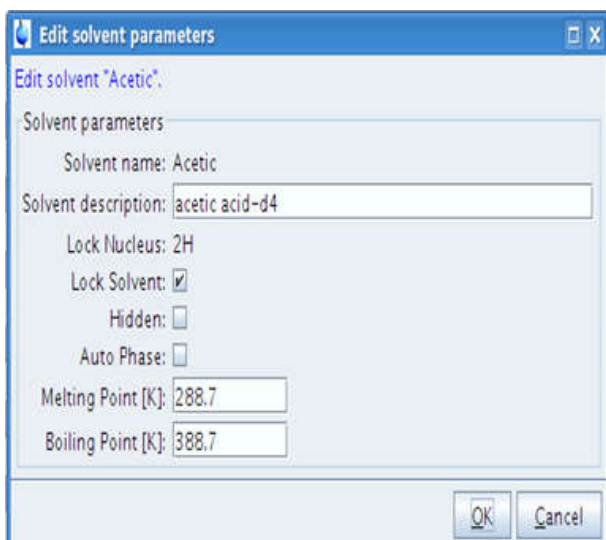


Figure 4.2: Edit Solvent Parameters.

In the submenu Lock window the solvents are listed, along with the probe for which the definition is valid (generic or specific probe), as well as other lock parameters (lock power, lock regulation triplet etc.).

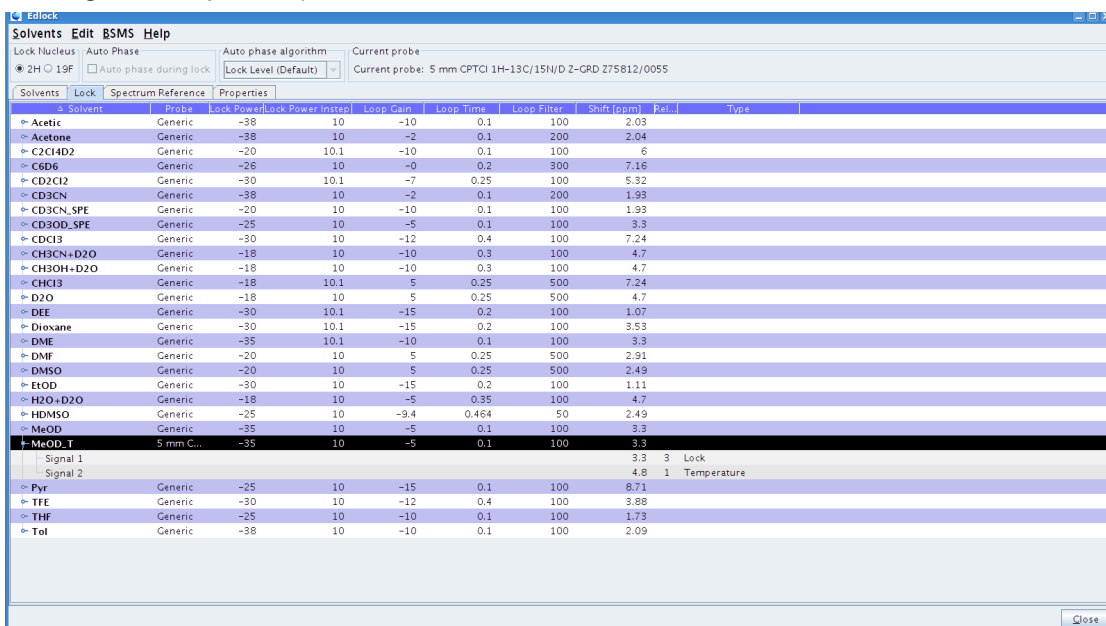


Figure 4.3: The Submenu Lock.

By executing a right mouse click on a solvent entry, a pull-down menu opens with the option for editing the lock parameters:

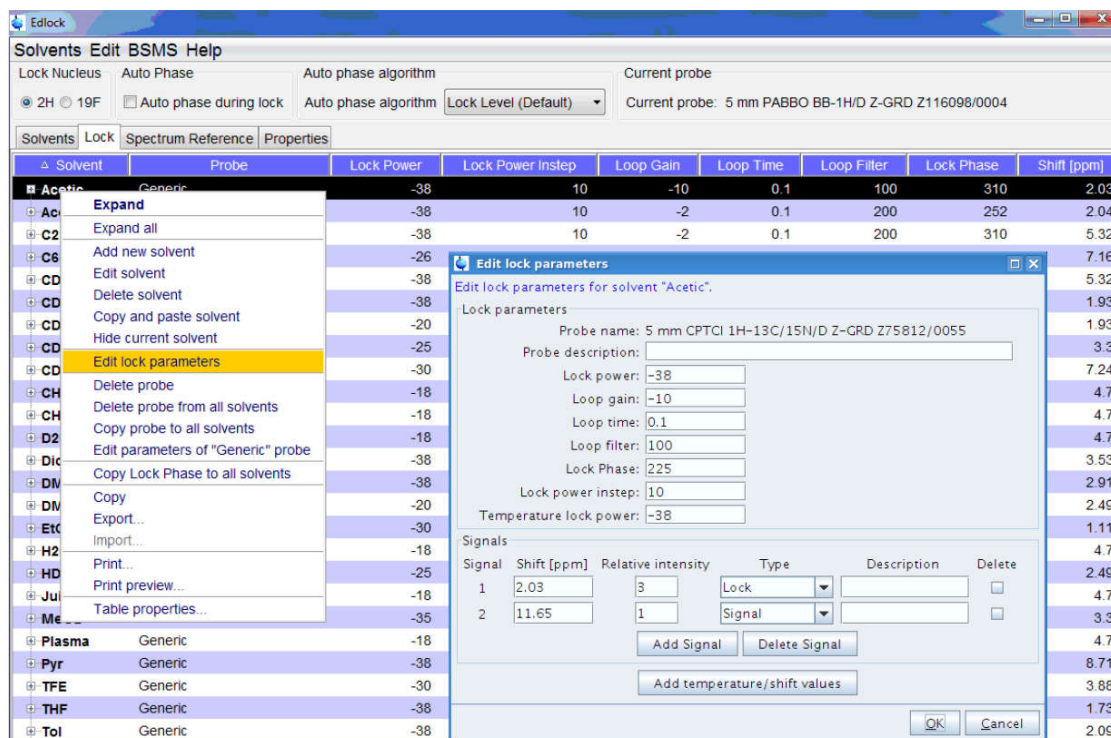


Figure 4.4: Edit Lock Parameters.

Every solvent entry contains information about every ^2H signal for that compound.

It is possible to define all the signals for solvents with more than one ^2H signal either as Signal, as Lock (signal used for field lock) or as Temperature (signal used for the NMR Thermometer).

For any NMR Thermometer substance, temperature and shift values can be added or imported (see below). The NMR Thermometer has its own lock power (figure above). For methanol (NMR Thermometer, standard sample: 99.8% deuterated methanol) a default solvent with corresponding lock parameters and temperature-shift value has already been defined.

4.2 Setting up the NMR System for the NMR Thermometer

Since the NMR Thermometer is observing a ^2H signal, the system has to be properly set up (lock, shimming, optimal lock parameters for the field lock and NMR Thermometer), similar to any other NMR measurement. This means that you should perform an automatic tuning and matching (*atma*), lock-in, and shimming. The lock-in procedure works as usual (type **lock** on the command line and select the solvent). If you lock directly on a solvent dedicated for the NMR Thermometer (methanol or another solvent), a temperature value is immediately shown, either in the monitoring mode or regulation mode (figure below). After that you can perform *topshim*. A decent line shape (no unusual line splitting) is mandatory for an exact temperature measurement inside the sample.

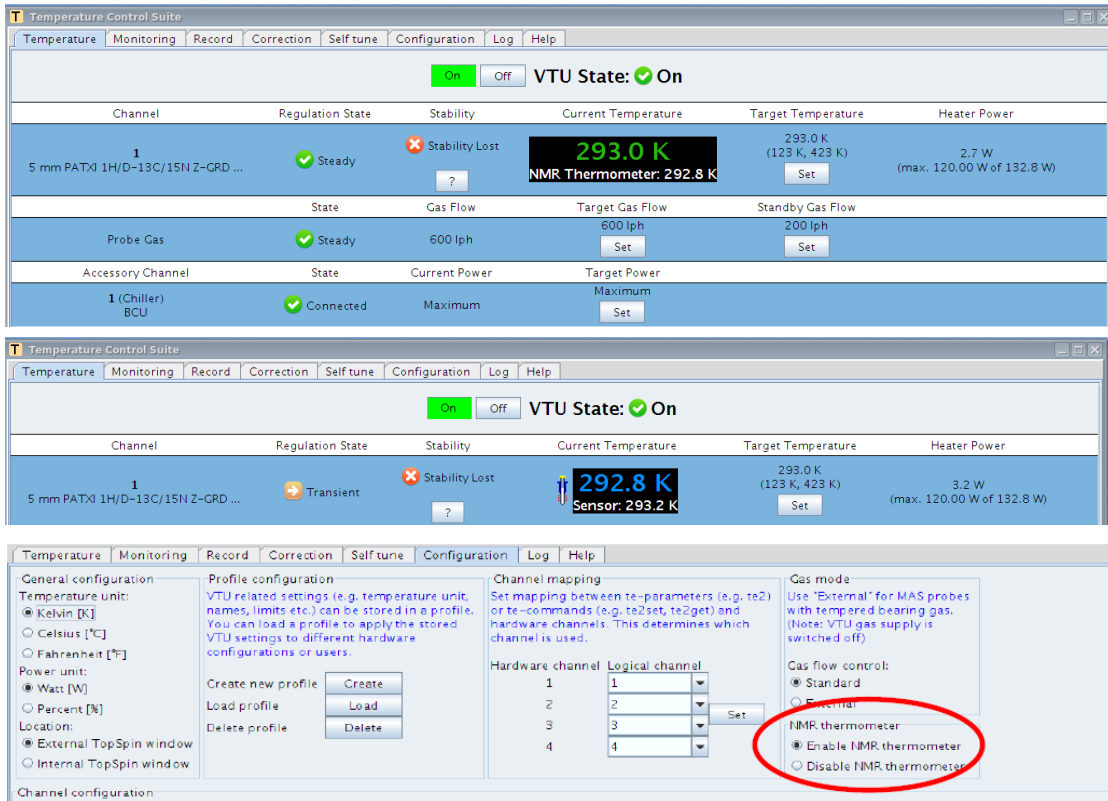


Figure 4.5: Edte Window: NMR Thermometer monitoring mode (disabled, top), regulation mode (enabled, middle) and the selection of both modes in the Configuration menu of the edte window (bottom).

Another important parameter is the *lockphase*, which can be optimized automatically by starting *autophase* (BSMS display). The procedure used for *autophase* is selected in the edlock window (Lock Level Default, Spectrum, and Enhanced Lock Level in the figure below).

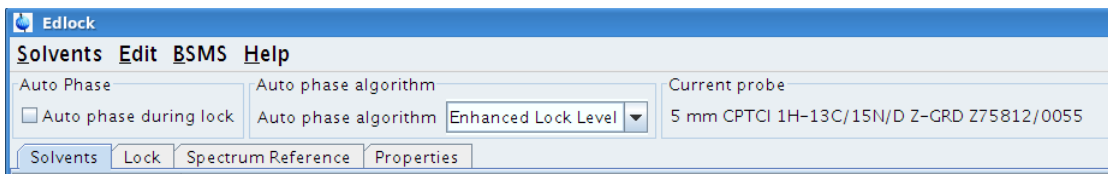


Figure 4.6: Selection of the Auto Phase Algorithm.

To prevent saturation, the *Lockpower* and *lockgain* for both lock channels should be optimized as well.

Since the NMR Thermometer contains both the lock and the temperature regulation components, the PID values for the temperature regulation need to be adjusted for each by using *selftune*. You will be notified by the system if a *selftune* is recommended:

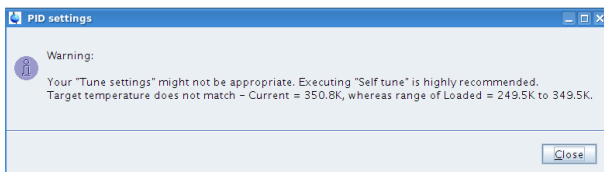


Figure 4.7: Selftune warning about PID parameters misfit.

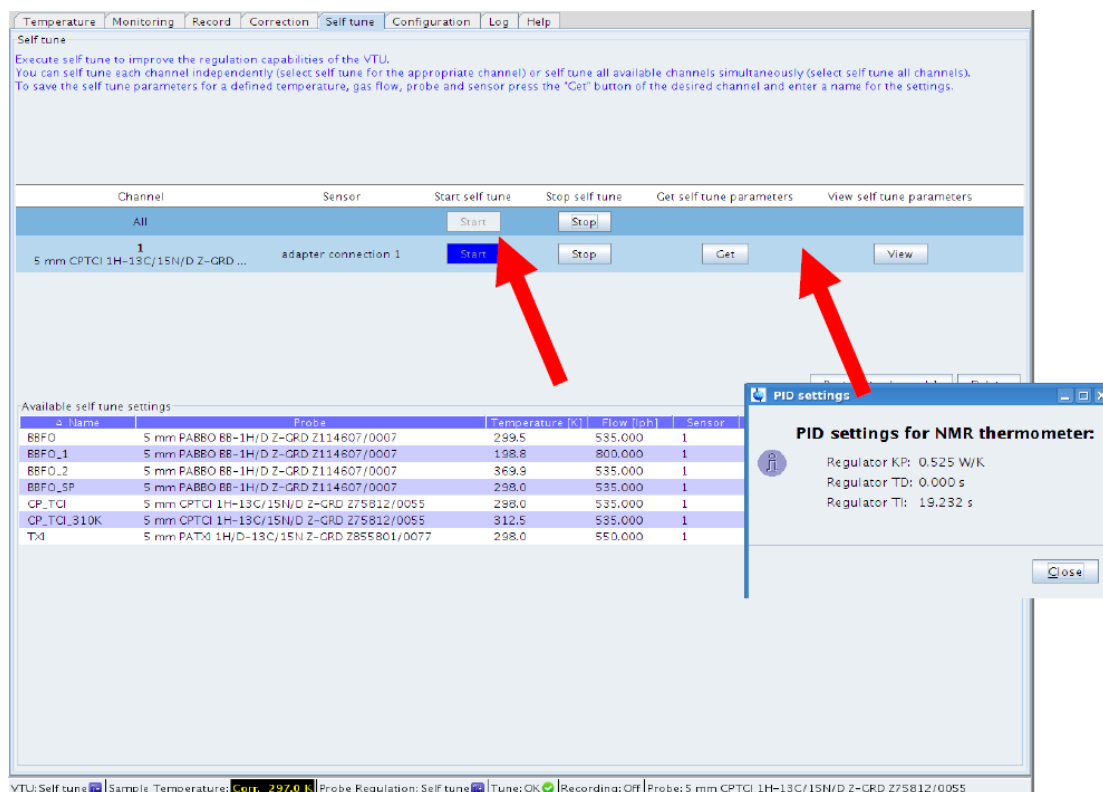


Figure 4.8: Starting the selftune procedure from the edte window.

The *selftune* should be carried out on both temperature channels (All).

4.3 Examples

An easy example to begin with is to use the methanol sample (standard sample: 99.8% deuterated).

Since this sample is the reference for the NMR Thermometer, the solvent entry in *edlock* containing the temperature and shift values is already predefined.

As mentioned, the NMR Thermometer is running in two different modes: monitoring and regulation mode (see above).

Assume that the lock parameters are already optimized.

4.3.1 Monitoring Mode

The following steps need to be performed:

- Insert sample.
- Tune and match by using the *atma* command.
- Lock-in – temperature appears.
- Topshim.
- Selftune.



Figure 4.9: Selftune in Progress.

- *Edte* – disable NMR Thermometer (if not already done).
- Change target temperature (*edte* for instance, a 10K temperature jump).

As an exercise we will perform a temperature jump (e.g. 10K) and follow the different temperature and other BSMS values (flow/heater) in the monitoring window (*edte* - Monitoring, as well as, activate NMR Sensor Temperature, NMR Thermometer, Target Temperature, Current Power etc.).



Figure 4.10: Monitoring several BSMS values during a temperature jump.

4.3.2 Regulation Mode

The following steps need to be performed:

- Insert sample.
- Tune and match – *atma*.
- Lock-in – temperature value appears.
- Topshim.
- Selftune (maybe already done in example 1).
- Edte – enable NMR Thermometer.
- Perform an experiment with internal RF heating (e.g. TOCSY).

After setting up the system (lock, shim) and enabling the NMR Thermometer (*edte*) the sample temperature is used for temperature regulation, which is visible on the status bar:

Sample 	Shim Coil Temperature 306 K	POWCHK 	Sample Temperature 298.2 K On Reg. State:	Temperature Channel 1 300.0 K	Spooler queued: 0 delayed: 0 cron: 0	BSMS status message Δ Z 2 Autoshim Locked Error	Time 16:53:21 May 04
------------	--------------------------------	------------	--	----------------------------------	---	--	----------------------------

Figure 4.11: TopSpin status bar with NMR Thermometer enabled.

To test the performance of the NMR Thermometer start a TOCSY experiment and follow the temperature values (sensor, NMR Thermometer) in the monitoring window of the VTU display (*edte*). One can nicely see that the temperature inside the sample increases over a certain

period of time and that the system immediately reacts to that by reducing the heater power (figure below - lower part), and hence the sensor temperature (figure below - upper part, white line).



Figure 4.12: Monitoring of the sensor and the sample temperature in edte during a TOCSY experiment

In the figure above, the upper part shows the temperature of the sensor and the NMR Thermometers; the lower part shows the heater power.

5 Advanced Operation

5.1 Define a New Solvent

To work with your own NMR Thermometer substances, you first have to define a new solvent in the *edlock* table.

Select a solvent which is similar to your mixture (similar lock parameters) and click the right mouse button. In the resulting pull-down menu you can add the new solvent:

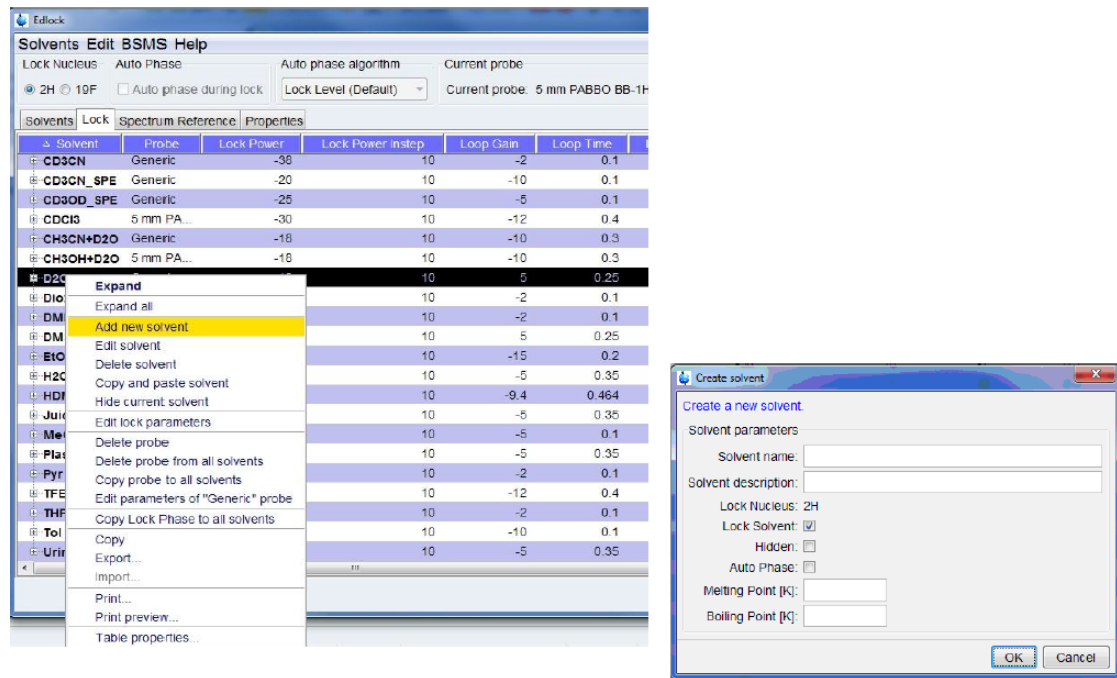


Figure 5.1: Adding New Solvents.

In the edit lock parameters window the signals can be defined (Shift) and assigned (Type) as Lock, Signal or Temperature:

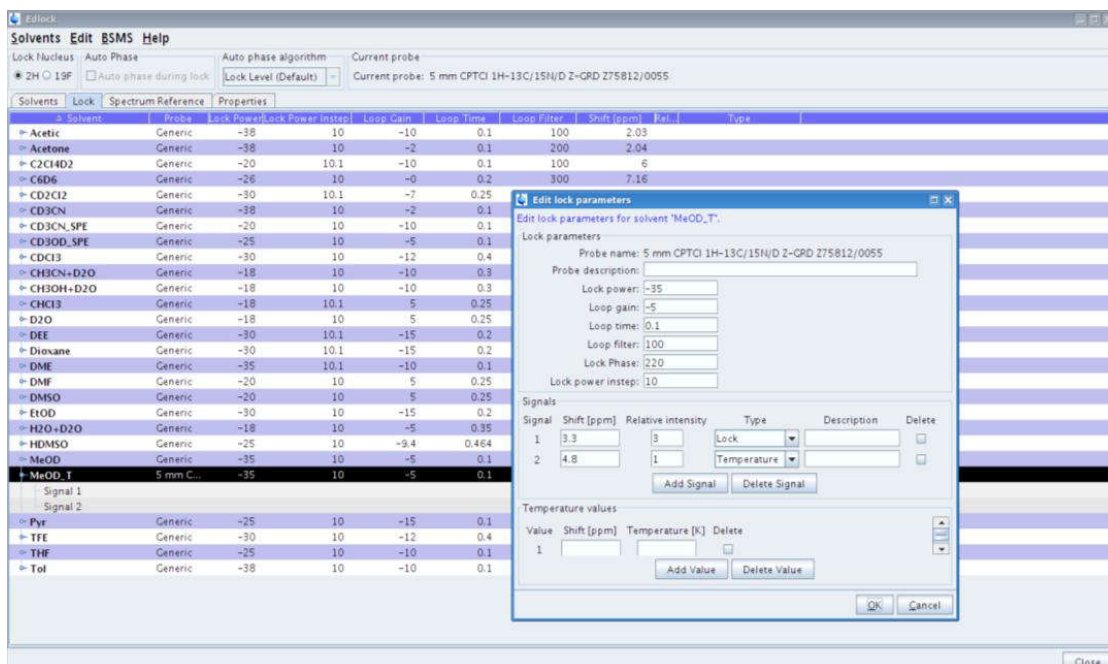


Figure 5.2: Assign Signals to Type of Signal (Signal, Lock, Temperature).

In the next step you should import the shift-temperature values (.csv or .xml format) or fill in the values manually. You can also create your own .xml (see below).

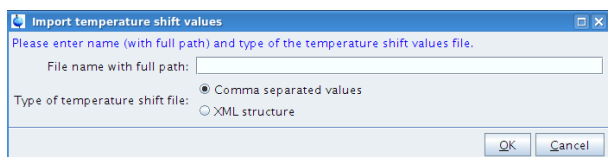


Figure 5.3: Dialog for importing temperature shift value files.

The .xml file (figure below) contains the name of the solvent (identical with the solvent name) and two shift values, one for the field shift and one for the temperature shift. The values have to be identical with the values defined for the particular solvent in *edlock*.

```
<table substance="Naac50">
  <field shift="4.7" />
  <temperature shift="1.8" />
  <point delta-shift="2.4766" temperature="346.1600" />
  <point delta-shift="2.5198" temperature="341.3700" />
  <point delta-shift="2.5634" temperature="336.6200" />
  <point delta-shift="2.6080" temperature="331.8000" />
  <point delta-shift="2.6532" temperature="327.0000" />
  <point delta-shift="2.6990" temperature="322.2400" />
  <point delta-shift="2.7458" temperature="317.4400" />
  <point delta-shift="2.7933" temperature="312.6200" />
  <point delta-shift="2.8422" temperature="307.8100" />
  <point delta-shift="2.8915" temperature="302.9800" />
  <point delta-shift="2.9422" temperature="298.1800" />
  <point delta-shift="2.9937" temperature="293.3600" />
  <point delta-shift="3.0464" temperature="288.5500" />
  <point delta-shift="3.1004" temperature="283.7200" />
  <point delta-shift="3.1566" temperature="278.9200" />
</table>
```

Figure 5.4: Example of a shift-temperature file in .xml format.

After the import the shift and temperature value are filled in the *edlock* table for the selected solvent and used afterwards for the lock-in process. Create a new XML file:

Edit lock parameters for solvent "MeOD_T".

Lock parameters

Probe name: 5 mm CPTCI 1H-13C/15N/D Z-GRD Z75812/0055

Probe description:

Lock power:

Loop gain:

Loop time:

Loop filter:

Lock Phase:

Lock power instep:

Temperature lock power:

Signals

Signal	Shift [ppm]	Relative intensity	Type	Description	Delete
1	<input type="text" value="3.3"/>	<input type="text" value="3"/>	Lock	<input type="text"/>	<input type="checkbox"/>
2	<input type="text" value="4.8"/>	<input type="text" value="1"/>	Temperature	<input type="text"/>	<input type="checkbox"/>

Temperature shift values

19	<input type="text" value="1.3449"/>	<input type="text" value="316.07"/>	<input type="checkbox"/>
20	<input type="text" value="1.3546"/>	<input type="text" value="315.1"/>	<input type="checkbox"/>
21	<input type="text" value="1.3644"/>	<input type="text" value="314.16"/>	<input type="checkbox"/>
22	<input type="text" value="1.3742"/>	<input type="text" value="313.19"/>	<input type="checkbox"/>
23	<input type="text" value="1.3839"/>	<input type="text" value="312.23"/>	<input type="checkbox"/>

Figure 5.5: Temperature shift values filled in.

For many NMR applications (for instance, Bio-NMR in aqueous solutions) an estimate of the slope (be aware of the fact that the slope can have negative or positive sign!) of the temperature dependency and the knowledge of the offset (absolute temperature correction) are sufficient (the necessary temperature range might be rather small) to create an .xml file. It is of course also possible to determine the correct slope automatically.

5.2 Selection of NMR Thermometer Compounds

The simplest compound for the NMR Thermometer is fully deuterated methanol used as a NMR solvent. While the deuterium signal of the methyl group is used for the field lock, the deuterium signal of the hydroxyl group is used for the NMR Thermometer. The deuterium chemical shifts of commonly used organic NMR solvents like CDCl₃, acetone-d₆ and DMSO-d₆ are virtually independent on the sample temperature. In that case, deuterated methanol could be placed in a capillary or mixed with the solvent.

The chemical shift of water strongly depends on the temperature. For samples in aqueous solutions, a deuterated organic small molecule can be added as a thermometer compound. Details will be described in the following section.

The compounds listed in the table below have been tested and can be considered.

Name	CAS number	Availability	Number of D	Chemical shift [ppm]	Minimum Concentration [CryoProbe TCI 5mm 600MHz]
Methanol-d4	811-98-3	yes	3	3.3 (CD3) 4.8 (OD)	
Acetic acid-d4	1186-52-3	yes	3	2.03	20-50mM
Sodium acetate-d3	39230-37-0	yes	3	1.8	20-50mM
DMSO-d6	2206-27-1	yes	6	2.49	20mM
Acetone-d6	666-52-4	yes	6	2.04	20mM
Acetonitril-d3	2206-26-0	yes	3	1.93	20-50mM
Tetramethylammoniumchlorid-d12	23789-03-9	yes	12	3.2	10mM
TRIS-d11	n.a.	yes	6	3.5	20mM
TSP-d14	n.a.	on request	9	0	n.a.
DSS-d16	n.a.	on request	9	0	n.a.
Pivalic acid	42983-07-3	yes	9	1	10-20mM

Figure 5.6: Suggested compounds for the NMR Thermometer and samples in aqueous solution (D2O 5 – 100%), their properties and estimated concentration. Fully deuterated DSS and TSP are currently not available but will be synthesized on Bruker's request at small amounts only for internal tests.

For the selection of a suitable compound the following points should be considered:

- Large number of deuterons: The larger the number of chemically equivalent deuterons, the lower the required concentration. Tetramethylammonium chloride-d₁₂ is in this respect, the preferred compound, whereas sodium acetate is less favorable.
- Possibly one additional ²H signal.
- Moderate or no salt effect: Salts and acids in aqueous solution will increase the conductivity of the sample, which might cause a loss of sensitivity and an increase of the pulse length.
- The compound should not interact with the sample or change essential structural properties of the sample.
- Acceptance. Difference compounds are commonly added to protein solutions. These can be buffers like TRIS or stabilizers like EDTA. These compounds can be used if deuterated and available at a sufficiently high concentration.
- Price and availability.

5.2.1 Predefined Solvents for the NMR Thermometer

The lock table contains solvents which are setup for usage with the NMR Thermometer. For these solvents a temperature calibration of the chemical shift difference of the two ²H signals has been performed and is included. These are the following solvents:

T_MeOD: methanol-d₄

T_H2O+D2O+NaAc: sodiumacetate-d₃ in 90% H₂O, 10% D₂O

T_H2O+D2O+Me4NCl: tetramethylammoniumchlorid-d₁₂ in 90%, H₂O, 10% D₂O

T_H2O+D2O+Pivalate: pivalic acid-d₉ sodium salt in 90%, H₂O, 10% D₂O

Depending on the probe, the lock power may need to be adjusted.

5.2.2 Technical Considerations

When adding a thermometer substance to the sample, there are a few technical points to consider.

Depending on your NMR system (room temperature or cryogenically cooled probe, spectrometer frequency etc.) you need to add a sufficient amount of the compound in order to obtain a decent lock signal. Another concern is the distance from the ^2H signal which is used for the field lock (figure below). In addition, the dynamic range, which is the intensity ratio of the ^2H signal used for frequency lock and the NMR Thermometer, plays an important role when a highly or fully deuterated solvent (e.g. D_2O) is used as a field lock solvent.

The requirements for obtaining reliable lock-in and lock regulation performance include:

- S/N (^2H) > about 200:1 (signal > ~ 3 ppm distant from the main signal)
- Intensity ratio < 500:1 (field lock signal: frequency lock signal; $\Delta \sim 3$ ppm)

The smaller the distance to the second signal, the higher the concentration of the compound used for the NMR Thermometer should be. As a rule of thumb: reducing the distance of the two lock signals by a factor of 2 increases the necessary concentration (decreases the possible intensity ratio of the two signals) of the NMR Thermometer compound by the same factor. Examples:

- The D of the NMR Thermometer signal to the field lock signals is about 3 ppm. The required signal to noise ratio of the NMR Thermometer compound has to be at least 200:1 (I-ratio < 500:1).
- The D of the NMR Thermometer signal to the field lock signals is about 1.5 ppm. The required signal to noise ratio of the NMR Thermometer compound has to be at least 400:1 (I-ratio < 250:1).
- If the field lock signal is very large (for example using pure D_2O) the concentration of the NMR Thermometer compound has to be adjusted accordingly (see figure below).

Thermometer substances with a signal closer than 1 ppm to the field lock signal are critical and should not be selected as NMR Thermometer signals.

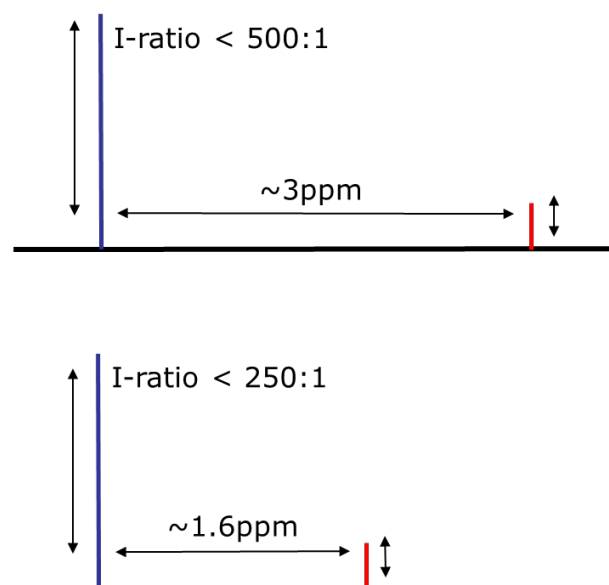


Figure 5.7: Approximate dynamic range of field lock (blue) vs. the NMR Thermometer signal (red) if the S/N of the frequency lock signal is sufficient (> 200:1).

The best performance (with low concentrations of the thermometer substance) can be expected using cryogenically cooled probes at high magnetic field strengths (600 MHz and higher), because the ^2H sensitivity will be the highest. Therefore, check the ^2H sensitivity of your system beforehand. Nevertheless, the NMR Thermometer also works at lower fields (e.g. 400 MHz) with room temperature probes.

The ^2H sensitivity of a room temperature probe, such as an inverse or broad band observe probe, is about a factor 5-10 times lower compared to a cryogenically cooled probe. The concentration of the NMR Thermometer compound therefore needs to be higher, as listed in the following table:

NMR Thermometer compound	Solvent	Chemical shift [ppm]	Minimum Concentration [CryoProbe TCI 5mm 600MHz]	Minimum Concentration [BBFO 5mm 500MHz]
Methanol-d4	methanol-d4	3.3 (CD3) 4.8 (OD)		
Acetic acid-d4	D2O, 5-100% in H2O	2.03	20-50mM	200-500mM
Sodium acetate-d3	D2O, 5-100% in H2O	1.8	20-50mM	200-500mM
DMSO-d6	D2O, 5-100% in H2O	2.49	20mM	200mM
Acetone-d6	D2O, 5-100% in H2O	2.04	20mM	200mM
Acetonitril-d3	D2O, 5-100% in H2O	1.93	20-50mM	200-500mM
Tetramethylammoniumchlorid-d12	D2O, 5-100% in H2O	3.2	10mM	100mM
TRIS-d11	D2O, 5-100% in H2O	3.5	20mM	200mM
TSP-d14	D2O, 5-100% in H2O	0	n.a.	n.a.
DSS-d16	D2O, 5-100% in H2O	0	n.a.	n.a.
Pivalic acid	D2O, 5-100% in H2O	1	10-20mM	100-200mM

Figure 5.8: Typical concentrations required for the NMR Thermometer compound. If the NMR Thermometer compound is used with pure D2O as solvent, the higher concentration of the thermometer compound is required due to the high dynamic range of the intensity of both lock compounds.

5.2.3 Considerations for Shimming with Topshim

Depending on the solvent, topshim will use ^1H or ^2H as a shim nucleus. For aqueous solutions, e.g. 5% D₂O in H₂O, ^1H is the shim nucleus, therefore the shim routine for Topshim does not need to be changed. For deuterated organic solvents ^2H is used as the shim nucleus. If the solvent contains more than one ^2H signal, like methanol-d₄ and pyridine-d₅, Topshim will use a selective ^2H pulse in order not to excite additional signals.

In the case where the thermometer compound is used together with an organic solvent, either in a mixture with the solvent or as an external capillary, the shim routine has to be adapted. In the following the procedure for the setup of the shimming routine for a new solvent is described.

Two steps should be performed:


- Define a new solvent with *edlock*. The procedure is described in the section [Define a New Solvent](#) [19] of this manual.
- Define the shimming routine for Topshim.

The command for defining the Topshim shim parameters is:


topshim solvcal solvent= <new solvent name as in edlock> where <new solvent name as in edlock> is exactly the name of the solvent as defined in the lock/solvent table.

The following example will show the setup for a solvent called *New*.


1. Enter the command **topshim solvcal solvent=New**. A window pops up allowing the selection of a solvent. Here it is of no importance which solvent is selected:

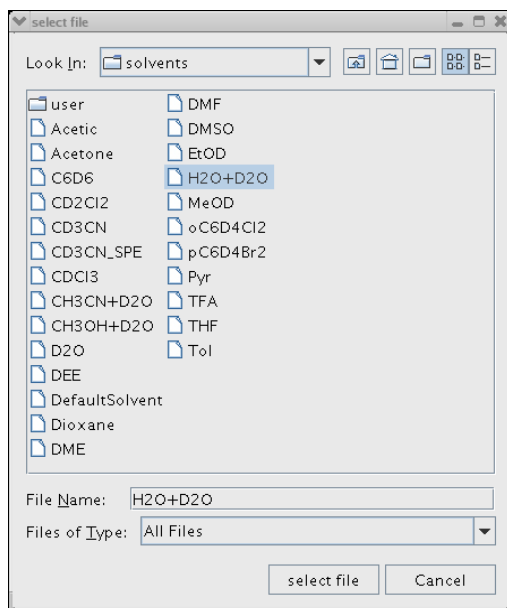


2. Select the solvent of your choice and press the **select file** button. A new window will open, select the **OK** button to modify the shim parameters:

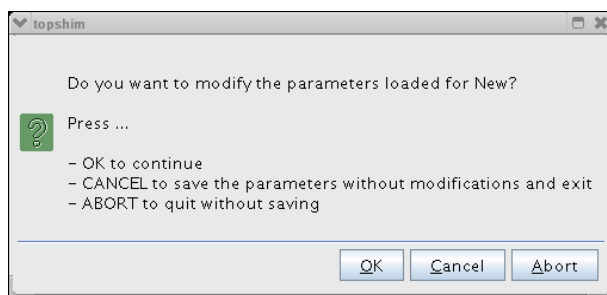


3. A new window will open and allow the selection of the shim nucleus. Enter **2** for ^2H as the shim nucleus:

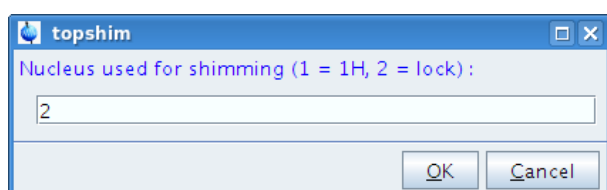


4. Now the frequency of the signal has to be defined. As default, the frequency of the field lock solvent, which is the so-called *lockshift*, will be selected. Enter **1**:

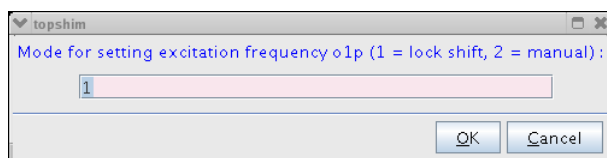


5. Enter **2** to activate the selective excitation:

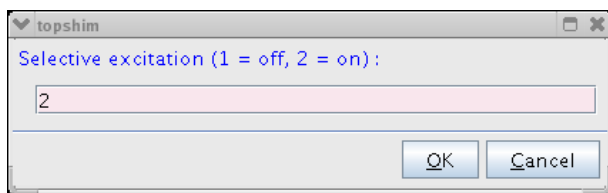

6. The definition of the selectivity for selective excitation is entered next. This depends on the difference of the chemical shift of the field lock solvent to the next closest signal of the thermometer compound. As an example, a selectivity of 0.5 ppm is sufficient for chemical shift difference of 1ppm. For larger shift differences a selectivity of 1 ppm shall be used:

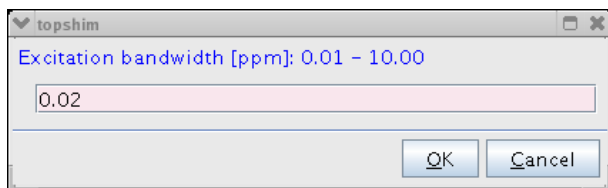


7. The T_1 relaxation time of the shim nucleus is used to define the repetition time for the shim procedure. Typical values for the T_1 relaxation time are about 1 sec. for D_2O , 3 sec for C_6D_6 and 5 sec. for acetone- d_6 .

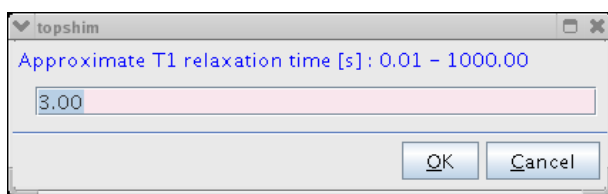

8. In a final step the optimization parameter has to be defined. There are three options:

- **ss**: Shims will be optimized for solvent suppression
- **ls**: Shims will be optimized for the line shape
- **Ishump**: Shims will be optimized for narrow hump.

In the example shown here, solvents suppression has been selected:

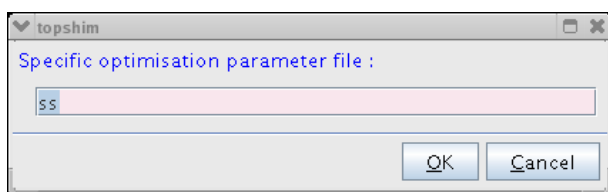

Further details of the setup of a new solvent for shimming can be found in the Topshim reference manual which can be accessed with the command **help topshim** in TopSpin.

6 Applications

6.1 Identical Chemical Shifts at Different Spectrometers

One of the most important applications of the NMR Thermometer is to use it as internal temperature reference and correct the temperature offset between sample and the temperature sensor of the probe. This is in general comparable to the conventional temperature correction (described elsewhere) using the methanol sample (or other temperature calibration samples) derived from two ^1H spectra measured at two different temperatures.

In the case of the NMR Thermometer you just insert the methanol sample (99.8% deuterated) and after setting up the system (tuning/matching/shimming) the sample temperature is displayed immediately. After enabling the NMR Thermometer the sample temperature can be used as target temperature (edte). If the temperature is stable the sample is replaced by, for example, 2 mM sucrose sample in 9:1 $\text{H}_2\text{O}/\text{D}_2\text{O}$ and a ^1H spectrum is acquired. Repeating the same procedure at a second spectrometer leads to a very small shift difference corresponding to a temperature difference of about 40 mK:

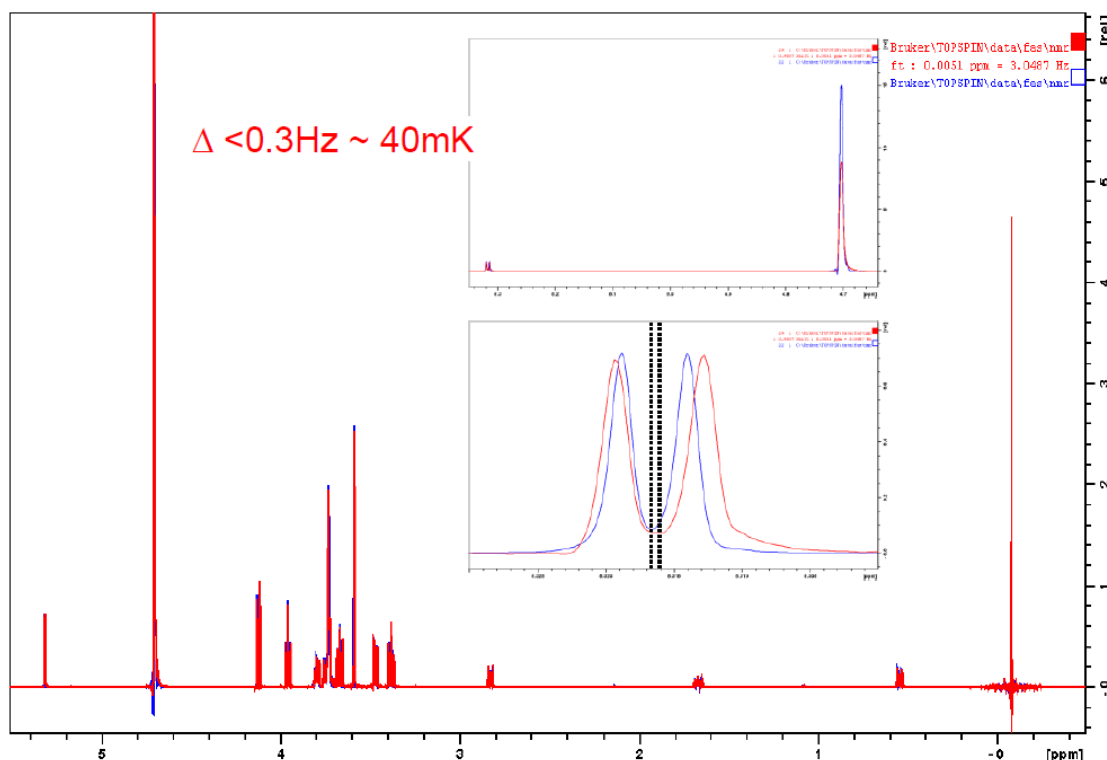


Figure 6.1: Overlay of a ^1H spectrum of 2 mM sucrose (standard sample) measured at 600 MHz (TXI probe) and 800 MHz (TCI CryoProbe).

Changing the sample to a 0.5 mM ubiquitin sample in 9:1 $\text{H}_2\text{O}/\text{D}_2\text{O}$ (figure below) nicely shows the precision of the temperature correction obtained by the NMR Thermometer using the methanol sample (99.8% deuterated) as temperature reference.

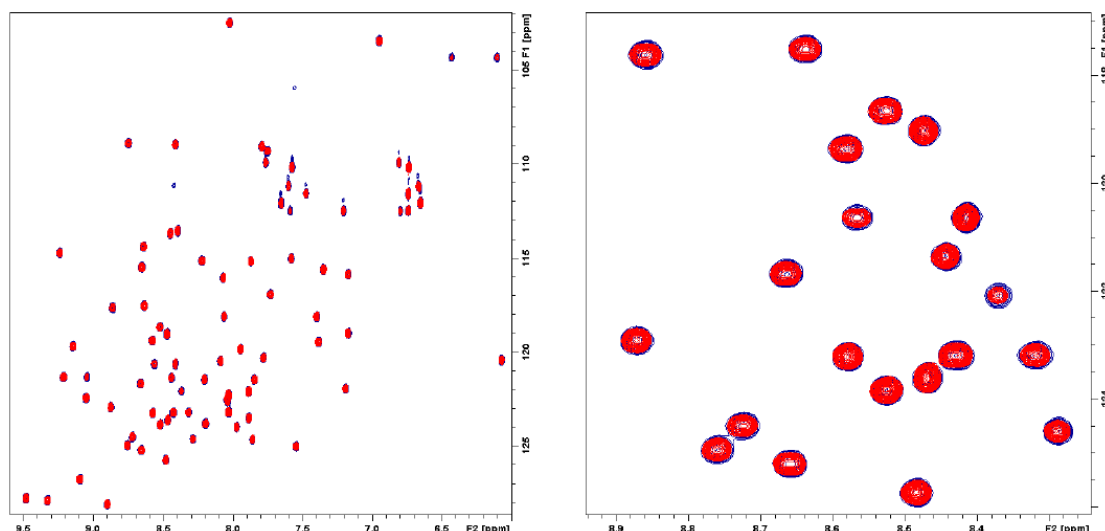


Figure 6.2: Overlay of ^{15}N HSQC spectra (overview: left part) of 0.5 mM ubiquitin 9:1 $\text{H}_2\text{O}/\text{D}_2\text{O}$.

6.2 Identical Chemical Shifts for Experiments with Different Heating due to Experimental Conditions (Temperature Compensation)

The superior feature of the NMR Thermometer is the ability to compensate for sample heating inside the sample due to different sources (RF heating, spinning speed, HR MAS).

6.2.1 RF Heating

RF heating inside the sample is caused by, for instance, a spinlock sequence as used in TOCSY-type of pulse sequences, or decoupling as well as CPMG sequences. Such experiments are widely used in biomolecular NMR. As an example we show a 0.5 mM ^{13}C - and ^{15}N -labeled ubiquitin in $\text{H}_2\text{O}/\text{D}_2\text{O}$ with sodium acetate- d_3 added as NMR Thermometer substance. In the example the 2D HSQC-planes of 3D NOESY-HSQC, TOCSY-HSQC and CPMG-HSQC experiments are compared. Reference planes were measured without temperature compensation (figure below, left part) and the others with NMR Thermometer in the regulation mode (figure below, right part).

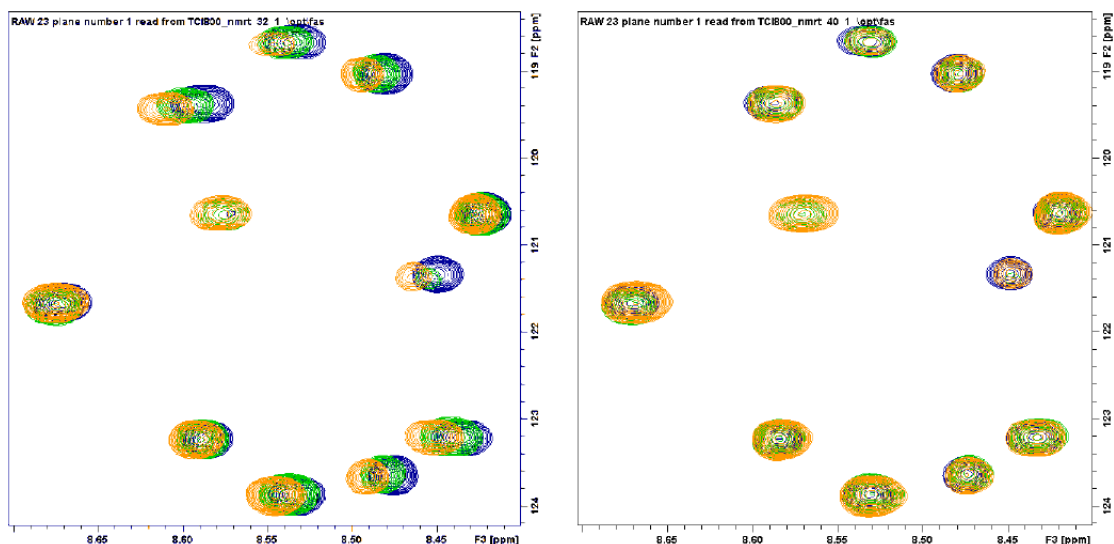


Figure 6.3: Overlay of NOESY-HSQC (blue), TOCSY-HSQC (orange) and CPMG-HSQC (green) spectra of 0.5mM ubiquitin in 95:5% H₂O/D₂O measured at 800MHz TCI CP.

In the figure above the left side shows the NMR Thermometer disabled (monitoring mode) and the right side the NMR Thermometer enabled (regulation mode).

6.2.2 Heating caused by Spinning Speed (HR-MAS)

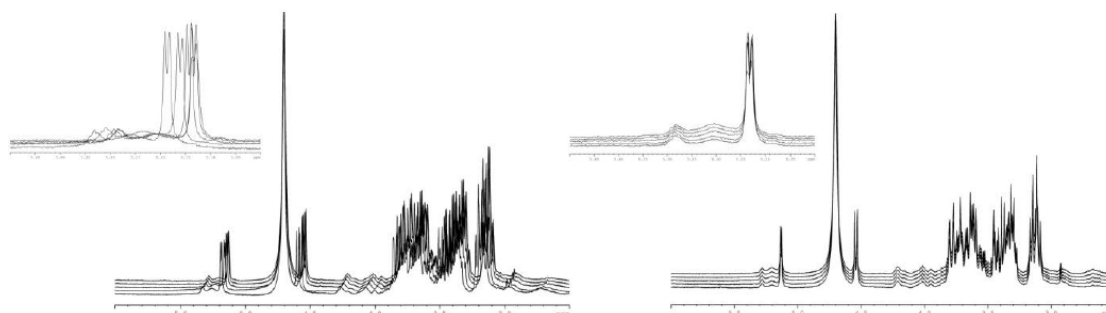


Figure 6.4: ¹H NMR spectra on a liver sample with sodium acetate added.

In the figure above, different spinning speeds were used (1, 2, 4, 6 and 8 kHz). Left part: NMR Thermometer disabled (monitoring mode); Right part: NMR Thermometer enabled (regulation mode).

An interesting application of the NMR Thermometer is high resolution MAS on biological material. Depending on the spinning speed used in HR-MAS (1 to 8 kHz) the frictional heating in the rotor is different and hence the temperature varies. To compensate for the heating is not only of interest for the comparability of the spectra, but it is also important in order to preserve sensitive sample material (like tissue material). The temperature difference between sample and probe sensor can be as much as 5K in the case of 8 kHz spinning speed.

Applying a spin lock sequence (TOCSY) on top of that could increase the temperature even further. Both heating effects can be compensated for by the NMR Thermometer:

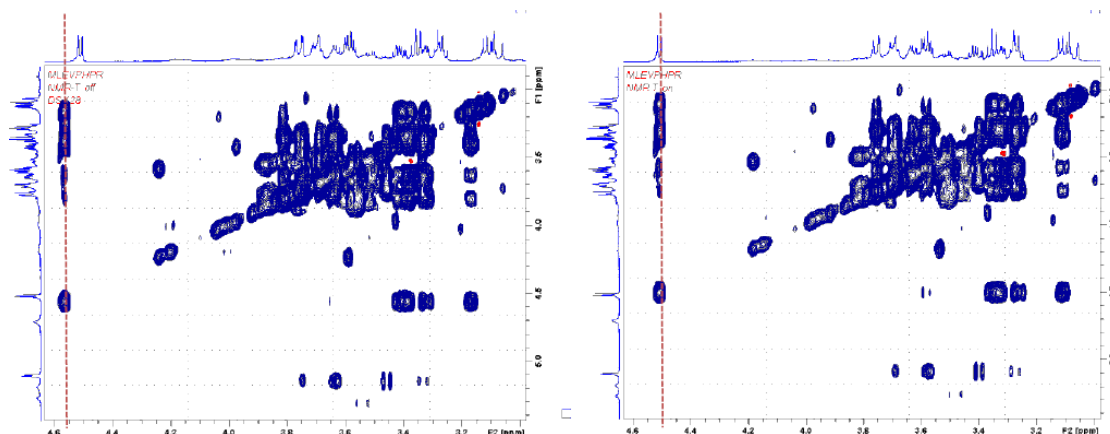


Figure 6.5: Comparison of TOCSY experiments of a liver sample with sodium acetate added measured with HR-MAS at 4 kHz spinning speed.

In the figure above is a comparison of TOCSY experiments of a liver sample with sodium acetate added measured with HR-MAS at 4 kHz spinning speed. The reference ^1H spectrum is plotted as projection. Left part: NMR Thermometer disabled (monitoring mode); Right part: NMR Thermometer enabled (regulation mode).

7 Frequently Asked Questions (FAQ)

The lock procedure works well with the solvent, but not with the frequency lock.

1. *The wrong solvent has been selected and thus the frequency of the compound used for the NMR Thermometer is out of range.*
2. *The concentration of the compound used for the NMR Thermometer is too low. Please check table 1 and 2 which give typical concentrations required.*
3. *The lock power for the thermometer compound is too high or too low. Change the value "Temperature Lock power" in the edlock table and repeat the lock procedure. The value of the temperature lock power typically is a few dB lower than for the field lock power.*
4. *The chemical shift of the compound used for the NMR Thermometer is too closed to the chemical shift of the solvent. Here it can happen that the lock procedure for both, the field and the frequency lock, is working fine on a 700 MHz spectrometer, while it fails on a lower field spectrometer due to the reduced shift difference (in Hz) at a lower field. In that case a different thermometer compound has to be used.*

The lock procedure worked well for field and frequency lock, but the temperature regulation of the NMR Thermometer is instable.

- *The lock power for the thermometer compound is too high. Reduce the value "Temperature Lock power" in the edlock table and repeat the lock procedure. The value for the temperature power typically is a few dB lower than for the field lock power.*

Does the NMR Thermometer work when experiments with pulsed field gradients are performed?

- *Yes, like the field lock the NMR Thermometer can be used together with pulsed field gradients. Both, the field and the frequency lock regulation are triggered with lock hold commands of the pulse program. Lock hold commands are standard in all standard pulse programs using pulsed field gradients.*

After enabling the NMR Thermometer the temperature starts to increase/decrease or during an experiment with RF heating (see above) temperature is decreasing?

- *The chemical shift – temperature values in the edlock table for the particular solvent are probably wrong for instance after creating an .xml file and using a positive sign for the slope instead of a negative one.*

8 Contact

Manufacturer:

Bruker BioSpin NMR
Silberstreifen
D-76287 Rheinstetten
Germany
Phone: +49 721-5161-6155
<http://www.bruker.com>
WEEE DE43181702

NMR Hotlines

Contact our NMR service centers.

Bruker BioSpin NMR provide dedicated hotlines and service centers, so that our specialists can respond as quickly as possible to all your service requests, applications questions, software or technical needs.

Please select the NMR service center or hotline you wish to contact from our list available at:

<http://www.bruker.com/service/information-communication/helpdesk.html>

List of Figures

Figure 2.1	Principle of the NMR Thermometer.....	7
Figure 4.1	The Submenu Bar: Solvent.....	11
Figure 4.2	Edit Solvent Parameters.....	12
Figure 4.3	The Submenu Lock.....	12
Figure 4.4	Edit Lock Parameters.....	13
Figure 4.5	Edte Window: NMR Thermometer monitoring mode (disabled, top), regulation mode (enabled, middle) and the selection of both modes in the Configuration menu of the edte window (bottom).....	14
Figure 4.6	Selection of the Auto Phase Algorithm.....	14
Figure 4.7	Selftune warning about PID parameters misfit.....	14
Figure 4.8	Starting the selftune procedure from the edte window.....	15
Figure 4.9	Selftune in Progress.....	16
Figure 4.10	Monitoring several BSMS values during a temperature jump.....	17
Figure 4.11	TopSpin status bar with NMR Thermometer enabled.....	17
Figure 4.12	Monitoring of the sensor and the sample temperature in edte during a TOCSY experiment.....	18
Figure 5.1	Adding New Solvents.....	19
Figure 5.2	Assign Signals to Type of Signal (Signal, Lock, Temperature).....	20
Figure 5.3	Dialog for importing temperature shift value files.....	20
Figure 5.4	Example of a shift-temperature file in .xml format.....	20
Figure 5.5	Temperature shift values filled in.....	21
Figure 5.6	Suggested compounds for the NMR Thermometer and samples in aqueous solution (D2O 5 – 100%), their properties and estimated concentration. Fully deuterated DSS and TSP are currently not available but will be synthesized on Bruker's request at small amounts only for internal tests.....	21
Figure 5.7	Approximate dynamic range of field lock (blue) vs. the NMR Thermometer signal (red) if the S/N of the frequency lock signal is sufficient (> 200:1).....	23
Figure 5.8	Typical concentrations required for the NMR Thermometer compound. If the NMR Thermometer compound is used with pure D2O as solvent, the higher concentration of the thermometer compound is required due to the high dynamic range of the intensity of both lock compounds.....	24
Figure 6.1	Overlay of a 1H spectrum of 2 mM sucrose (standard sample) measured at 600 MHz (TXI probe) and 800 MHz (TCI CryoProbe).....	27
Figure 6.2	Overlay of 15N HSQC spectra (overview: left part) of 0.5 mM ubiquitin 9:1 H2O/D2O.....	28
Figure 6.3	Overlay of NOESY-HSQC (blue), TOCSY-HSQC (orange) and CPMG-HSQC (green) spectra of 0.5mM ubiquitin in 95:5% H2O/D2O measured at 800MHz TCI CP.....	29
Figure 6.4	1H NMR spectra on a liver sample with sodium acetate added.....	29
Figure 6.5	Comparison of TOCSY experiments of a liver sample with sodium acetate added measured with HR-MAS at 4 kHz spinning speed.....	30

Index

A

Absolute temperature correction	21
Aqueous solutions	21
autophase	14

B

Bio-NMR	21
Buffers	22

C

Chemical shift distance	23
Chemical shift of water	21
Chemical shifts	5
Compounds	22

D

Digilock 2G	9
Dynamic range	23

E

edlock	11, 20
edsolv	11
exercise	16

I

import shift and temperature values	20
---	----

L

Lock parameters	12
lock phase	14
LTRX	7

M

Monitoring	16
------------------	----

N

New solvent	19
NMR Thermometer compounds	21

O

Overview	11
----------------	----

P

PID values	14
Predefined Solvents	22
Principle	7
pulsed field gradients	31

R

RF heating	5
------------------	---

S

Salt effect	22
Selecting a suitable compound	22
selftune	14, 15
Slope estimation	21
SmartVT	9
Stabilizers	22

T

Thermometer substance	7
Topshim	13, 24

V

VT control	11
------------------	----

X

xml files	20
-----------------	----





Bruker Corporation

info@bruker.com
www.bruker.com

Order No: Z33085